

Cloud Setup Assistant



Create a cloud account at
thermofisher.com/connect



ThermoFisher
SCIENTIFIC



Turn on your instrument,
click **“Get Started,”**
and then click **“Connect”**



ThermoFisher
SCIENTIFIC

Is your instrument
connected to the internet?

YES



NO

Click one of the options above to continue



ThermoFisher
SCIENTIFIC



Your device is connected to the internet

1

Select "Instrument"

2

Enter your Thermo Fisher Connect
username and password

3


Select "link account"



ThermoFisher
SCIENTIFIC



Your device is not connected to the internet

- 1** Go to thermofisher.com/connect on your computer and enter your username and password
 - 2** On your cloud dashboard, select the instrument icon on the left sidebar
 - 3** Select "Add Instrument" in the upper-right corner
 - 4** Select your instrument from the drop-down list and click "Next"
 - 5** On your instrument screen, select "PC" to connect. A linking code will be displayed.
 - 6** Enter the linking code into your cloud dashboard on your computer.
- 



ThermoFisher
SCIENTIFIC

Your instrument is now connected

You're now ready to send instrument data to the platform using integrated data analysis apps and collaboration tools.



Want to restart the helper?

Click the icon above to restart from the beginning, or click on one of the steps below to jump to a specific place.



Having installation issues?

[Contact customer support for assistance](#)



ThermoFisher
SCIENTIFIC