



AppliedBiosystems™ RapidLINK™ Software v1.1.5

RapidLINK upgrade installer workflow


1. Download the RapidLINK installer and save it on S:\ drive of the laptop where the earlier version of RapidLINK is already installed
2. After successful download, close the RapidLINK application
3. Navigate to the S:\ drive where the 'RapidLINKsetup.exe' is saved and right click on the .exe to run it as an Administrator
4. Click on 'Yes' to continue with Installation

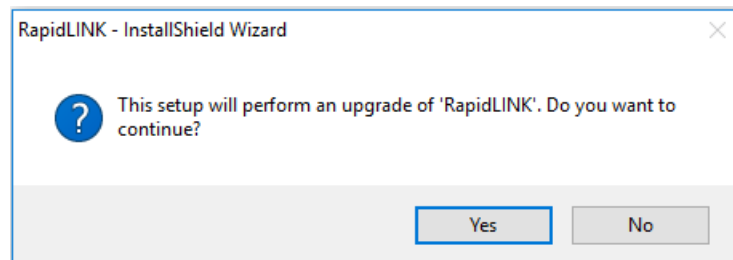


5. If promoted to Accept the EULA, select the 'I accept the terms in the license agreement' option and click 'Next'

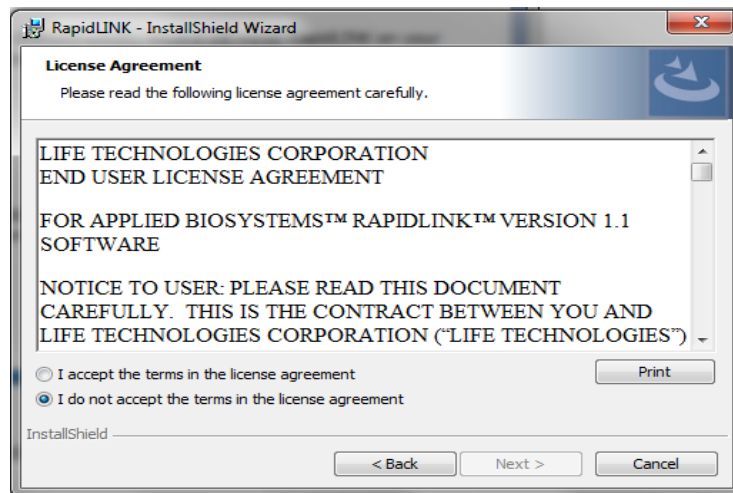

- Click on 'Next' to proceed with the install

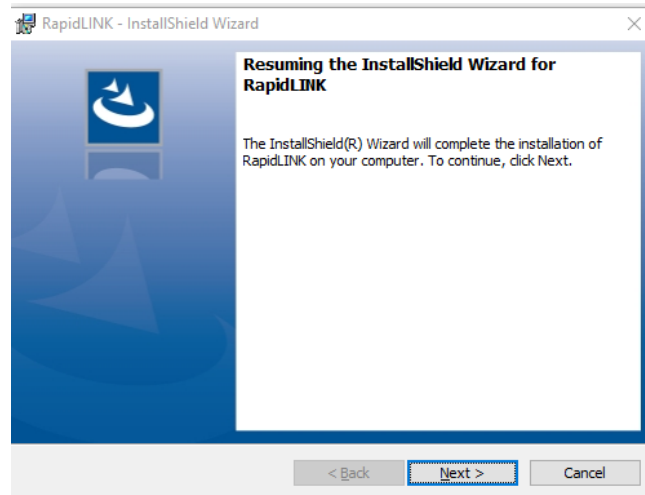


- Click on 'Finish' when installation is complete

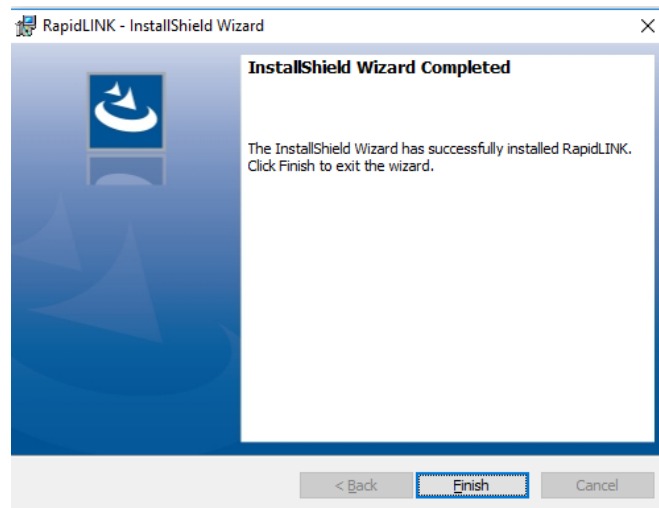

- Launch RapidLINK from the desktop shortcut
RapidLINK can also be launched by clicking on RapidLINK.exe from C:\RapidLINK

Steps to verify successful upgrade

1. Launch RapidLINK, click on Settings icon on the left side of the screen.
2. The build version should be 1.1.5

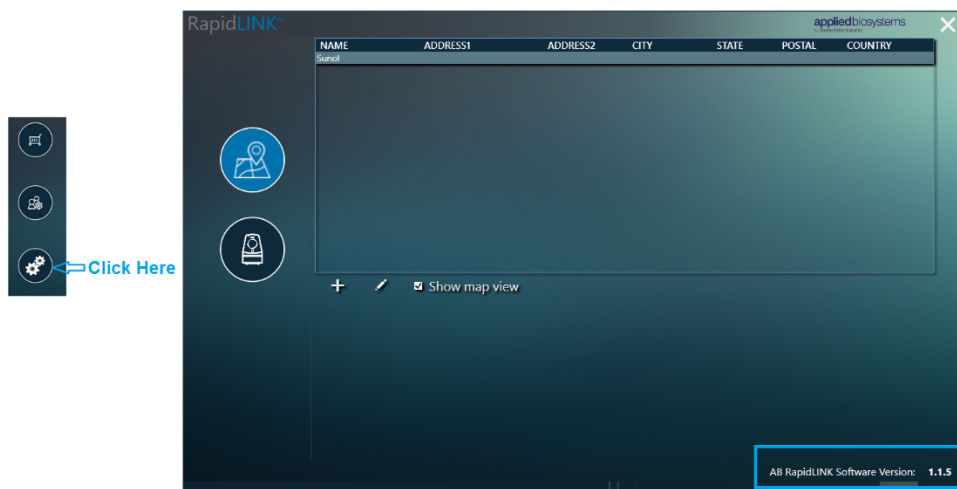

LEGAL STATEMENTS

For Forensics, Human Identification or Paternity/Kinship Use Only. Not for use in diagnostic or therapeutic applications

NOTICE TO PURCHASE: DISCLAIMER OF LICENSE

Purchase of this software product alone does not imply any license under any process, instrument or other apparatus, system, composition, reagent or kit rights under patent claims owned or otherwise controlled by Thermo Fisher Scientific, either expressly, or by estoppel.

DISCLAIMER

LIFE TECHNOLOGIES CORPORATION, PART OF THERMO FISHER SCIENTIFIC INC., AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES, PART OF THERMO FISHER SCIENTIFIC INC., AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

The trademarks mentioned herein are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

© 2019 Thermo Fisher Scientific. All rights reserved.