

# Protocol for Sanger sequencing of SARS-CoV-2 B.1.1.7 and B.1.351 strain lineages

The global emergency of SARS-CoV-2 has caused significant disruptions to economic and personal activities. Recently, two new lineages that appear to have increased infectivity have been identified. The B.1.1.7 lineage, first identified in England and subsequently in the US, contains 23 mutations scattered across the SARS-CoV-2 genome [1]. The other lineage, B.1.351, first identified in South Africa, contains a set of mutations clustering in the SARS-CoV-2 spike (S) gene [2]. Since both lineages are thought to have increased transmissibility relative to other lineages, it is important to determine whether samples contain the mutations identified in the two lineages [2,3].

We developed a Research Use Only protocol for detecting these lineages by Sanger sequencing. The primer sequences used here are based on those published by the Centers for Disease Control and Prevention (CDC) [4]. Briefly, cDNA synthesis is performed on a sample containing viral RNA. Next, the cDNA is used in specific regions of target amplification using primers that have been selected to cover the novel mutations. For this, the Applied Biosystems™ BigDye™ Direct Cycle Sequencing Kit and M13 sequence–tagged primer sets are used. The amplified sequences are then subjected

to cycle sequencing using either M13-forward or M13-reverse primers provided in the BigDye Direct Cycle Sequencing Kit. Unincorporated nucleotides and primers are next removed using the Applied Biosystems™ BigDye XTerminator™ Purification Kit, and the sequences are read by standard capillary electrophoresis (CE). The sequences obtained can be read by any sequencing program, such as SeqA or Geneious™ software, and compared with known or expected SARS-CoV-2 sequences (Figure 1).

Some of the sequences generated by this method will produce CE traces that may be difficult to interpret. To determine whether a sequencing trace was useful, we employed quality control metrics generated by Applied Biosystems™ Sequence Scanner Software v2.0. These metrics include trace score (average of basecaller quality values for bases in the clear range), contiguous read length (CRL), and QV20+ (total number of bases in the entire trace that have a basecaller quality value of ≥20). Guidelines for using these metrics for QC and analysis of results are given at the end of the protocol. However, standard analysis of sequencing traces is often sufficient to determine whether a novel sequence is present.



**Figure 1. Workflow for detection of SARS-CoV-2 lineages using Sanger sequencing.** RNA is purified from samples using standard techniques. cDNA is synthesized from the RNA, and specific M13 sequence–tagged amplicons are generated by PCR. The amplicons are sequenced in the forward and reverse directions using universal M13 primers and the BigDye Direct Cycle Sequencing Kit. The sequencing reactions are cleaned using the BigDye XTerminator kit and subjected to CE. The resulting sequencing traces can be analyzed and compared to reference SARS-CoV-2 sequences to determine if the lineages are present.

**This protocol and the reagents described within are for Research Use Only. Not for use in diagnostic procedures.**

**IMPORTANT:** This protocol is very sensitive; therefore, utmost care must be taken to prepare the stock solutions and set up the amplification reactions in an amplicon-free environment.

## 1. Materials needed

### 1.1 Equipment

Product	Supplier	Cat. No.
Veriti 96-Well Fast Thermal Cycler, ProFlex 96-Well PCR System, or similar thermal cycler	Thermo Fisher Scientific	4375305 or 4484075
MicroMixer E-36 for 96-well plates	Taitec	0027765-000
Single-channel and multichannel micropipettes of various sizes capable of pipetting volumes from 1.00 $\mu$ L to 1,000.0 $\mu$ L	MLS	Any
Cold block or ice	MLS	Any
Plate centrifuge	MLS	Any
Microcentrifuge or mini centrifuge	MLS	Any
Vortex mixer	MLS	Any

### 1.2 Reagents, kits, and consumables

Product	Supplier	Cat. No.
SuperScript™ IV VILO™ Master Mix	Thermo Fisher Scientific	117565500
Nuclease-Free Water	Thermo Fisher Scientific	AM9937 or equivalent
BigDye Direct Cycle Sequencing Kit	Thermo Fisher Scientific	4458688 or equivalent
BigDye XTerminator Purification Kit	Thermo Fisher Scientific	4376486 or equivalent
MicroAmp Fast Optical 96-Well Reaction Plate with Barcode, 0.1 mL	Thermo Fisher Scientific	4346906 or 4366932
MicroAmp Optical Adhesive Film	Thermo Fisher Scientific	4311971, 4313663, or 4360954
Nonstick, RNase-Free Microcentrifuge Tubes, 1.5 mL	Thermo Fisher Scientific	AM12450 or equivalent
5 mL tube, PCR clean	MLS	Any
Sterilized aerosol barrier (filter) pipette tips	MLS	Any

### 1.3 Primers

- Primer sequences are given in Table 1.

**Table 1. Sequences of M13-tagged primers for detecting mutations found in the lineage B.1.1.7.** A subset of primer pairs that focus on specific regions of the SARS-CoV-2 genome can be chosen according to researchers' needs; a complete list that covers all regions mutated in the B.1.1.7 and South African lineages is provided here. Mutations that are synonymous and produce no amino acid change have a :syn suffix (e.g., C15279T:syn). Mutations that are associated with the South African B.1.351 lineage are indicated with (SA). Mutations that are common to both lineages are indicated with (common). The M13 sequence tags are highlighted in red.

Mutation covered by primer pair	Forward primer name	Forward primer sequence	Reverse primer name	Reverse primer sequence
G174T:nc (SA)	SC2M1-1_ LEFT_M13	TGTAAACGACGGCCAGT TTTGCCTCAACTTGAACAGCCC	SC2M1-1_ LEFT_M13	CAGGAAACAGCTATGACCTGTCTCACCCTAGGACCGTAC
C913T C1059T:T265I (SA)	SC2M1-3_ LEFT_M13	TGTAAACGACGGCCAGT AACAACTTCTGTGCCCTGATG	SC2M1-3_ RIGHT_M13	CAGGAAACAGCTATGACCTCTGAATTGTGACATGCTGGACA
C3267T:T1001I	SC2M1-8_ LEFT_M13	TGTAAACGACGGCCAGT ACTTACACCCTGGCATTGATT	SC2M1-8_ RIGHT_M13	CAGGAAACAGCTATGACCTGCAACACCTCCTCCATGTTT
G5320T:K1655N (SA)	SC2M1-13_ LEFT_M13	TGTAAACGACGGCCAGT TCCTACCACATCCACCTAGATGG	SC2M1-13_ RIGHT_M13	CAGGAAACAGCTATGACCGCACAAAAGTTAGCAGCTTCACC
C5388A:A1708D	SC2M1-14_ LEFT_M13	TGTAAACGACGGCCAGT ACTTCTATTAATGGGAGATAACAAGTGT	SC2M1-14_ RIGHT_M13	CAGGAAACAGCTATGACCGCAGCACCCTCTGCAATACAAAGT
T6954C:I2230T	SC2M1-18_ LEFT_M13	TGTAAACGACGGCCAGT AACCGTGTTTGTACTAATATATGCCTT	SC2M1-18_ RIGHT_M13	CAGGAAACAGCTATGACCTGCCAAAACCACTCTGCAACT
A10323G:K3353R (SA)	SC2M1-26_ LEFT_M13	TGTAAACGACGGCCAGT AGTACAAGTATTTAGTGGAGCAATGGA	SC2M1-26_ RIGHT_M13	CAGGAAACAGCTATGACCTGGGCCTCATAGCACATTGGTA
del11288-11296i: delSGF3675-3677 (common)	SC2M1-29_ LEFT_M13	TGTAAACGACGGCCAGT AGTCCAGAGTACTCAATGGTCTTTGT	SC2M1-29_ RIGHT_M13	CAGGAAACAGCTATGACCAATACCTCTGGCCAAAACATGA
C14408T:P4715L (common)	SC2M1-37_ LEFT_M13	TGTAAACGACGGCCAGT TTTTCATACAAACCACGCCAGGT	SC2M1-37_ RIGHT_M13	CAGGAAACAGCTATGACCGTGCAGCTACTGAAAAGCACGT
C15279T:syn	SC2M1-39_ LEFT_M13	TGTAAACGACGGCCAGT ACGATGGTGGCTGTATTAATGCT	SC2M1-39_ RIGHT_M13	CAGGAAACAGCTATGACCGGTGTGACAAGCTACAACACGT
T16176C:syn	SC2M1-42_ LEFT_M13	TGTAAACGACGGCCAGT GGAGTATGCTGATGCTTTCATTGTAC	SC2M1-42_ RIGHT_M13	CAGGAAACAGCTATGACCGCGTTTCTGCTGCAAAAAGCTT
del21765-21770: delHV69-70 A21801C_S:D80A (SA)	SC2M1-55_ LEFT_M13	TGTAAACGACGGCCAGT AGGGGTACTGCTGTTATGCTTTAAA	SC2M1-55_ RIGHT_M13	CAGGAAACAGCTATGACCAAGTAGGGACTGGGTCTTCGAA
del21991-21993:delY144 A22206G_S:D215G (SA) del22281-22289: del241/243 (SA)	SC2M1-56_ LEFT_M13	TGTAAACGACGGCCAGT TGGGACCAATGGTACTAAGAGGT	SC2M1-56_ RIGHT_M13	CAGGAAACAGCTATGACCAACAGCTGTCCAACCTGAAGAA
G22813T:K417N (SA)	SC2M1-58_ LEFT_M13	TGTAAACGACGGCCAGT ACTTGTGCCCTTTGGTGAAGT	SC2M1-58_ RIGHT_M13	CAGGAAACAGCTATGACCTGCTGGTGCATGTAGAAGTTCA
C23271A:A570D G23012A_S:E484K (SA) A23063T_S:N501Y (common) AA23403GC:D614G (SA) C23604A:P681H C23709T:T716I C23664T_S:A701V (SA)	SC2M1-59_ LEFT_M13	TGTAAACGACGGCCAGT CCGGTAGCACACCTTGTAAATGG	SC2M1-59_ RIGHT_M13	CAGGAAACAGCTATGACCCCTATTAAACAGCCTGCACG
G24914C:D1118H	SC2M1-60_ LEFT_M13	TGTAAACGACGGCCAGT ACCAGTGTGCTTCTTATCAGG	SC2M1-60_ RIGHT_M13	CAGGAAACAGCTATGACCGAGTATTCCAGTTAAAGCACGGT
G625563TC:Q57H (SA)	SC2M1-63_ LEFT_M13	TGTAAACGACGGCCAGT AAATGATATCCTTTCAGCTTGACAAA	SC2M1-63_ RIGHT_M13	CAGGAAACAGCTATGACCTGAGTCTAATTGAGTTGCAAGGA
G625563TC:Q57H (SA)	SC2M1-65_ LEFT_M13	TGTAAACGACGGCCAGT CTAGGTTTTATAGCTGGCTTGATTGC	SC2M1-65_ RIGHT_M13	CAGGAAACAGCTATGACCCATTTCCAGCAAAGCCAAAGCC
C25904T:S171L (SA)	SC2M1-66_ LEFT_M13	TGTAAACGACGGCCAGT CTCACACCTTTTGTCTGTTGCT	SC2M1-66_ RIGHT_M13	CAGGAAACAGCTATGACCGTGTCTACAAAGGCACGCTAGT
C26456T:T205I (SA)	SC2M1-67b_ LEFT_M13	TGTAAACGACGGCCAGT TCACACAATCGAGGTTTATCC	SC2M1-67b_ RIGHT_M13	CAGGAAACAGCTATGACCGTACCCTTGGAAATCTGCCATGG
C27972T:Q27stop G28048T:R521 A28111G:Y73C	SC2M1-71_ LEFT_M13	TGTAAACGACGGCCAGT TGTTTCATCAGACAAGAGGAAGTTCA	SC2M1-71_ RIGHT_M13	CAGGAAACAGCTATGACCGCAACAACGCACTACAAGACT
GATG28280CTAT:D3L	SC2M1-72b_ LEFT_M13	TGTAAACGACGGCCAGT GTAGGAGCTAGAAAATCAGCACC	SC2M1-72b_ RIGHT_M13	CAGGAAACAGCTATGACCCATATGATGCCGCTTTGTTAGC
RG203-204KR C28977T:S235F	SC2M1-73_ LEFT_M13	TGTAAACGACGGCCAGT TGGCTACTACCGAAGAGCTACC	SC2M1-73_ RIGHT_M13	CAGGAAACAGCTATGACCGCTTCTAGAAGCCTCAGCAGC

- Primers can be ordered from our **custom oligo ordering web page** (<https://www.thermofisher.com/order/custom-standard-oligo>).

– 25 nmol of dried and desalted primers can be ordered, but order can be scaled up as needed.

- Resuspend dried oligos to final concentration of 100  $\mu$ M with TE.

#### 1.4. Amplification mixes of primers

- Prepare the target-specific amplification primer mixes:

– Label clean microcentrifuge tubes for each primer pair (e.g., SC2M1-3). Add 492  $\mu$ L TE to each tube.

– Add 4  $\mu$ L each of both the left and right oligos of a pair to the appropriate tube (i.e., SC2M1-3\_LEFT\_M13 and SC2M1-3\_RIGHT\_M13 in one tube, SC2M1-8\_LEFT\_M13 and SC2M1-8\_RIGHT\_M13 to the next, etc.).

– These will be the 10X sequencing amplification primer mixes, with each oligo at 0.8  $\mu$ M, that will be used in step 3.1–3.2.

## 2. cDNA synthesis

2.1. For each sample, combine:

Reagent	Final volume
	50 $\mu$ L
5X SuperScript IV VILO Master Mix	10 $\mu$ L
Sample	1–15 $\mu$ L
Water	To final 50 $\mu$ L

2.2. Vortex for 2–3 seconds, then centrifuge briefly (5–10 seconds) at 1,000 x *g*.

**Note:** Sample input volume can be adjusted for sensitivity. For example, up to 15  $\mu$ L of a sample that is expected to have low titer may be used.

2.3. Reverse transcription

2.3.1. Program a thermal cycler with the following profile:

Parameter	Stage/step			
	Annealing	Polymerase extension	Polymerase inactivation	Hold
Temperature	25°C	50°C	80°C	4°C
Time	10 min	15 min	10 min	Indefinitely

2.3.2. Put samples in the thermal cycler and run the program.

**Note:** Samples can be held at 4°C or on ice for up to 8 hours; for longer storage, freeze at –20°C.

### 3. PCR amplifications of targets

3.1. For each sample, a forward and reverse reaction will be run. The initial PCR amplification, therefore, requires two identical reactions to be set up. An example 96-well plate setup for two samples is shown below:

	1	2	3	4	5	6	7	8	9	10	11	12
A	SC2M1-3 primers	SC2M1-3 primers	SC2M1-42 primers	SC2M1-42 primers	SC2M1-71 primers	SC2M1-71 primers	SC2M1-3 primers	SC2M1-3 primers	SC2M1-42 primers	SC2M1-42 primers	SC2M1-71 primers	SC2M1-71 primers
B	SC2M1-8 primers	SC2M1-8 primers	SC2M1-55 primers	SC2M1-55 primers	SC2M1-72 primers	SC2M1-72 primers	SC2M1-8 primers	SC2M1-8 primers	SC2M1-55 primers	SC2M1-55 primers	SC2M1-72 primers	SC2M1-72 primers
C	SC2M1-14 primers	SC2M1-14 primers	SC2M1-56 primers	SC2M1-56 primers	SC2M1-74 primers	SC2M1-74 primers	SC2M1-14 primers	SC2M1-14 primers	SC2M1-56 primers	SC2M1-56 primers	SC2M1-74 primers	SC2M1-74 primers
D	SC2M1-15 primers	SC2M1-15 primers	SC2M1-59 primers	SC2M1-59 primers	No primers	No primers	SC2M1-15 primers	SC2M1-15 primers	SC2M1-59 primers	SC2M1-59 primers	No primers	No primers
E	SC2M1-18 primers	SC2M1-18 primers	SC2M1-60 primers	SC2M1-60 primers	NTC	NTC	SC2M1-18 primers	SC2M1-18 primers	SC2M1-60 primers	SC2M1-60 primers	NTC	NTC
F	SC2M1-29 primers	SC2M1-29 primers	SC2M1-62b primers	SC2M1-62b primers	Positive control	Positive control	SC2M1-29 primers	SC2M1-29 primers	SC2M1-62b primers	SC2M1-62b primers	Positive control	Positive control
G	SC2M1-38 primers	SC2M1-38 primers	SC2M1-64 primers	SC2M1-64 primers	Blank	Blank	SC2M1-38 primers	SC2M1-38 primers	SC2M1-64 primers	SC2M1-64 primers	Blank	Blank
H	SC2M1-39 primers	SC2M1-39 primers	SC2M1-68 primers	SC2M1-68 primers	Blank	Blank	SC2M1-39 primers	SC2M1-39 primers	SC2M1-68 primers	SC2M1-68 primers	Blank	Blank

**Note:** Reactions using the same cDNA sample have identical color coding.

**Note:** The layout above is for querying all sites known to be mutated in the B.1.1.7 strain. If only a subset of amplicons is to be analyzed, the layout can be adjusted accordingly.

**Note:** Wells reserved for controls are suggested in the layout. For the “No primers” control, add sample but do not add amplification primers. For the no-template control (NTC), perform a cDNA synthesis with no added sample, and amplify using any primer pair. For the positive control, use a sample that is known to have SARS-CoV-2 present, and any primer pair.

3.2. In each well of a 96-well PCR plate, combine:

- 1.5 µL of 10X sequencing amplification primer mix in duplicate (as suggested in the table above)
- 5 µL of 2X BigDye Direct PCR Master Mix (supplied in kit)
- 1 µL of cDNA sample from completed step 2.3
  - Leftover cDNA sample can be frozen at –20°C.
- Water to 10 µL total volume

3.3. Seal the plate; vortex for 2–3 seconds, then centrifuge briefly (5–10 seconds) at 1,000 x g.

3.4. Place the plate into a thermal cycler and run the following program:

Parameter	Stage/step				
	Polymerase activation	Cycling (40 cycles)			Hold
		Denaturation	Annealing	Extension	
Temperature	95°C	96°C	62°C	68°C	4°C
Time	10 min	3 sec	15 sec	30 sec	Indefinitely

**Note:** Samples can be held at 4°C or on ice for up to 8 hours; for longer storage, freeze at –20°C.

**Note:** Do not exceed 10 µL total reaction volume. The cycle sequencing and BigDye Xterminator steps have been optimized for 10 µL input volumes.

#### 4. Cycle sequencing

4.1. Once the PCR in step 3.4 is complete, the plate can be used directly for cycle sequencing.

4.2. Remove the seal from the plate.

4.3. To each well of the plate, add:

- 2  $\mu$ L of BigDye Direct Sequencing Master Mix (supplied in kit)
- 1  $\mu$ L of BigDye Direct M13 Forward or M13 Reverse primer (supplied in kit)

**Note:** It is important to add the M13 Forward primer to one of the duplicate PCR reactions, and the M13 Reverse primer to the other reaction. An example based on the plate setup is shown below.

	1	2	3	4	5	6	7	8	9	10	11	12
A	SC2M1-3 M13 For	SC2M1-3 M13 Rev	SC2M1-42 M13 For	SC2M1-42 M13 Rev	SC2M1-71 M13 For	SC2M1-71 M13 Rev	SC2M1-3 M13 For	SC2M1-3 M13 Rev	SC2M1-42 M13 For	SC2M1-42 M13 Rev	SC2M1-71 M13 For	SC2M1-71 M13 Rev
B	SC2M1-8 M13 For	SC2M1-8 M13 Rev	SC2M1-55 M13 For	SC2M1-55 M13 Rev	SC2M1-72 M13 For	SC2M1-72 M13 Rev	SC2M1-8 M13 For	SC2M1-8 M13 Rev	SC2M1-55 M13 For	SC2M1-55 M13 Rev	SC2M1-72 M13 For	SC2M1-72 M13 Rev
C	SC2M1-14 M13 For	SC2M1-14 M13 Rev	SC2M1-56 M13 For	SC2M1-56 M13 Rev	SC2M1-74 M13 For	SC2M1-74 M13 Rev	SC2M1-14 M13 For	SC2M1-14 M13 Rev	SC2M1-56 M13 For	SC2M1-56 M13 Rev	SC2M1-74 M13 For	SC2M1-74 M13 Rev
D	SC2M1-15 M13 For	SC2M1-15 M13 Rev	SC2M1-59 M13 For	SC2M1-59 M13 Rev	No primers M13 For	No primers M13 Rev	SC2M1-15 M13 For	SC2M1-15 M13 Rev	SC2M1-59 M13 For	SC2M1-59 M13 Rev	No primers M13 For	No primers M13 Rev
E	SC2M1-18 M13 For	SC2M1-18 M13 Rev	SC2M1-60 M13 For	SC2M1-60 M13 Rev	NTC M13 For	NTC M13 Rev	SC2M1-18 M13 For	SC2M1-18 M13 Rev	SC2M1-60 M13 For	SC2M1-60 M13 Rev	NTC M13 For	NTC M13 Rev
F	SC2M1-29 M13 For	SC2M1-29 M13 Rev	SC2M1-62b M13 For	SC2M1-62b M13 Rev	Positive control M13 For	Positive control M13 Rev	SC2M1-29 M13 For	SC2M1-29 M13 Rev	SC2M1-62b M13 For	SC2M1-62b M13 Rev	Positive control M13 For	Positive control M13 Rev
G	SC2M1-38 M13 For	SC2M1-38 M13 Rev	SC2M1-64 M13 For	SC2M1-64 M13 Rev	Blank	Blank	SC2M1-38 M13 For	SC2M1-38 M13 Rev	SC2M1-64 M13 For	SC2M1-64 M13 Rev	Blank	Blank
H	SC2M1-39 M13 For	SC2M1-39 M13 Rev	SC2M1-68 M13 For	SC2M1-68 M13 Rev	Blank	Blank	SC2M1-39 M13 For	SC2M1-39 M13 Rev	SC2M1-68 M13 For	SC2M1-68 M13 Rev	Blank	Blank

4.4. Seal the plate. Vortex for 2–3 seconds, then centrifuge briefly (5–10 seconds) at 1,000 x g.

4.5. Place the plate into a thermal cycler and run the following program:

Parameter	Stage/step						
	Post PCR cleanup	Post PCR inactivation	Polymerase activation	Cycling (25 cycles)			Hold
				Denaturation	Annealing	Extension	
Temperature	37°C	80°C	96°C	96°C	50°C	60°C	4°C
Time	15 min	2 min	1 min	10 sec	5 sec	75 sec	Indefinitely

## 5. Sequencing cleanup

5.1. Spin the reaction plate at 1,000 x *g* for 1 minute, then remove the seal.

5.2. Prepare a mix with SAM Solution and BigDye XTerminator™ Solution in an appropriately sized tube.

5.2.1. Calculate the amount of SAM Solution and XTerminator Solution needed for all samples. You will need 45 µL of SAM Solution and 10 µL of XTerminator Solution per well.

5.2.2. Add the calculated volume of SAM Solution to a new tube using a conventional pipette tip.

**Note:** Make sure there are no particulates in the SAM Solution before pipetting. If there are particulates, heat the SAM Solution to 37°C and mix to dissolve. Cool to room temperature before using.

5.2.3. Vortex the bulk container of XTerminator Solution at maximum speed for at least 10 seconds, until the solution is homogeneous.

5.2.4. Using a wide-bore pipette tip, add the calculated volume of XTerminator Solution to the tube.

**IMPORTANT:** Avoid pipetting from the top of the liquid.

5.2.5. Mix the tube of combined reagents until homogeneous.

5.3. Add 55 µL of the SAM Solution/XTerminator Solution mix to each well.

**IMPORTANT:** Avoid pipetting from the top of the liquid. When aliquoting into the plate, re-vortex the SAM Solution/XTerminator Solution mix every 8–10 wells to homogenize the bead mixture.

5.4. Seal the plate with Applied Biosystems™ MicroAmp™ Optical Adhesive Film. Make sure the plate is sealed well.

5.5. Vortex the reaction plate for 40 minutes.

5.6. In a swinging-bucket centrifuge, spin the plate at 1,000 x *g* for 2 minutes.

## 6. Collect data

6.1. Make sure the instrument is calibrated with the correct sequencing standard (Z-dye set matrix and sequencing standard)

- For details, see the Applied Biosystems™ 3500/3500xL Genetic Analyzer User Guide or SeqStudio™ Genetic Analyzer Getting Started Guide.

6.2. Remove the MicroAmp film and replace it with a 96-well plate septum.

6.3. Load plates into the genetic analyzer.

6.4. Select or create an appropriate run module according to your capillary length, number of capillaries, and polymer type on your instrument. The recommended default run modules are listed below:

- For 3500xL instruments with 50 cm capillaries:

- Instrument protocol: BDxFastSeq50\_POP7xl\_Z

**Note:** Replace 50 with 36 in the instrument protocol name if you have a 36 cm capillary installed.

- Analysis Module: BDTv3.1\_PA\_Protocol-POP7

- For SeqStudio instruments:

- MedSeqBDX



## 7. Analyze results using a sequencing program

Sequence Scanner v2.0 is free software for viewing electropherograms. It provides an easy way to perform a high-level sequencing data quality check or general data review that includes summary tables and electropherograms as well as a general raw or analyzed data view for .ab1 files.

7.1. To obtain the software, go to [thermofisher.com/pages/WE28396/](https://www.thermofisher.com/pages/WE28396/)

7.2. Using Sequence Scanner Software v2.0, generate a QC report. For each sequencing trace, determine the trace score, CRL, and the QV20+ score.

7.3. Suggested acceptance criteria:

- A sequencing trace is acceptable as positive if two of the three thresholds are met:
  - Trace score greater than 31
  - CRL greater than 50
  - QV20+ greater than 50
- A sequencing trace is acceptable as negative if two of the three thresholds are met:
  - Trace score less than 14
  - CRL less than 24
  - QV20+ less than 24
- Sequencing traces that do not fit the above criteria are indeterminate and should be repeated.

7.4. Using BLAST™ alignment or another sequence alignment tool, align positive traces to the SARS-CoV-2 genome.

- Alignments greater than 85% over read length are considered homologous to the SARS-CoV-2 genome
- Discard any sequences that are not homologous to SARS-CoV-2

7.5. For lineage analysis in any of the amplicons, these criteria should be met:

- Positive (passable) traces in both directions (7.3)
- Homology to the SARS-CoV-2 genome (7.4) in regions outside the putative lineage
- Negative traces in no-template control reactions (7.3)

7.6. Test runs that fail for reasons not attributable to system performance, such as equipment malfunction, operator error, or other demonstrable cause, will be designated as invalid runs. Invalid runs will be retested and documented in the study report(s).

## References

1. Wise J (2020) Covid-19: New coronavirus variant is identified in UK. *BMJ* Dec 16;371:m4857. doi: [10.1136/bmj.m4857](https://doi.org/10.1136/bmj.m4857)
2. Tegally H et al. (2020) Emergence and rapid spread of a new severe acute respiratory syndrome-related coronavirus 2 (SARS-CoV-2) lineage with multiple spike mutations in South Africa. *medRxiv* <https://doi.org/10.1101/2020.12.21.20248640>
3. Public Health England (2020) Investigation of novel SARS-COV-2 variant: Variant of Concern 202012/01. Technical briefing 1. [https://assets.publishing.service.gov.uk/government/uploads/system/uploads/attachment\\_data/file/947048/Technical\\_Briefing\\_VOC\\_SH\\_NJL2\\_SH2.pdf](https://assets.publishing.service.gov.uk/government/uploads/system/uploads/attachment_data/file/947048/Technical_Briefing_VOC_SH_NJL2_SH2.pdf)
4. Paden CR et al. (2020) Rapid, sensitive, full-genome sequencing of severe acute respiratory syndrome coronavirus 2. *Emerg Infect Dis.* 26(10):2401-2405. doi.org/10.3201/eid2610.201800

Find out more at [thermofisher.com/sangercoronavirus](https://thermofisher.com/sangercoronavirus)

**ThermoFisher**  
SCIENTIFIC