Maxima H Minus Reverse Transcriptase

Engineered to impress

Thermo Scientific™ Maxima™ H Minus Reverse Transcriptase enables consistent success in cDNA synthesis in RT-PCR and RT-qPCR applications. Developed through molecular evolution, with the introduction of multiple favorable mutations into wild-type Moloney murine leukemia virus (M-MuLV) reverse transcriptase, the enzyme delivers impressive thermostability, processivity, and cDNA synthesis rates.

Why use Maxima H Minus Reverse Transcriptase?

Maxima H Minus Reverse
Transcriptase is an enzyme with
significantly increased thermostability,
enhanced processivity, and
diminished RNase H activity. Due
to these improved features, the
enzyme consistently produces high
yields of full-length cDNA even at
elevated temperatures. The enzyme
is available as a stand-alone enzyme
or formulated for convenient kits to
support different needs for first-strand
cDNA synthesis.

Features

- High yields of cDNA over a wide temperature range (up to 65°C) (Figure 1)
- Efficient RT-PCR even with extremely long amplification targets (up to 20 kb) (Figure 2)
- Enhanced linearity, sensitivity, and early C, in two-step RT-qPCR (Figure 3)
- Multiple formats, from stand-alone enzyme to first-strand cDNA synthesis kit, double-stranded cDNA synthesis kit, and one-tube reverse transcription (RT) master mix

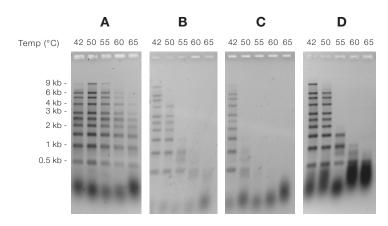


Figure 1. Maxima H Minus Reverse Transcriptase generates full-length cDNA at a wider range of reaction temperatures than reverse transcriptases from other suppliers. cDNA synthesis using 1 µg of Invitrogen™ Millennium™ RNA Markers (poly(A)-tailed) with an oligo(dT)₁₈ primer was performed with (A) Maxima H Minus Reverse Transcriptase; and reverse transcriptases from other suppliers: (B) TaKaRa™ PrimeScript™ Reverse Transcriptase, (C) Promega™ GoScript™ Reverse Transcriptase, at different temperatures per suppliers' instructions. The optimal temperatures are 50°C for Maxima H Minus Reverse Transcriptase and 42°C for reverse transcriptases from other suppliers. cDNA products were detected by alkaline gel electrophoresis.



Technical details

- Supports a wide range of starting total RNA amounts (1 pg–2.5 μg)
- Can synthesize RNA at elevated temperatures (42–65°C)
- RT reaction can be completed in only 15–30 minutes
- Up to 50x higher processivity compared with wild-type M-MuLV reverse transcriptase enzyme
- Demonstrates templateswitching activity
- Capable of incorporating modified nucleotides

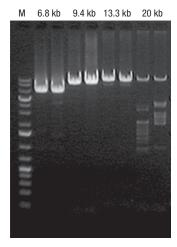


Figure 2. Two-step long RT-PCR from RNA transcripts up to 20 kb. Total RNA was extracted from human and mouse cells and used to reverse-transcribe 6.8 kb and 9.4 kb transcripts from the human RNA, and 13.3 kb and 20 kb transcripts from the mouse RNA, using 1 µg of total RNA per reaction and Maxima H Minus Reverse Transcriptase. The synthesized cDNAs were used as templates for PCR, and the PCR products were visualized by gel electrophoresis. M: Thermo Scientific™ GeneRuler™ 1 kb Plus DNA Ladder.

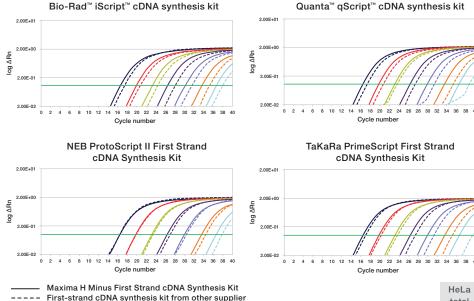


Figure 3. The Thermo Scientific™ Maxima™ H Minus First Strand cDNA Synthesis Kit demonstrates consistently better RT efficiency than that obtained with other suppliers' kits. RT-qPCR of the human β₂ macroglobulin gene was performed on 10-fold serial dilution of 1 μg–1 pg of total RNA from HeLa cells. First-strand cDNA was generated using the Maxima H Minus First Strand cDNA Synthesis Kit and four other commercial first-strand cDNA synthesis kits. cDNA was amplified using Applied Biosystems™ TaqMan® Universal Master Mix II with uracil N-glycosylase (UNG), on the Applied Biosystems™ ViiA™ 7 Real-Time PCR System.

total
RNA
1 µg
100 ng
10 ng
1 ng
100 pg
1 pg
1 pg

Ordering information

Product	Quantity	Cat. No.
Maxima H Minus Reverse Transcriptase	2,000 U	EP0751
	10,000 U	EP0752
	4 x 10,000 U	EP0753
Maxima H Minus First Strand cDNA Synthesis Kit	20 rxns/100 rxns	K1651/K1652
Maxima H Minus First Strand cDNA Synthesis Kit, with dsDNase	20 rxns/100 rxns	K1681/K1682
Maxima H Minus Double-Stranded cDNA Synthesis Kit	10 rxns	K2561
Maxima H Minus cDNA Synthesis Master Mix	50 rxns/200 rxns	M1661/M1662
Maxima H Minus cDNA Synthesis Master Mix, with dsDNase	50 rxns/200 rxns	M1681/M1682

Find out more at thermofisher.com/maxima



In the United States:

For customer service, call 1-800-766-7000 To fax an order, use 1-800-926-1166 To order online: fishersci.com

In Canada

For customer service, call 1-800-234-7437 To fax an order, use 1-800-463-2996 To order online: fishersci.ca



For Research Use Only. Not for use in diagnostic procedures. © 2017 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. Promega and GoScript are trademarks of Promega Corporation. NEB and ProtoScript are trademarks of New England Biolabs. TaKaRa and PrimeScript are trademarks of Takara Holdings Inc. TaqMan is a registered trademark of Roche Molecular Systems, Inc., used under permission and license. Bio-Rad and iScript are trademarks of Bio-Rad Laboratories, Inc. Quanta and qScript are trademarks of Quanta BioSciences, Inc. BN0605133 0817