

TrueDesign Genome Editor

Workflow guide for creating a gene knockout using insertion of stop codons


The Invitrogen™ TrueDesign™ Genome Editor is easy-to-use, free online software for designing and ordering the reagents needed for precise genome editing by homology-directed repair with RNA-guided nucleases and single-stranded DNA donors.

This workflow guide will walk through the steps for creating a gene knockout by insertion of stop codons. To accomplish gene knockout using imperfect repair following a double-strand break, see our workflow guide for creating a gene knockout by causing frameshift indels.


Please note that algorithm-designed CRISPR gRNAs optimized for functional knockout are available for all human and mouse genes and may be ordered directly from [thermofisher.com/truedesign](https://www.thermofisher.com/truedesign).

Step 1:


Go to [thermofisher.com/truedesign](https://www.thermofisher.com/truedesign). Select one of the links to launch the software.

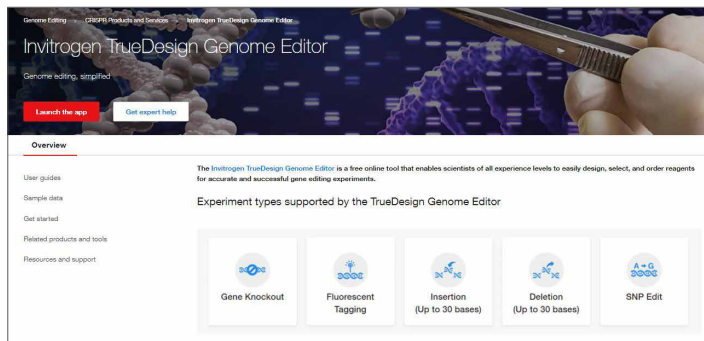

You may be prompted to sign in. Use your existing credentials, or simply provide an email address to register as a new user.

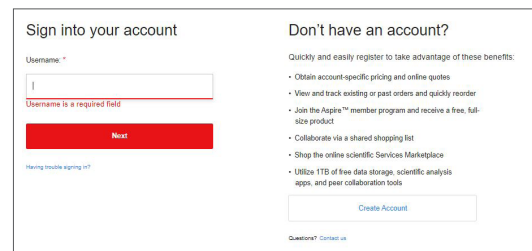


Alternatively, go directly to the **Thermo Fisher™ Connect Platform** and navigate to the TrueDesign Genome Editor.

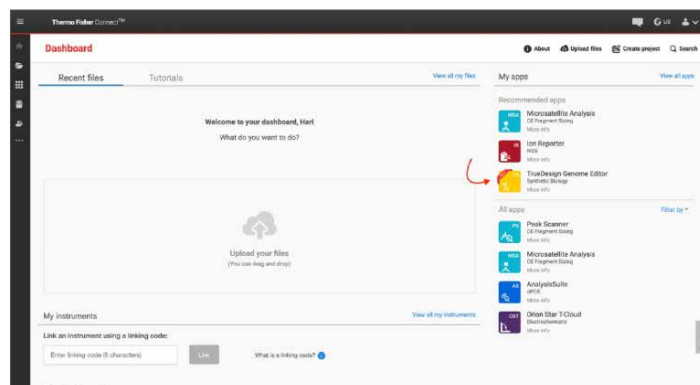

A “Terms of Use” window may pop up. Read the content, scroll to the bottom of the screen, and click “Accept”.


Step 2:

In the TrueDesign software, choose **Gene Knockout** as your experiment type and click “Next”.

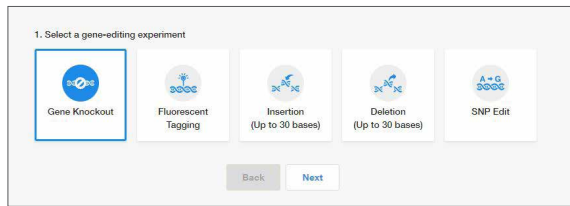

Step 3:


Select the host organism for your knockout experiment and click “Next”.

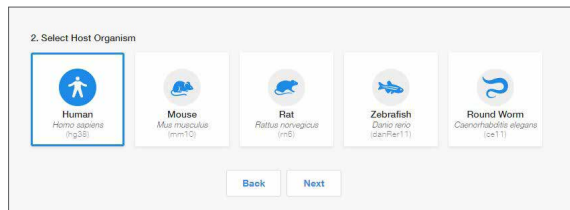

Step 4:


Select **Gene Symbol/Entrez ID** to identify your gene of interest. You may also navigate to a transcript of interest by entering a DNA sequence or chromosome location.

Click “Next”.

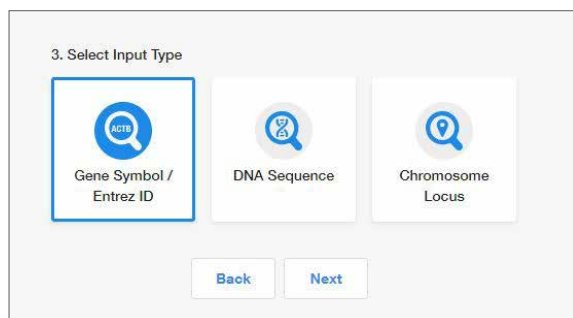

Step 5:


Begin typing the gene symbol or Entrez ID in the gene identifier box. A filtered drop-down list will appear. Select your gene of interest and click “Search Gene”.

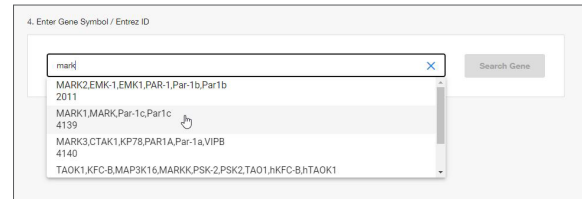

Step 6:

All of the protein-coding transcripts for your selected gene will be displayed. If there is more than one protein-coding transcript and you are unsure of which one to select, click the transcript ID hyperlink to be taken to the NCBI website, where you can better visualize the transcript maps.

After you make a selection, click “Edit”.

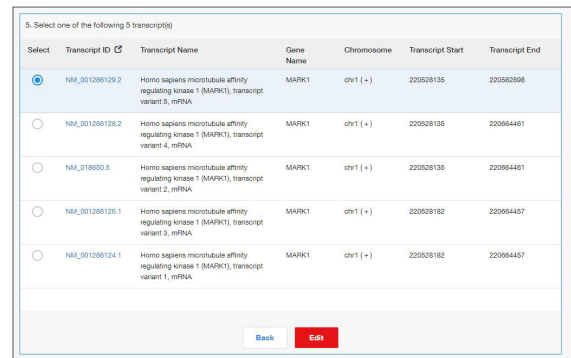

Select	Transcript ID	Transcript Name	Gene Name	Chromosome	Transcript Start	Transcript End
<input checked="" type="radio"/>	NM_001286128.2	Homo sapiens microtubule affinity regulating kinase 1 (MARK1), transcript variant 5, mRNA	MARK1	chr1 (-)	220528135	220582698
<input type="radio"/>	NM_001286128.2	Homo sapiens microtubule affinity regulating kinase 1 (MARK1), transcript variant 4, mRNA	MARK1	chr1 (+)	220528135	22064461
<input type="radio"/>	NM_018500.5	Homo sapiens microtubule affinity regulating kinase 1 (MARK1), transcript variant 2, mRNA	MARK1	chr1 (+)	220528135	22064461
<input type="radio"/>	NM_001286126.1	Homo sapiens microtubule affinity regulating kinase 1 (MARK1), transcript variant 3, mRNA	MARK1	chr1 (+)	220528182	22064457
<input type="radio"/>	NM_001286124.1	Homo sapiens microtubule affinity regulating kinase 1 (MARK1), transcript variant 1, mRNA	MARK1	chr1 (+)	220528182	22064457

Step 7:

The next screen will display the entire transcript’s topology along the top of the screen, with a zoomed-in sequence-level view below.


You will be prompted to choose **Frameshift Indels** or **Insert STOP Codon**.

- **Frameshift indel:** The most common approach for gene knockout, this process relies on imperfect repair following a double-stranded DNA break. You will be required to select a region on the transcript to search for custom CRISPR and TALEN™ designs. Pre-designed gRNAs for human and mouse are available with the **Invitrogen™ TrueGuide™ Synthetic gRNA** product line.

Step 7 (continued)

- **Insert stop codon:** A more precise method for knockout of gene expression, this process uses a DNA donor to insert multiple stop codons to prevent translation. By default, this insertion will be designed adjacent to the transcriptional start site.


For this workflow example, select **Insert STOP Codon**. Please refer to the workflow guide for causing frameshift indels for details on that experiment type.

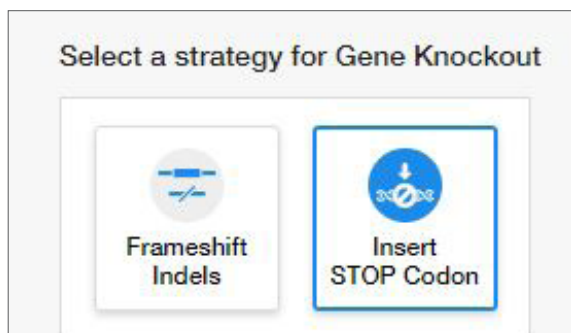

Step 8:


The view will update with a selection box. Click on the stop codon you wish to insert. The relative frequency of each codon's usage within your selected species will be displayed.

Next, choose to add either 3 or 4 stop codons from the drop-down list. Click "Add".

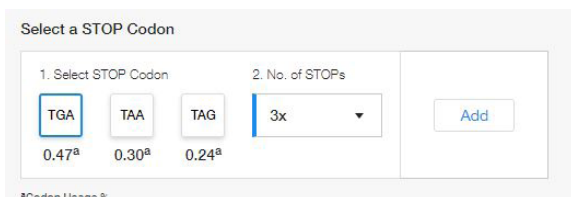

Step 9:

The sequence editor will update to display the location of the inserted stop codons with the inserted DNA highlighted in blue.

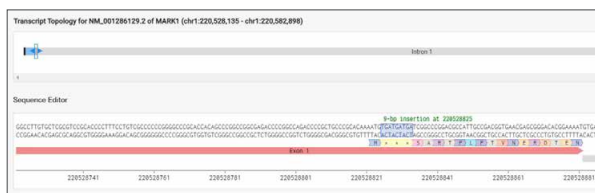

An edit list will appear, displaying the type and location of the edit. You can undo this edit by clicking the blue "undo" arrow in the edit list.


When you are satisfied with your changes, click "Design". This will initiate the design process for the software to find and analyze available TALEN pairs and CRISPR gRNA target regions and check them for specificity.

Step 10:

When complete, you will see a table with two tabs: CRISPR and TALEN targets. View the target location on the sequence editor display by clicking anywhere in the row.

Each row in the CRISPR table displays:

- The CRISPR gRNA design; green checkmarks indicate recommended gRNAs due to score and proximity to the insertion site
- The gRNA PAM site
- The gRNA score, which is a weighted score for efficiency and specificity
- The number of predicted off-targets; click the link to open a pop-up window that displays the location and mismatch information for each potential off-target
- The edit site's distance from the cut site
- The DNA oligo designed as a donor template for the insertion of the stop codons

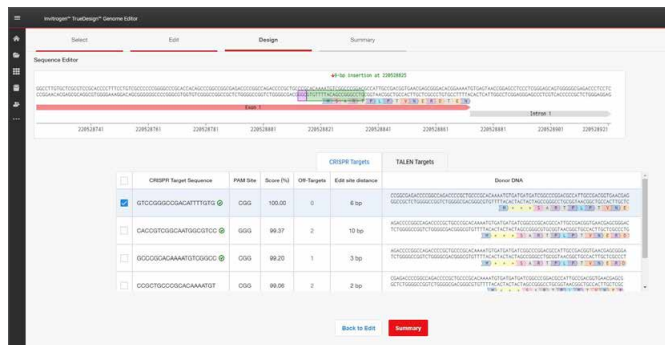

To view the TALEN targets, click on the TALEN tab of the table and similar information will be displayed for each TALEN pair. TALEN pairs are recommended when there are no PAM sites within 10 bp of the knock-in site, or if the efficiency and specificity of the gRNAs are not optimal. Green checkmarks in the design results table will indicate the recommended technology. Learn more about TALEN technology at thermofisher.com/tal.

Step 11:


To select one or more CRISPR gRNAs or TALEN pairs to add to your experiment, use the checkboxes in the table and click “Summary”.

Step 12:

The summary page will display all the reagents needed for your knockout experiment and give you the opportunity to add additional products to complete your workflow.

For a stop codon insertion experiment using CRISPR technology, the gRNAs and donor oligos selected in the design step are automatically included. The tool will also add Invitrogen™ TrueCut™ Cas9 Protein v2 and Invitrogen™ Lipofectamine™ CRISPRMAX™ Cas9 Transfection Reagent.

Products may be deselected from the product summary area, or different quantities may be selected from the drop-down lists.

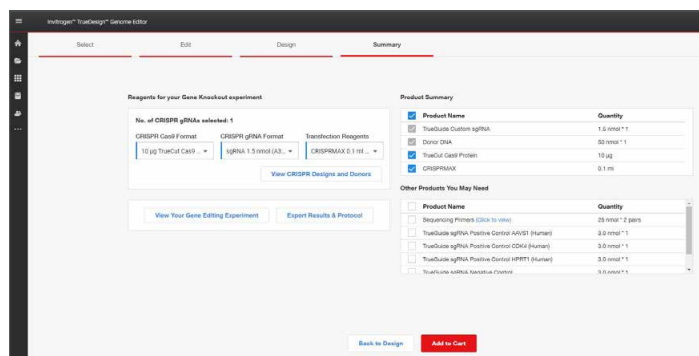

Step 13:

You may use the checkboxes to add additional items such as sequencing primers or positive and negative experimental controls.

When you have completed your product selections, click “Add to Cart” for easy one-step ordering of all selected reagents.

If “Add to Cart” is not enabled in your region or you want to send the list of reagents to your purchasing agent, you can download and save a detailed report of your experiment by clicking “Export Results & Protocol”. The resulting Microsoft™ Excel™ file contains multiple tabs that include all the designs generated by the software, plus all of the gene-specific experimental details and ordering information.



Get started at thermofisher.com/truedesign