

Updating the Bead Lot 175946 expiration date in the Attune™ Acoustic Focusing Cytometer

Each lot of Attune™ Performance Tracking beads has an associated bead lot file that can be downloaded from this website and then imported into the Attune™ software version 1.2.5 or later. The available bead lot files can be downloaded from the Attune™ Performance Tracking Beads (Cat. 4449754) web page located within the Thermo Fisher website:

<https://www.thermofisher.com/order/catalog/product/4449754>

To download and install the bead lot file with updated expiration date, please follow the steps below:

1. Download and save the new bead lot file:
 - a. Visit the Attune™ Performance Tracking Beads (Cat. 4449754) web page located within the Thermo Fisher website.
 - b. Under the Documents section of the webpage, download and save the new bead lot file for lot 1759476 (called Attune (not NxT) Performance Tracking Bead Lot File – Jan 2019 – Lot 1759476). Note that bead lot files are .csv files.
 - c. On your Attune™ Acoustic Focusing Cytometer computer, save the updated .csv file (4449754 Lot 1759476_UPDATED.csv)
2. Remove the old bead lot file:
 - a. On your Attune™ Acoustic Focusing Cytometer computer, navigate to X:\Users\Public\Documents\Applied Biosystems\Attune 2.1\Performance Tracking\Beadlots\ 4449754 Lot 1759476.csv.
 - b. Delete the .csv file (4449754 Lot 1759476.csv)
3. Update the bead lot file in the software using the new file that was saved to the computer.
 - a. In the Attune™ cytometric software under performance test, click Get New Lot # to navigate to and select the bead lot file. It will automatically populate the drop-down list and update the expiration date.



4. Go to the Performance Test Module.
5. Re-run baseline and performance test using the new bead lot file.

For other questions or concerns, please contact your local technical support team.