

Release Notes

Qubit™ 4 Fluorometer Firmware v2.10

Contents

OVERVIEW

FEATURES

DOWNLOAD INSTRUCTIONS

UPGRADE INSTRUCTIONS

LEGAL STATEMENTS

Overview

Firmware v2.10 for the Qubit™ 4 Fluorometer

This firmware release is for Qubit4™ Fluorometer only:

Q33226 Qubit™ 4 Fluorometer

New Features/Changes of version 2.10

1. Add support for Wi-Fi network using [Thermo Fisher High-Power USB Wi-Fi module](#)
2. Enable Thermo Fisher Cloud connectivity
3. Add Verification Assay support

Downloading the Firmware File

Download the attached file onto a USB memory stick (shown in .pak format).

1. Save/Transfer the .pak file (not in any folder) on your USB stick.
2. Make sure that no other .pak files are present on the memory stick.

Updating your Qubit™ instrument

1. Power on the instrument.
2. Insert the USB memory stick containing the firmware update into one of the USB ports in the Qubit™ instrument.
3. On the Home screen, touch 'Settings'.
4. In the Settings menu, touch 'Update Software'.
5. If you have downloaded the firmware properly, the instrument should locate the file.
6. Press 'Upgrade' at the bottom of the screen to confirm that you want to upgrade.
IMPORTANT! Ensure that you do not turn off the instrument when the upgrade is in progress.
7. When the upgrade process is complete, the instrument will restart automatically. When it has rebooted, you can remove your USB stick.

LEGAL STATEMENTS

Important licensing information:

These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

The trademarks mentioned herein are the property of Life Technologies Corporation and/or its affiliate(s) or their respective owners.

LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

For Research Use Only. Not for use in diagnostic procedures.

© Copyright 2018 Life Technologies Corporation. All rights reserved.