


# CaptureSelect LambdaXP Affinity Matrix – New Product Introduction


New product introduction – August 2021

 The world leader in serving science

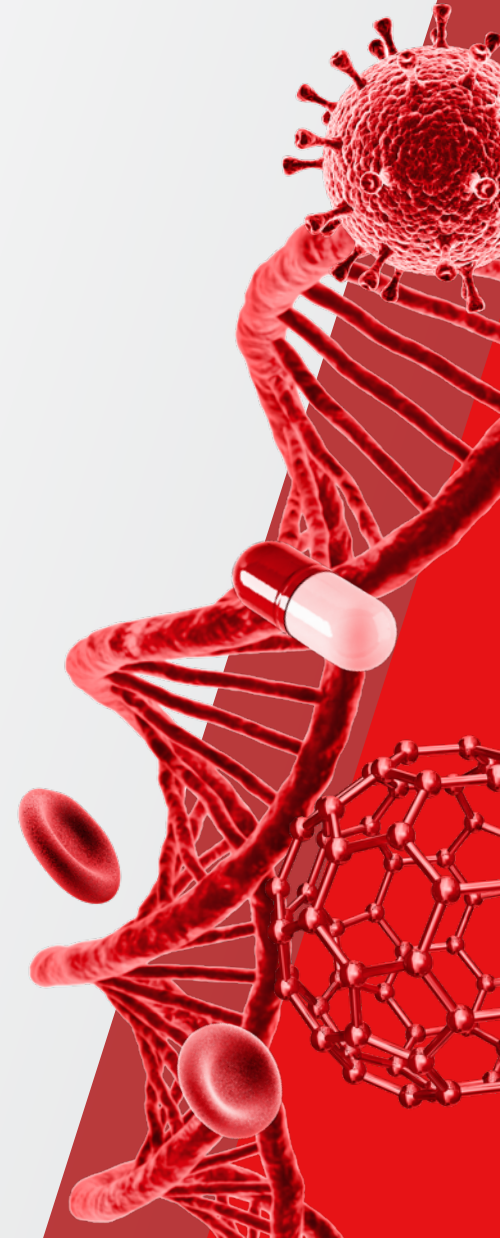


# Thermo Scientific CaptureSelect LambdaXP Resin

## Next-generation lambda light chain binder

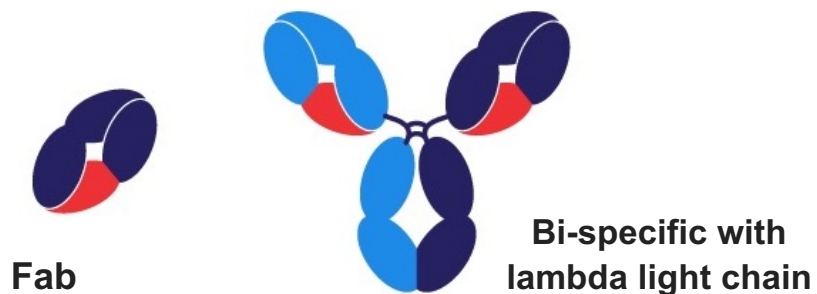


- ✓ Generic: 100% lambda subtype coverage for all immunoglobins containing a lambda light chain
- ✓ Specific to human, does not bind to bovine antibodies
- ✓ High binding capacity: >35 g/L IgG
- ✓ Small elution pool volume at pH 3.5 – 4
- ✓ Excellent scalability Animal derived free product

## CaptureSelect™ resin and antibody selectivity

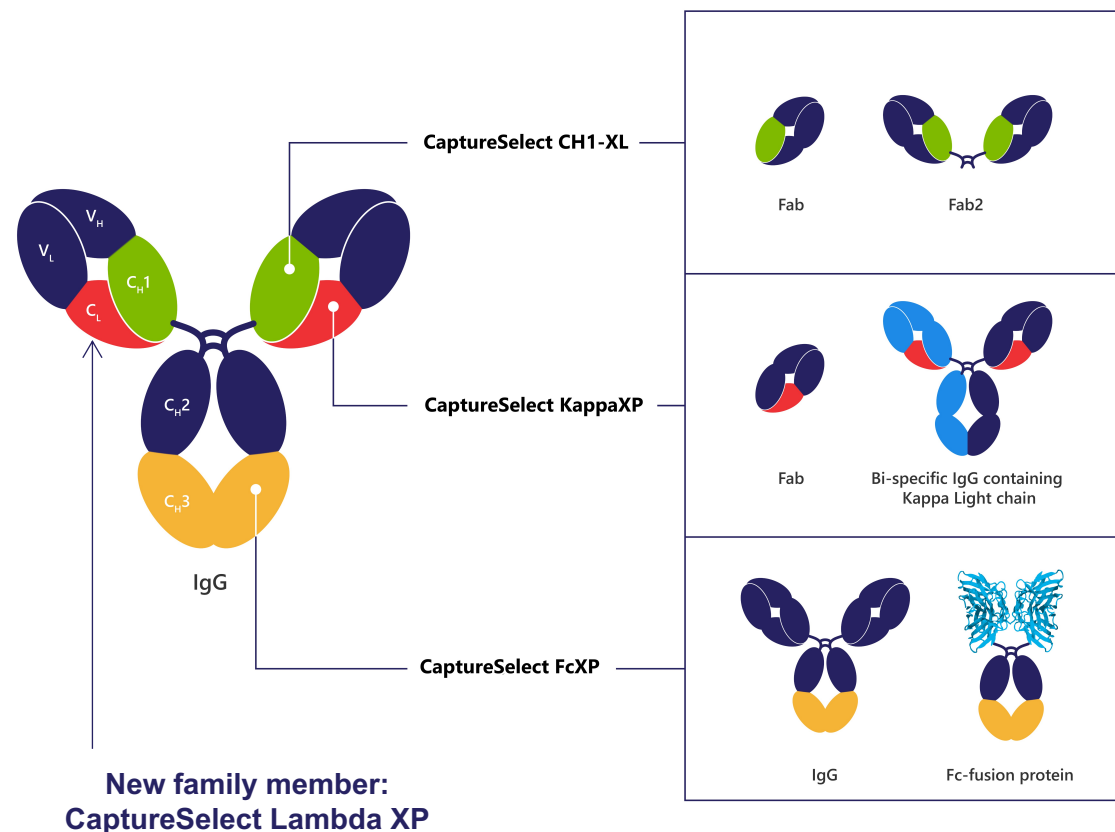


# Bi-specific antibody purification using LambdaXP resin

## Dynamic Binding Capacity at 4 min residence time

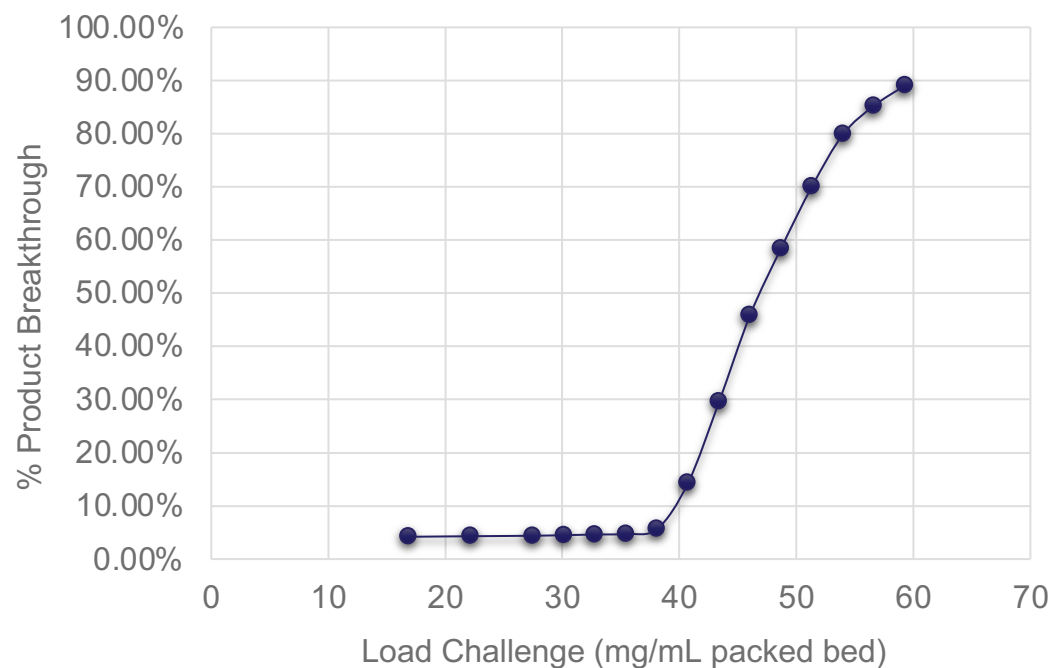

✓ High Dynamic Binding Capacity at 10% breakthrough ~40 mg/mL


## Elution performance – 25 mM sodium acetate

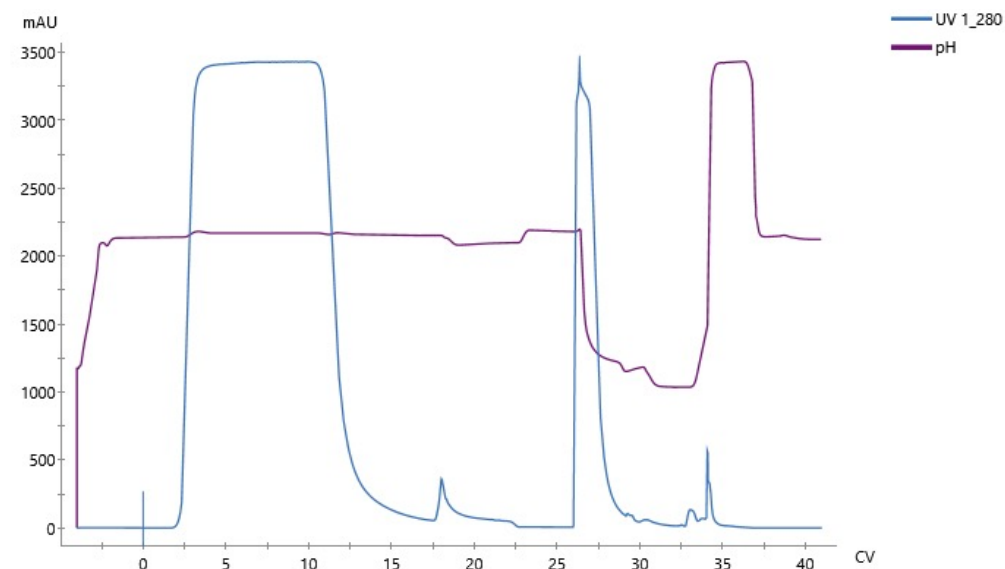

✓ Efficient elution (3 CV) of Bi-specific antibody at pH 3.6 using a load concentration of 32 mg/mL

# CaptureSelect LambdaXP - superior performance compared to alternative affinity resin


## Breakthrough analysis - comparison

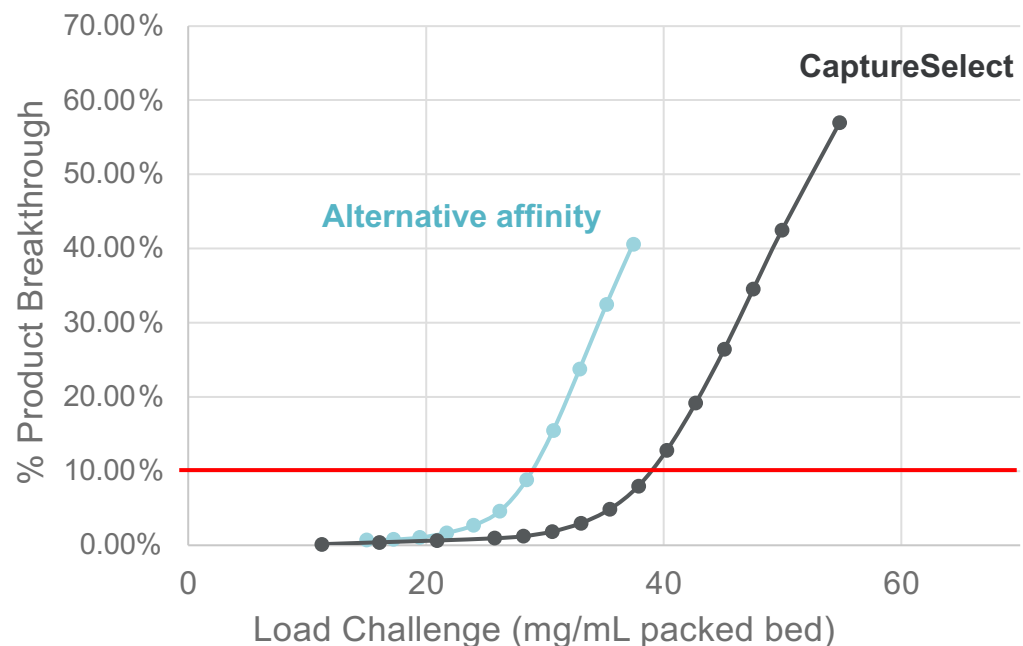

The CaptureSelect™ affinity matrix shows improved binding capacity compared to the alternative affinity resin

Load: Bi-specific antibody – 4 min residence time

## Elution performance – comparison

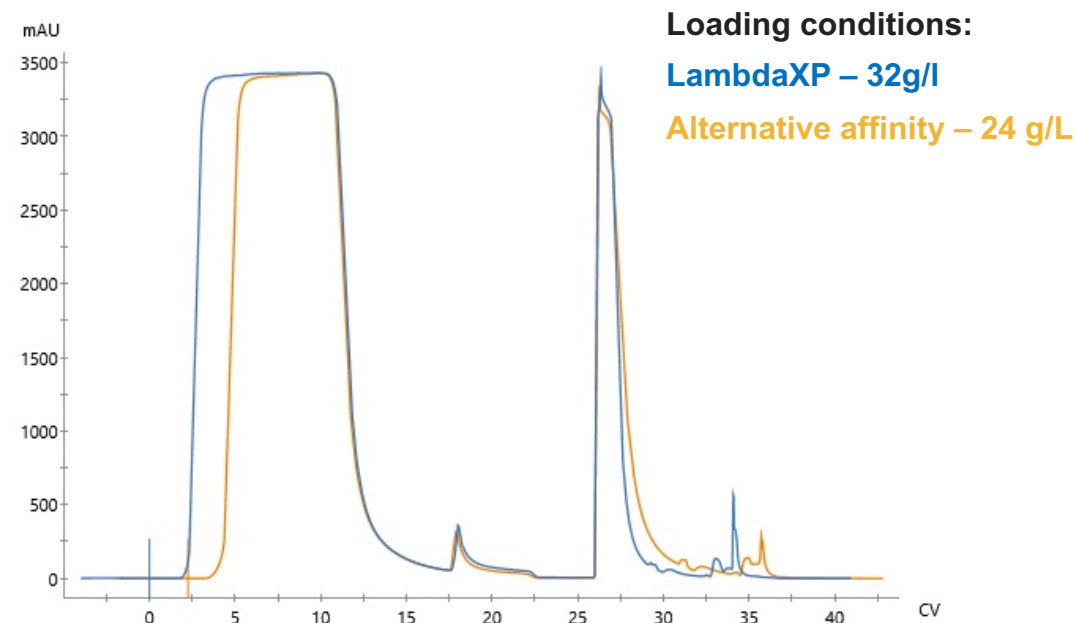

The CaptureSelect™ affinity matrix shows improved elution efficiency compared to the alternative affinity resin

Load: Bi-specific antibody – elution: 25mM sodium acetate, pH 3.6

# Process performance compared to alternative affinity resin



Media	Load Challenge (g/L)	Yield (%)	Aggregate by HP-SEC (%)	Fragment by HP-SEC (%)	HCP (ng/mg)	DNA (ng/mg)	0.5-0.5 OD Elution CV
Alternative affinity resin	24	89	1.2	1.0	1951	< 5.6 E-5	4.5
CaptureSelect™ Lambda XP	32	89	1.3	1.4	1662	< 1.2 E-4	3.0

- Elution with 25mM sodium acetate, pH 3.6

The CaptureSelect LambdaXP resin shows superior performance and impurity clearance compared to the alternative affinity resins

# CaptureSelect LambdaXP advised cleaning strategy



## Acidic strip after every run

- 0.1-0.5 M Citric Acid pH 2 (has chelating properties)
- When preferred, other acids can be used



## Additional cleaning/sanitization steps (when needed)

- Process optimization mainly depending on type of feed
  - Cleaning after every run, or after 2-5 runs with 20 mM NaOH
- Sanitization with PAB (phosphoric acid, acetic acid, benzyl alcohol) as alternative for NaOH





# CaptureSelect LambdaXP resin characteristics and available products



*Purification of Ig's, Fab, and Fab2 fragments directly from complex source materials in a single step with high purity and yield.*

## MAIN RESIN CHARACTERISTICS

- Matrix:** agarose-based, epoxide activated
- Average particle size:** 65 ± 10 µm
- Ligand:** CaptureSelect™ LambdaXP affinity ligand
- Ligand immobilization method:** Epoxide immobilization of the ligand
- Binding capacity:** >35 g IgG per liter resin depending on flow rate, column height, and residence time
- Elution conditions:** 25 mM sodium acetate, pH 3.5
- Flow characteristics:** 150–300 cm/h (up to 2 bar)
- Formulation buffer:** 20%(v/v) ethanol



SKU	Product
2943752005	CS LambdaXP AFF MTR 5 ML
2943752010	CS LambdaXP AFF MTR 10 ML
2943752050	CS LambdaXP AFF MTR 50 ML
1943752250	CS LambdaXP AFF MTR 250 ML
194375201L	CS LambdaXP AFF MTR 1 L*
194375205L	CS LambdaXP AFF MTR 5 L*
810375201	CS LambdaXP Leakage Elisa -1 ASSAY**
810375210	CS LambdaXP Leakage Elisa -10 ASSAY**
5943752001	CS LambdaXP MiniChrom 1 ML***
5943752005	CS LambdaXP MiniChrom 5 ML***

\* Products come with full regulatory support (RSF) enabling use in commercial manufacturing  
\*\* Available from September  
\*\*\* Available from October

# Thank you

For Research Use or Further Manufacturing. Not for use in diagnostic procedures

© 2021 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

