

Thermo Fisher S C I E N T I F I C

Handle Food Samples with Care for Reliable Rheological Results

Dr. Klaus Oldörp

Overview

- Food and rheology
- Sample handling before the measurement
- The right measuring geometry
- The suitable measuring method



FOOD AND RHEOLOGY

How a Rheologist sees Food

- Natural Food
 - Has been grown
 - Can be tested for texture
 - Natural variations are commonly accepted
- Processed Food
 - Has undergone a production process
 - Can be tested for
 - Properties during production
 - Properties of the final product
 - Constant properties are linked to the brand
 - In spite of variations of raw materials
 - In spite of different production sites



Rheology and Food – Where is the Connection?

During Production

Pumping

Mixing

Cooking

Filling

Storage Stability

During Consumption?

Look

Mouthfeeling

Creaminess

Spreadability

• ...

Viscosity

Viscoelasticity

Yield Stress

Moduli

Temperature Dependence

Time Dependence

Elongational Viscosity



Texture – One of Food's Key Properties

- For the Consumer
 - Part of the pleasure when consuming food
 - Important criterion to accept or reject food
- For the Rheologist
 - Sth. to describe with rheological parameters
 - Sth. to bear in mind when
 - Handling the sample
 - Selecting the measuring geometry
 - Designing the test method
 - A lot of trouble if ignored!



SAMPLE HANDLING

Handling of Texturized Samples

- Visual assessment
 - Does it flow?
 - Does it appear elastic?
- Handle with care
 - No shaking
 - No stirring
- Use suitable tools
 - Spoon, spatula
 - No pipettes etc.
- Possible risks
 - Damage of texture
 - Inclusion of air bubbles











Preparing the Sample into the Rheometer

- Take sample from container carefully
 - Not squeezing it through a nozzle etc.
 - Best using spatula or spoon
- Carefully placing it onto the lower plate
 - Centred
 - Not squeezing or distributing it
- Closing the gap with reduced speed
- Allow suitable recovery time
 - Thermal equilibration
 - Remove mechanical stress

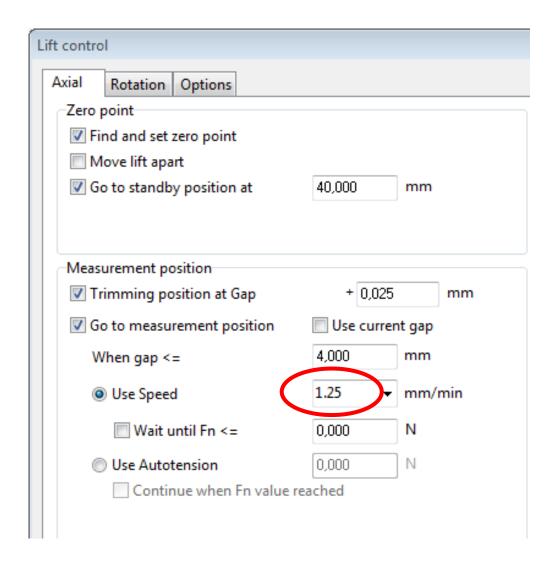

Closing the Gap with Reduced Speed


- Takes more time
- Causes less damage of sample structure
- Saves recovery time afterwards
- Reduces risk of including air bubbles
- Results are less affected by method

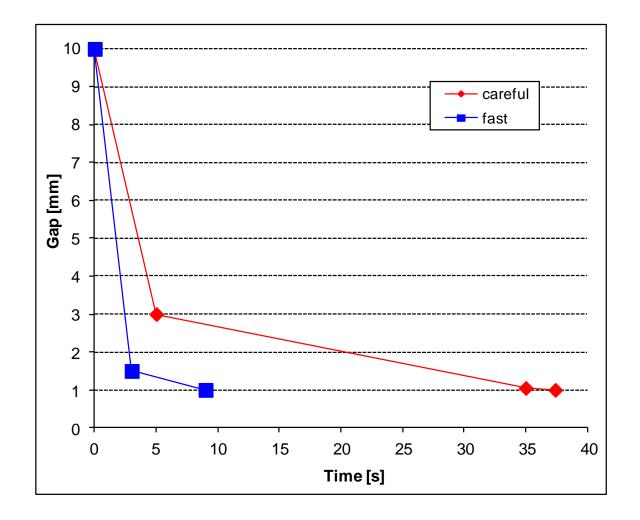

THE MEASURING GEOMETRY


Is There a Universal Measuring Geometry?

- Concentric cylinder geometry
 - Due to historical reasons
 - Large geometry surface increases sensitivity
 - Relatively simple instruments can be used
 - Used in several standards (e.g. ICA 76)
- Disadvantages
 - Difficult for higher viscosities or samples with higher yield stress
 - Significant preshearing when inserting bob
 - Fixed gap size
 - Difficult cleaning
 - Long thermal equilibration time (15 30 min)
 - Heating/cooling ramps very slow

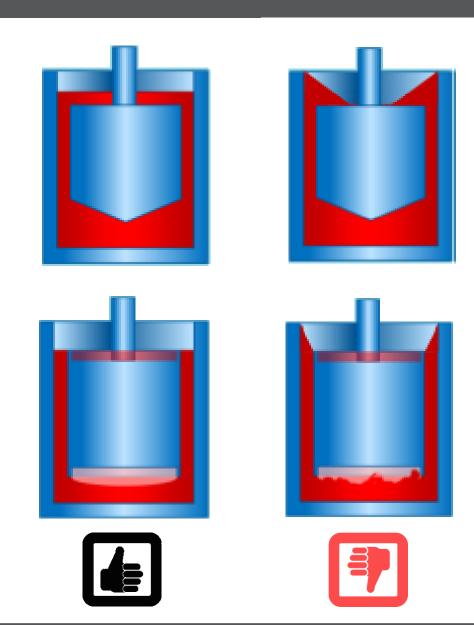

Choose the Right Geometry

- There is no such thing as the universal measuring geometry!
- Have a look at the sample
- Consider the test planned
- Select the size
 - Low viscosities → big geometries
 - High viscosities → small geometries
- When using special geometries
 - It's a relative test
 - Standardize the method

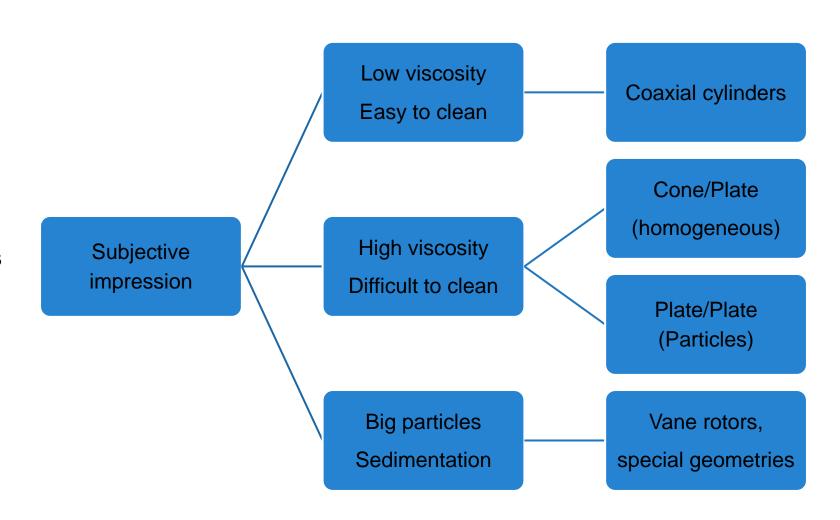


Cone/Plate and Plate/Plate Geometries

- Advantages
 - Smaller sample volume
 - Faster temperature equilibration
 - Much easier to clean
 - Using plate/plate, gap can be adapted
- Disadvantages
 - Drying effects can be more disturbing
 - Sample can leave the gap
 - At higher speeds
 - Due to elastic effects




Samples Containing Particles

- Cone/plate geometry
 - 35 mm / 1°
 - 52 µm gap (fixed)
 - Particles interlock under shear
 - Grinding and blocking
- Plate/plate geometry
 - 35 mm
 - 1 mm gap (flexible)
 - Gap can be adapted to avoid grinding and blocking
- Example
 - Smoothies with seeds

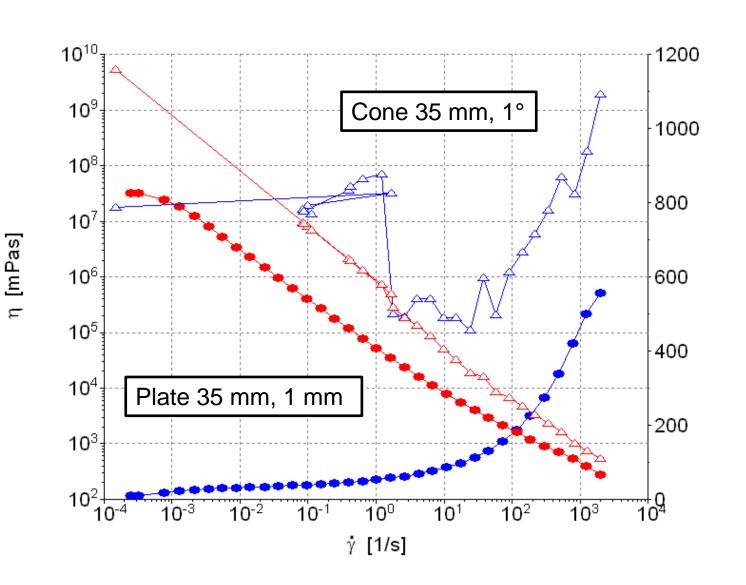


Particles in the Measuring Gap

- Keep in mind for
 - Cone/plate geometries
 - Plate/plate geometries
 - Concentric cylinder geometries
- Rule of thumb
 - Gap > 5 × max. particle diameter
 - Gap > 10 × average particle diameter
- Solution
 - Use bigger gap or cone angle
 - Use vane rotors for relative tests

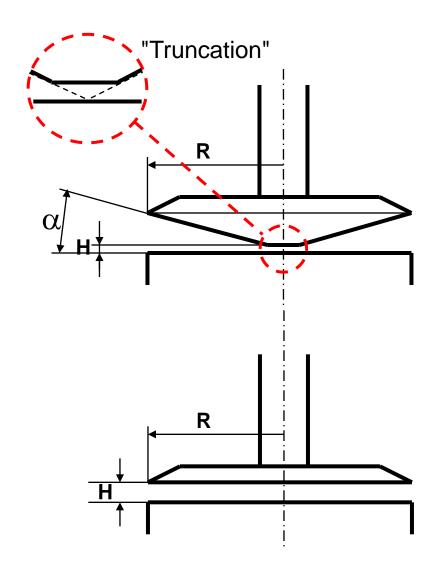


Use Right Amount of Sample

- PP- or CP-geometry: overfill and trim
- CC geometry: use pipette or balance
- Correct amount is given in RheoWin
- If correct amount not possible
 - Overfill with a constant amount
 - Standardize filling procedure for comparable results

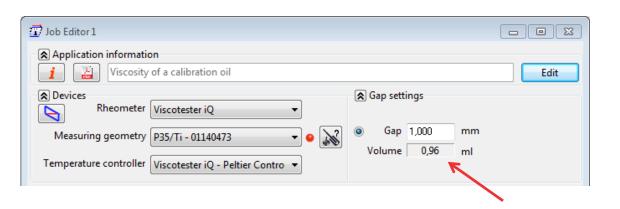


η [mPas]

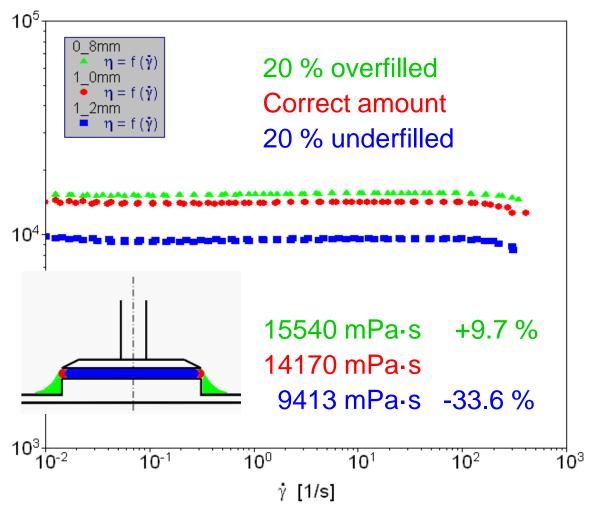

Effect of Underfilling Quantified


- Underfilling reduces contact surface
- Smaller surface results in smaller response
- Rheometer detects e.g. lower viscosity
- Small geometries are affected the most
- Carefully remove excess sample
- Use trimming position



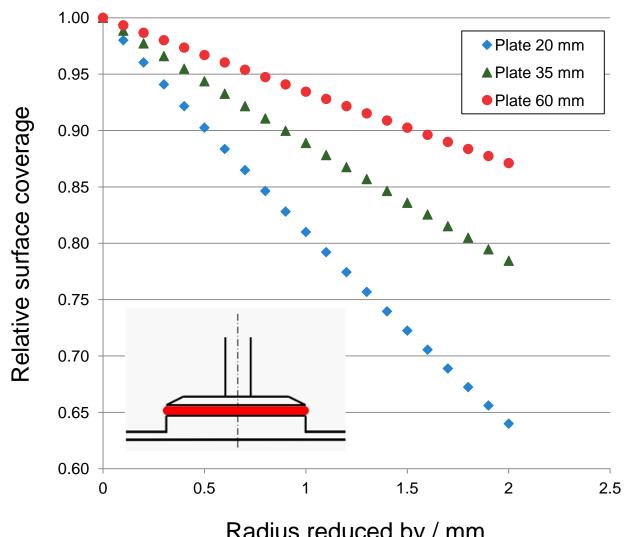

Radius reduced by / mm

- Happens earlier with increasing
 - Radius
 - Gap (or cone angle)
 - Elasticity of sample
- To reduce this effect use
 - Smaller radii
 - Smaller gaps or cone angles
 - CC-geometries
 - Oscillatory measurement

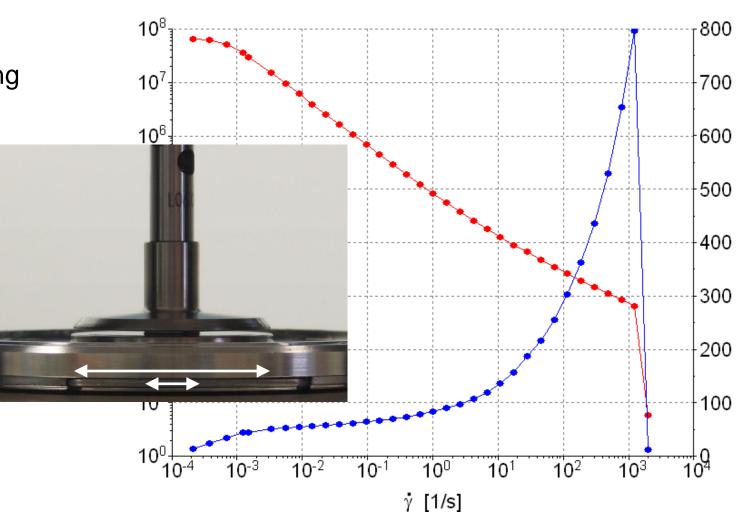



Drying Effects

- The majority of foods contains water
- Evaporation can occur during:
 - Long time measurements
 - Tests at higher temperatures
- Evaporation causes
 - Underfilled geometry
 - Film formation
 - Lost contact to part of the geometry
- Countermeasures:
 - Use solvent trap
 - Cover sample with low viscosity oil
 - Use CC-geometry (plus cover with oil)





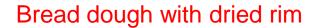


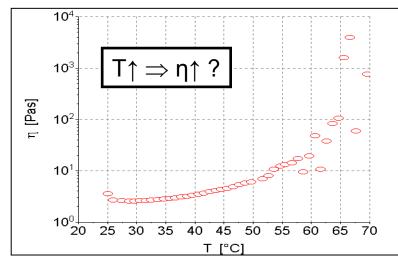



Indication of Drying Sample

- Rheological data looks unsuspicious
- Stress curve reveals reduced contact with upper geometry
- Sometimes F_N can also indicate drying/shrinking

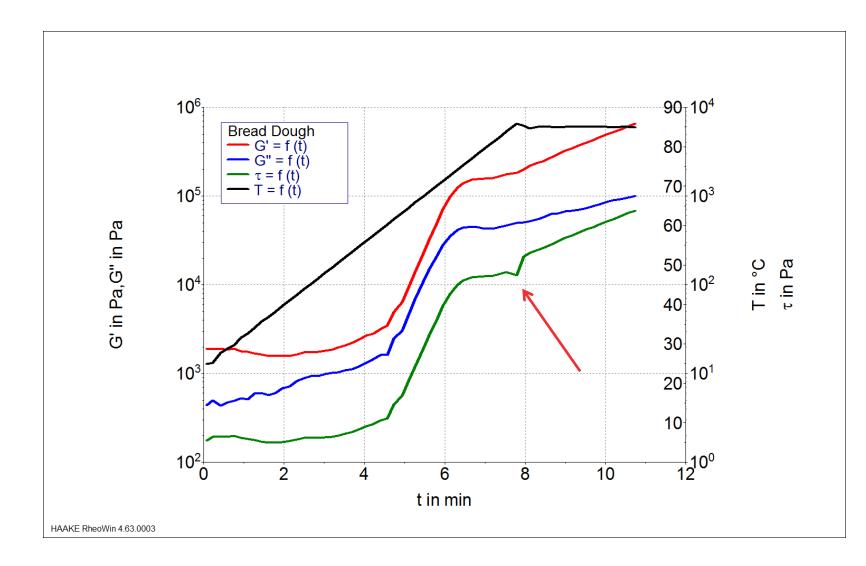

THE MEASURING METHOD


Viscosity

- Most frequently measured rheological parameter
- Can be measured with a variety of methods
 - Capillaries
 - Cups
 - Falling ball viscometer
 - Rotational viscometer
- Many historical methods are used
 - For simplicity
 - To be able to compare results
 - Due to loss of knowledge




Viscosity: The Single Point Measurement

- Popular test because it is fast → QC
 - Often done with constant shear rate
 - Shear rate should relate to application
 - Wait until values stays constant over time
 - Values are taken from plateau
- It's an absolute viscosity (if done like described above)
 - Thermal equilibrium
 - Constant laminar flow
- But it does not show the shear rate dependence
- Always measure viscosity over a suitable shear rate range

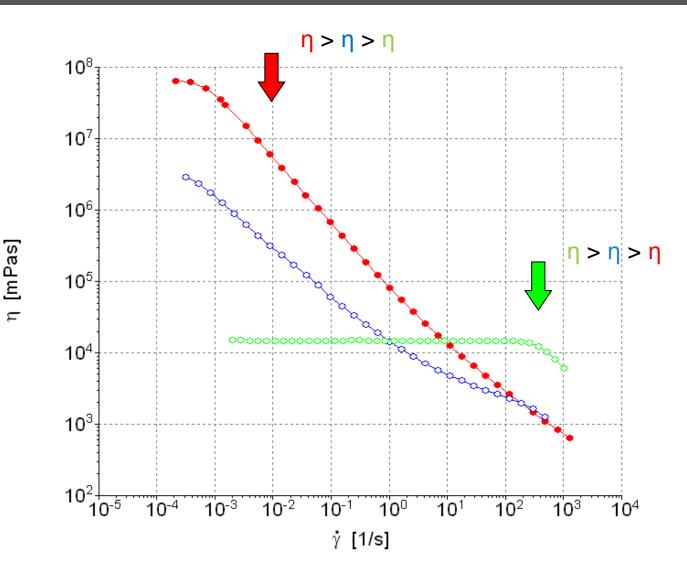


The Rotational Ramp

- Commonly used
- Fast, usually 1 2 min
- Reproducible
- QC method "fingerprint"
- Transient measurement
- Gives no absolute viscosity
- Relative test
- Only comparable using the same ramp

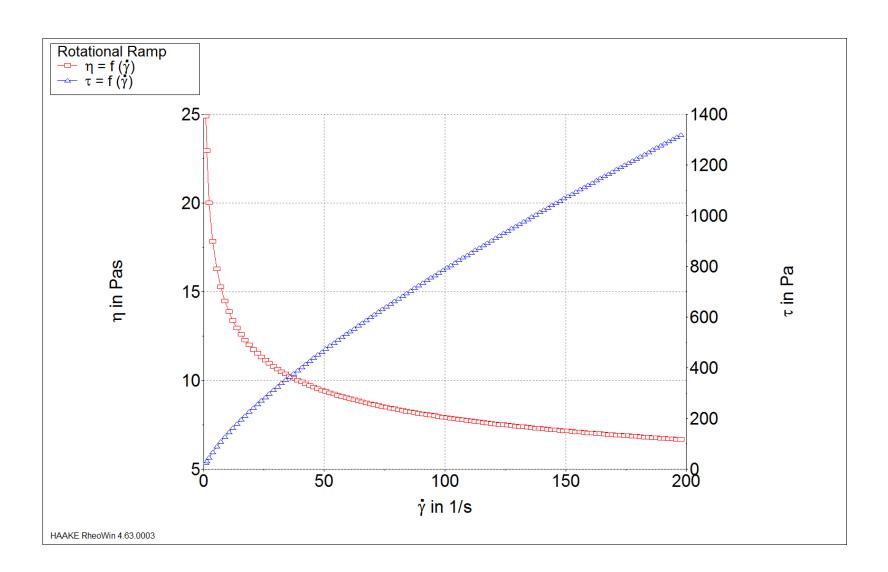


The Real Viscosity Curve

This viscosity curve looks good

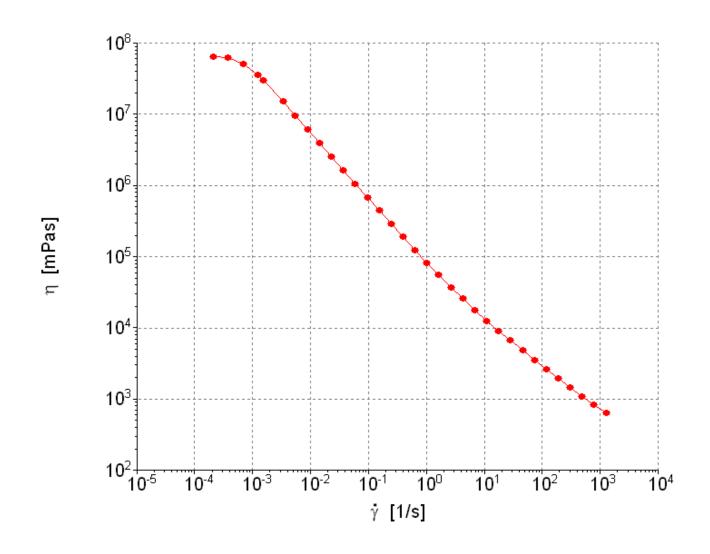

The Real Viscosity Curve

- This viscosity curve looks good
- But this one comes from the same product

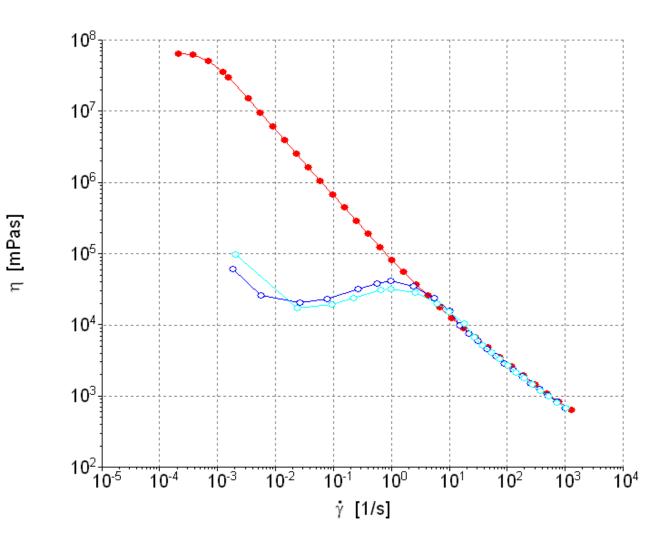


The Real Viscosity Curve

- This viscosity curve looks good
- But this one comes from the same product


[mPas]


- And this one as well.
- The difference lies in the method used
 - Step wise increase of shear rate
 - Linear shear rate ramp
 - Logarithmic shear rate ramp
- Big differences at low shear rates
- Identical results at high shear rates

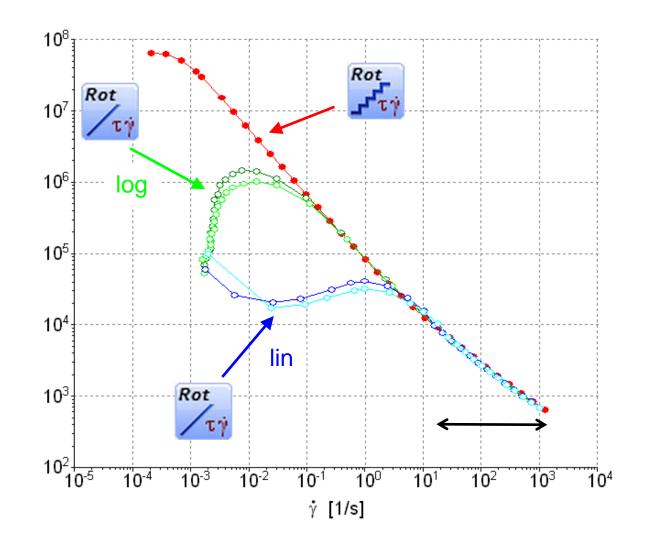

Shear Rate Dependent Equilibrium


- Instrument instantaneously at right speed
- Time to establish flow profile depends on
 - Sample viscosity
 - Sample elasticity
 - Shear rate

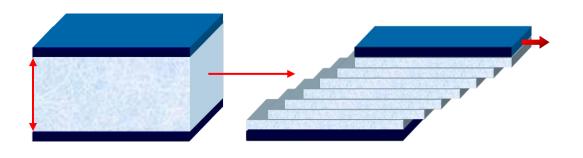


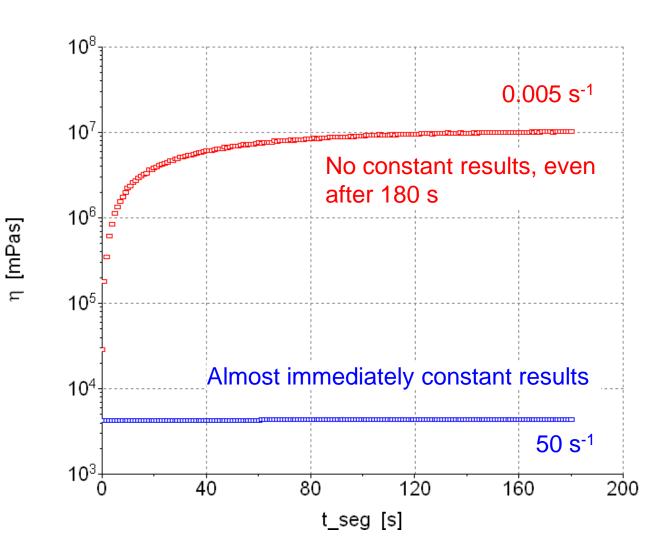

The Time Efficient Viscosity Curve

- Mayor criticism against shear rate step: it takes more time than a shear rate ramp
- This can be true, depending on
 - Shear rate range
 - Number of data points (→ interpolation)
 - Sample's properties

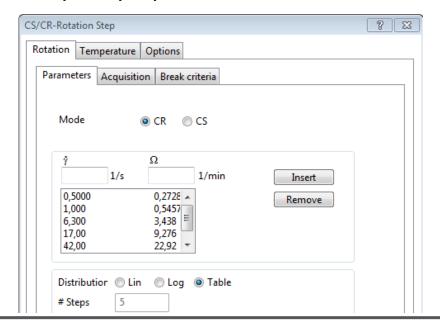

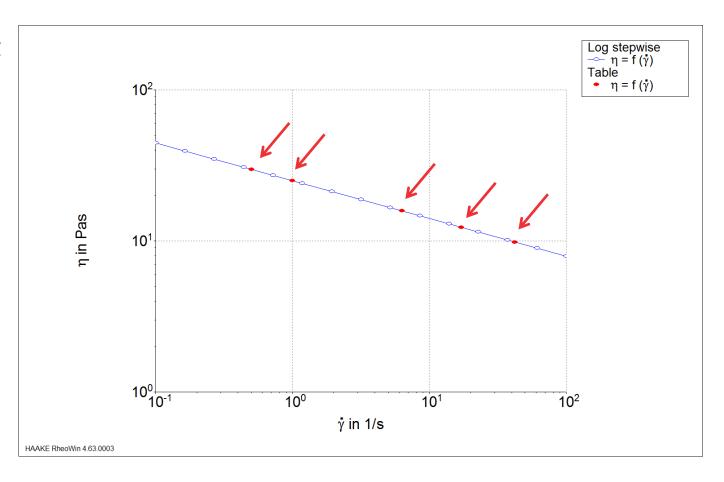

Using a shear rate table saves time but still gives more info!



Rheology Beyond Viscosity

- Viscosity is needed to quantify
 - Pumping (production)
 - Dosing (e.g. syrup from a dispenser)
 - Visual impression (sauces, syrups, molten cheese...)
- It is not sufficient to quantify
 - Long time stability (salad dressing)
 - Gel stability (pudding)
 - Mouth feeling (sausage-meat)
 - •








Oscillatory Measurements – A possible Alternative

- When sample comes out of the gap due to elastic effects
- When geometry loses contact with sample
 - CC geometry in peanut butter
 - CC geometry in mayonnaise
- When you are interested in the sample's structure
 - Emulsion
 - Gel




This is what rotation sees

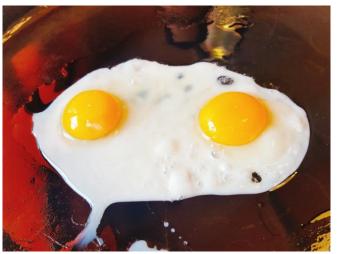

Oscillation sees the undamaged structure

Oscillatory Measurements – Stability of a Structure

- Stability against sedimentation
- Stability of a foam
- Stability of an emulsion
- Related to yield stress

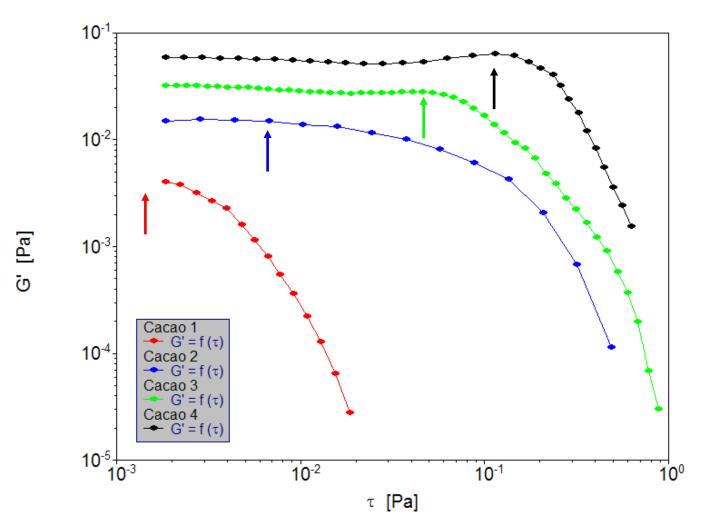


Different chocolate milks with different stabilizers

Oscillatory Measurements – Development of a Structure

- Gelling of protein
- Baking of dough
- Crystallization of fat
- Oscillation
 - does not affect the development of the structure

 Yields valuable information about the structure

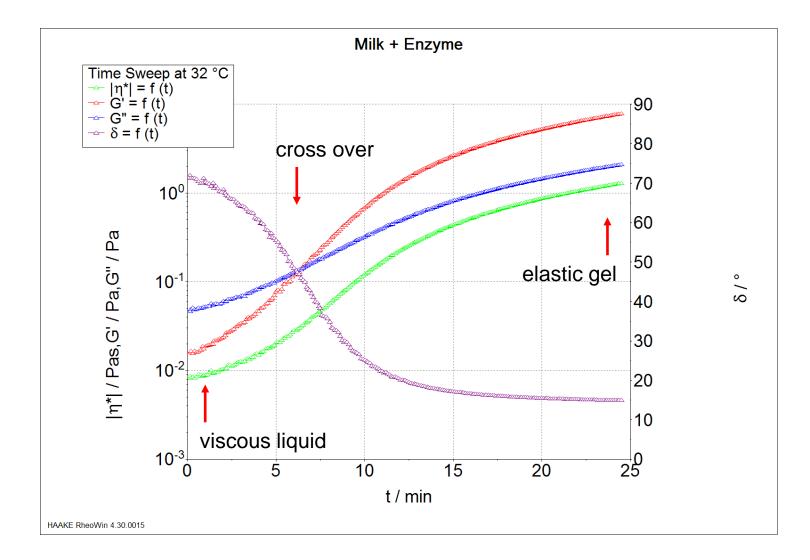

Summary

- Food samples can be tricky to handle
- But reliable results are easily achived when
 - The sample is handled carefully
 - The right geometry is used
 - The measuring methods fits to the sample
- There is no universal test method!
- The measuring method should be
 - More than one dot
 - But still time efficient
 - Sometimes an oscillatory measurement



Thank you for your attention!

Please let me have your questions now!