

Performance Features of an Extended Range Beamsplitter for Mid- and Near-IR Spectroscopy

Michael Bradley, Ph.D., Thermo Fisher Scientific, Madison, WI, USA

Key Words

- Smart NIR Integrating Sphere
- Smart UpDRIFT
- Fiber Optics
- FT-Raman
- Mid-Infrared
- Near-Infrared

The multitude of samples facing the modern analytical laboratory makes versatility in the instrumentation a critical concern. The typical laboratory FT-IR spectrometer is equipped with mid-infrared capabilities (MIR), operating from 4000 to 400 cm^{-1} . However, many applications utilizing the near-IR (NIR) spectral region, from about 10,000 to 4000 cm^{-1} are also critical, including raw material identification, moisture analysis and FT-Raman spectroscopy. The MIR and NIR ranges are typically covered by two instruments, or one instrument with a change in beam splitter. This technical note describes the use of a Thermo Scientific XT-KBr™ beamsplitter to collect both MIR and NIR data. Application data showing the trade-offs in performance is also presented.

The XT-KBr beamsplitter covers the spectral range from 11,000 to 375 cm^{-1} . The detectors and sources need to be selected correctly for optimal performance. The wide range eliminates the need to change the beamsplitter when studying a sample with important MIR and NIR features, or when swapping between MIR and NIR applications. Both of the Nicolet™ series spectrometers can be equipped with the XT-KBr beamsplitter, though the Thermo Scientific Nicolet 6700/8700 spectrometers are better equipped to deal with the multiple ranges through multiple-installed sources and detector, selectable by software.

The data shown were collected using a Nicolet 6700 spectrometer equipped with both the Thermo Scientific Ever-Glo™ ETC™ MIR source and the quartz-halogen NIR source. Raman data were collected using the Thermo Scientific Nicolet NXR FT-Raman Module with an InGaAs room temperature detector. Other conditions were as noted in the experiment description.

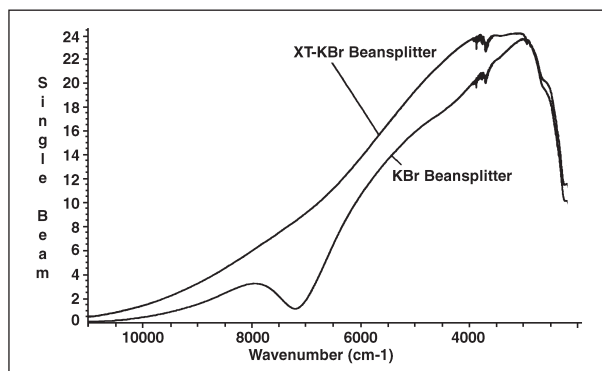


Figure 1

Figure 1 shows single beam (non-ratio) spectra of the XT-KBr and standard KBr beamsplitters using a common scale, collected using identical parameters at 4 cm^{-1} resolution, a DTGS detector and white light source. The XT-KBr beamsplitter shows significantly greater energy beyond 2000 cm^{-1} versus the KBr beamsplitter. The reflective germanium coating on the KBr beamsplitter absorbs strongly at about 7400 cm^{-1} , limiting its use in the NIR, whereas the XT-KBr beamsplitter provides a transparent energy window to 11,000 cm^{-1} .

NIR FT-Raman has become an important tool in providing additional structural information for a wide variety of laboratory samples. The technique has value to analytical labs because it permits analysis with several advantages:

- No sample preparation
- No atmospheric CO_2 and H_2O problems
- Structural elucidation complimentary to FT-IR
- Samples may be housed in glass or plastic
- Spectral data can be obtained down to 100 cm^{-1} (Raman shift)
- Low potential for fluorescence (relative to dispersive Raman)

Figure 2 shows FT-IR and FT-Raman spectra of a powdered laundry detergent. The FT-IR data was collected using a Thermo Scientific Smart™ Performer accessory with the sample pressed against the ZnSe crystal, and a DTGS detector. The sample for the FT-Raman spectrum was placed into an NMR tube. Both spectra were measured using the XT-KBr beamsplitter. No manual change of optical components was required for the measurements. Computerized selection of the FT-Raman accessory can be made from the OMNIC™ spectroscopy software, with changeover from FT-IR to FT-Raman occurring within seconds.

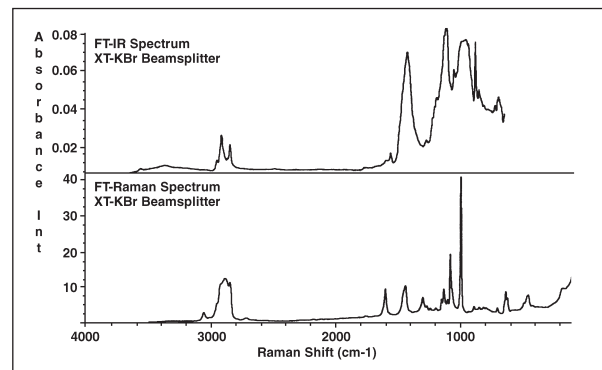


Figure 2

Normally, an NIR beamsplitter such as CaF₂ would have to be installed for the FT-Raman experiment. Comparison of the performance of the CaF₂ and XT-KBr beamsplitter options is shown in Figure 3. These single beam spectra were collected using a DTGS detector and white light source. The CaF₂ beamsplitter clearly provides higher NIR energy beyond 5000 cm⁻¹. Figure 4 shows a histogram of the relative signal-to-noise ratios (SNR) of CaF₂ vs. XT-KBr obtained from the 100% lines. This chart confirms the higher performance of CaF₂ above 5000 cm⁻¹.

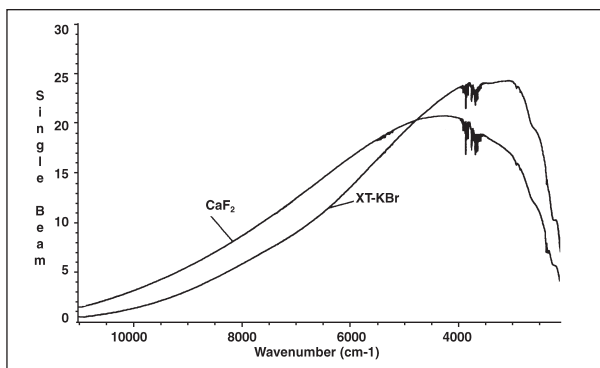


Figure 3

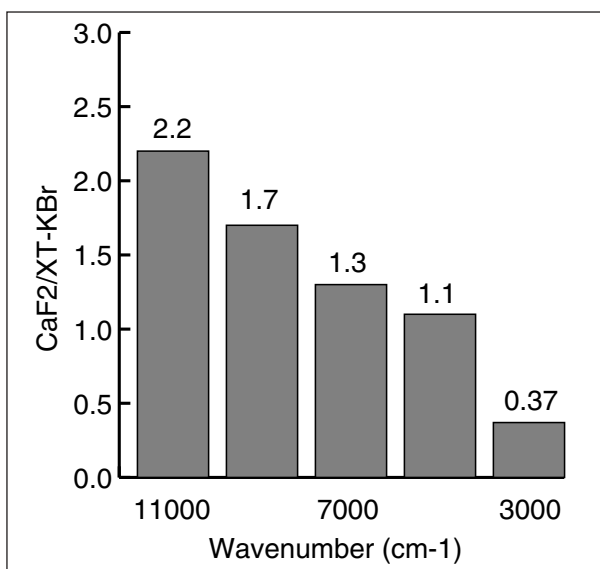


Figure 4

In the NIR range for the Raman shifted spectrum (9298-5668 cm⁻¹, representing a 3600-100 cm⁻¹ Raman shift), the signal-to-noise ratio for a cyclohexane spectrum is about 1.6 times better with the CaF₂ beamsplitter. However, the spectral quality shown in figure 5 is excellent for either choice of beamsplitter. The vast majority of FT-Raman sampling can easily be performed using the XT-KBr beamsplitter.

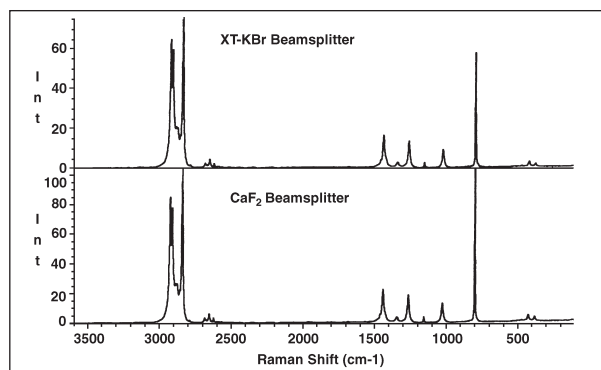


Figure 5

NIR fiber optic sampling is another technology for simplification of laboratory data collection. The sampling advantages include:

- No sample preparation
- Remote sampling
- Reflective samples may be housed in glass containers
- Very fast analysis times (10 seconds is typical)

NIR fiber optics sampling accessories are available both coupled to a sample compartment fiber-launcher or for mounting onto the face of the Nicolet 6700 FT-IR. There are also two accessories – the Smart NIR UpDRIFT and the Smart NIR Integrating Sphere – which fit into the sample compartment and provide enhanced NIR sampling capabilities.

The Thermo Scientific Nicolet SabIR™ NIR fiber optic sampling accessory was used to measure the spectrum of an analgesic tablet using the XT-KBr beamsplitter. The spectrum shown in figure 6 was produced in 20 seconds at 16 cm⁻¹ resolution.

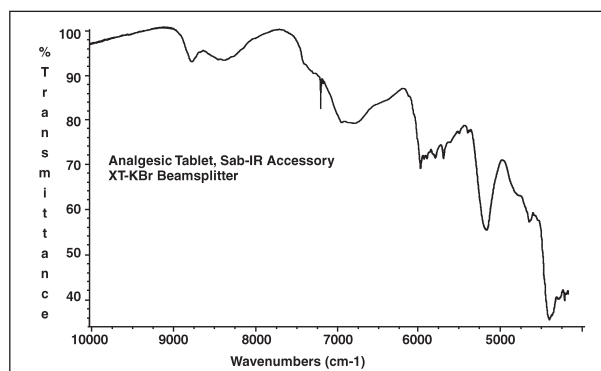


Figure 6

The XT-KBr beamsplitter also performs very well in the MIR. Figure 7 compares single beam spectra for the KBr and XT-KBr beamsplitters using the Ever-Glo ETC MIR source and DTGS detector (common scaled). These curves show that above 2600 cm^{-1} , the XT-KBr has greater throughput than KBr. Below 2600 cm^{-1} , the situation is reversed.

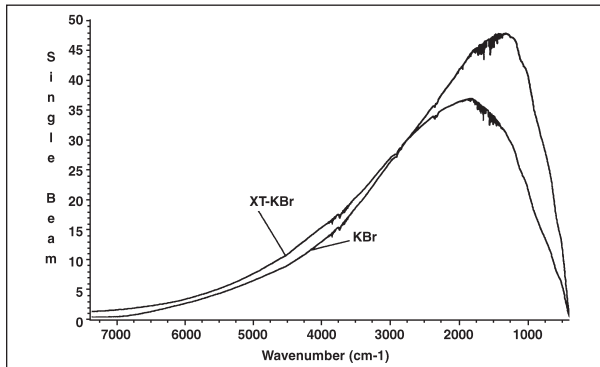


Figure 7

A numerical comparison of the MIR signal-to-noise ratios (SNR) for the two beamsplitters is shown in figure 8. At 2600 cm^{-1} , the SNR is identical. At 7000 cm^{-1} , the XT-KBr beamsplitter outperforms KBr by 4.6 times. At 450 cm^{-1} , the KBr beamsplitter outperforms the XT-KBr by about 2 times. The SNR of the Nicolet series of FT-IR spectrometers is significantly greater than most applications require, so the XT-KBr beamsplitter will provide reasonable performance for MIR spectroscopy over its full operating range.

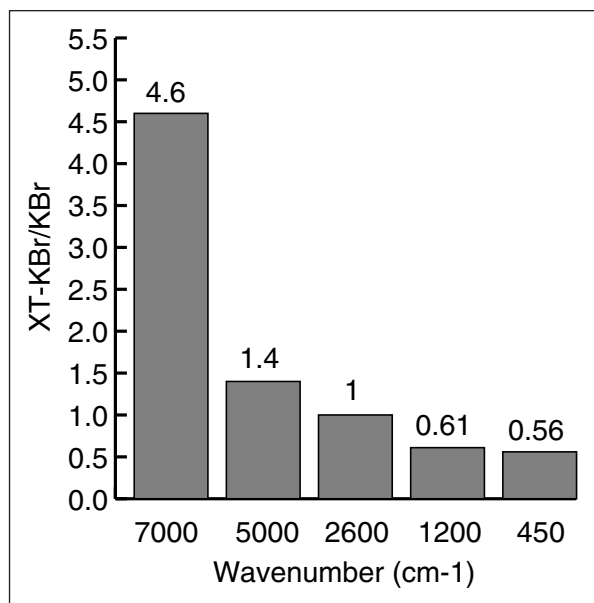


Figure 8

Conclusions

The XT-KBr beamsplitter provides the ability to collect mid and near-IR spectral data from 11,000 to 375 cm^{-1} . As with any spectroscopic solution, there are performance compromises over subsets of this range, but most routine mid-IR and near-IR applications can be easily done with the XT-KBr beamsplitter. The simplicity and time savings provided by the use of a single beamsplitter to address two spectral ranges is a significant advantage.

In addition to these offices, Thermo Fisher Scientific maintains a network of representative organizations throughout the world.

Africa
+43 1 333 5034 127

Australia
+61 2 8844 9500

Austria
+43 1 333 50340

Belgium
+32 2 482 30 30

Canada
+1 800 530 8447

China
+86 10 8419 3588

Denmark
+45 70 23 62 60

Europe-Other
+43 1 333 5034 127

France
+33 1 60 92 48 00

Germany
+49 6103 408 1014

India
+91 22 6742 9434

Italy
+39 02 950 591

Japan
+81 45 453 9100

Latin America
+1 608 276 5659

Middle East
+43 1 333 5034 127

Netherlands
+31 76 579 55 55

South Africa
+27 11 570 1840

Spain
+34 914 845 965

**Sweden/Norway/
Finland**
+46 8 556 468 00

Switzerland
+41 61 48784 00

UK
+44 1442 233555

USA
+1 800 532 4752

www.thermo.com



Thermo Electron Scientific Instruments LLC, Madison, WI USA is ISO Certified.

TN51432_E 04/08M

©2008 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.

Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.