### thermo scientific

## Step up

productivity, confidence, and compliance for pesticides analysis with a sample-to-result LC-MS/MS solution

Thermo Scientific Pesticide Explorer



Pesticide Explorer workflow: productivity, robustness, and regulatory compliance

The Thermo Scientific<sup>™</sup> Pesticide Explorer workflow is an easy to implement sample-to-result, LC-MS/MS analytical workflow, tailored to deliver robust, sensitive quantitation in compliance with global regulatory requirements. With a pre-tested and validated multi-residue pesticide method for over 400 analytes in different matrices, the workflow allows food-safety monitoring and testing laboratories to step up their productivity with reduced startup time and cost, easier compliance, and robust performance. Regardless of laboratory expertise, the Pesticide Explorer workflow is designed to simplify the process of obtaining reliable, unambiguous, high-quality LC-MS/MS results.

#### **PESTICIDE EXPLORER**



## Achieve immediate success with a comprehensive, single-provider workflow

The Pesticide Explorer workflow is an optimized single-provider solution, built by scientists for scientists working in food safety laboratories around the world. The comprehensive solution provides a reliable, high-quality LC-MS/MS workflow for analyzing hundreds of pesticide residues in a wide range of matrices.

#### Meet all requirements for robust multi-class pesticide residue analysis

Laboratories face numerous challenges in developing robust, sensitive quantitation workflows for multi-class pesticide residue analysis. Adding to the challenge is the difficultly of reducing cost per sample while obtaining high performance and certainty in results. The Pesticide Explorer workflow solves these challenges with a complete solution from sample to results, optimized to ensure successful analysis of pesticide residues in a variety of food matrices.

The workflow includes the TSQ Quantis triple-stage quadrupole mass spectrometer, renowned for excellent sensitivity, robustness, and speed, enabling laboratories to meet current and future regulations for multi-class pesticide residue analysis. The immediate benefits for laboratories are:

- Rapid start-up with a comprehensive solution—Thermo Scientific<sup>™</sup> HyperSep<sup>™</sup> Dispersive SPE (QuEChERS) sample preparation kit, quality control (QC) text mix, industry-leading TSQ Quantis triple-stage quadrupole mass spectrometer, state-of-the-art Vanquish Flex Binary UHPLC system, Thermo Scientific<sup>™</sup> Accucore<sup>™</sup> aQ columns, TraceFinder software methods, and detailed user guide—tailored to meet the most stringent pesticide residue regulations for food safety analyses
- **Productivity** with a ready-to-run workflow that empowers users of any expertise to analyze over 400 multi-class pesticide residues in a single 15-minute analysis
- **Confidence** with robust, pre-tested, and validated multi-residue pesticide data acquisition and processing methods and extensive pesticide database of over 600 compounds with selected reaction monitoring (SRM) transitions and retention times
- **Compliance** with European Union SANTE Guidelines for pesticide residues in food for ultimate confidence in results





Extracted ion chromatogram of over 200 pesticide residues in baby food at 10 ng/g shown in TraceFinder software



Vanquish Flex Binary UHPLC system with TSQ Quantis triple-stage quadrupole mass spectrometer

### Pesticide Explorer workflow components

- **QuEChERS** (Quick, Easy, Cheap, Effective, Rugged and Safe) sample preparation reagent slim pouch kit with full procedures
- Quality Control (QC) standards for performance check
- Vanquish Flex Binary UHPLC system
- Accucore aQ columns
- TSQ Quantis triple-stage quadrupole mass spectrometer
- **TraceFinder software** with pre-tested, validated data acquisition and processing methods, along with a comprehensive compound database
- Detailed user guide

#### Convenient, easy, and effective: QuEChERS sample preparation

The QuEChERS sample preparation kit provides a time-saving yet extremely effective approach for the extraction and cleanup of pesticide residues in food matrices. Tested sample preparation procedures are included for wine, olive oil, and baby food matrices. With the QuEChERS kit, users can easily extract and analyze samples with confidence.



QuEChERS sample preparation method: proven, simple, and easily transferrable

#### Benefits of the QuEChERS sample preparation method

Characteristic	QuEChERS method						
Variability	LOW: proven reproducibility for hundreds of pesticides across multiple matrices						
Lab portability	HIGH: one method can be easily implemented and transferred in multiple labs and geographies						
Ease of use	HIGH: easy-to-use configurations and methods allow for simple workflow implementation						
Cost per sample	LOW: economical, users with varied expertise achieve similar results						

#### Productivity and quality: Vanquish Flex Binary UHPLC system

With excellent retention time stability and maximum pressure up to 1000 bar, the Vanquish Flex Binary UHPLC systems increase productivity without compromising quality of results. The system shares all Vanquish system values such as a design focused on uptime, robustness, and quality, enabling laboratories to address any quantitation challenge.

#### Speed and resolution: Accucore aQ columns

The high-throughput, high-resolution Accucore aQ columns are ideally suited for pesticide-residue monitoring. The rugged solid-core column with polar end-capped C18 chemistry delivers excellent reproducibility for the polar and hydrophobic analytes typically found in multi-class, multi-residue pesticide analyses. When combined with the power of the Vanquish Flex Binary UHPLC system, users benefit from the extended pressure capabilities and robustness required for consistent, high-throughput separations.

#### Confident quantitation: TSQ Quantis mass spectrometer

With robust analysis of pesticide residues in complex matrices, the TSQ Quantis triple-stage quadrupole mass spectrometer enables food-safety testing labs to address regulatory requirements with confidence and ease. Users become productive quickly with built-in automated compound optimization and application-specific methods and reports. Outstanding scan speed permits simultaneous quantitation of hundreds of compounds at regulatory limits.



TSO QUANTIS

Matrix extracted standard (MES) of Azoxystrobin in wine matrix at a concentration of 1.0 ng/g, with the quantitation ion peak (left), qualitative confirming ions (middle), and calibration curve range from 0.5 to 100 ng/g (right)

#### Five clicks to success: TraceFinder software

TraceFinder software includes an extensive built-in compound database; the ability to screen and quantify analytes; streamlined data review; customized, editable report templates; and automatically generated reports—specifically tailored for food safety applications. From data acquisition to the final report, TraceFinder software makes workflows fast and simple.





Matrix-extracted standard (MES) of difenacoum at 0.5 ng/g in wine matrix showing the quantitation ion (left) and the overlay with the qualifier ions (right) in TraceFinder software



# Become more productive with a ready-to-run TraceFinder software workflow

Laboratories can substantially reduce time for high-confidence results using LC-MS/MS pesticide residue analysis methods that are included with the workflow. For ultimate ease of use, all necessary LC-MS/MS instrument parameters are specified. Users simply load and launch a method. If desired, laboratories can easily develop customized methods using the TraceFinder software compound database.

und Database PE2 Full C	MPD ISI															
Tang v V X	Peak View Pane			_		_							-		_	
Name od Al. Collacer Al.	# Compound Name	2 Peak Label 2	Peak Morkflow		lasociated larget Peak		- MS Order		Precursor m/2	- Product m/2	w(t	- Adduct		Polarity	- Oage	
Post 1	118 apineutluron	T1C2-425109->226	Contaming	• 1	2 425 109->13		ms2		425.309	226.970	425.109	None		Positive	* 1	
24-dimethylaniline 26-dichtorsbercamide	119 apinghos-ethyl	T1: 346.044 + 76.917	TargetPrais				m2		345.044	75.917	345.044	Nord		Positive	- 1	
2-methol 4-6-dinitrophen	120 scinphos-sthul	T1C1 345 544->131	Conferring		2 345 544-+75		ma2		345.044	131.917	345.044	None		Positive		
145-binethacalb	121 scinghos-ethyl	T1C2 345 644->260	Conferring		3 345 544->36		ms2		345.044	250.687	346.044	None		Positive		
6-chloro-4-hydroxy-3-phe abarrel fo	122 approtone	T2: 226-085-+124.92	TargetPark				ms2		225.066	124.929	225.085	None		Positive	- 1	
abarrectin acrphete	123 adjorothme	T1C1 225 585->155	Contemino	- 1	2 225 585->12	( - L	ma2	-	225.066	155.917	226.085	None	-	Positive		
acetamiprid acetochior	124 approxime	T1C2: 226.086->197.	Contaning	- 1	1 225.045->12		ms2		225.066	197,657	226.086	Nore		Positive	- 1	
aciterzolar-s-methyl	125 Appointing	T2:400195->178.06	TarpetPark				ms2		420.225	178.042	400195	None		Posting		
adonies according AHA	126 Assoyciation	T1C1-400.195->248	Conferring		1.400195->17		ms2		420.295	248.917	430195	None		Positive		
election albendapole	127 Azocyclotie	T1C2-400.195->274	Conferring		3 430195->17		ms2		430.285	274.687	430195	Nore		Postar		
aldica-b	128 appropriate	T1:404.124->328.65	TargetPask				ms2		404.124	328.657	404124	None		Positive	* 1	
aldcarb-sulfone-NH4+ aldcarb-sulfoxide	129 appropriate	T1C2-404.124-+348	Confirming	• 1	2-404124->32		ms2		404.124	343.929	404.124	None		Positive		
alettria	130 accostration	T1C2-404124->371	Conferring		2-454.124->22		m42		404.124	371,929	404124	Nace		Positive		
alldochlor Alloxidm sodio	131 beflubutered	T1: 256125->64.857	TargetPrais				ms2		256.126	64.857	256.126	None		Postar	- 1	
alpha-cypermethrin-NH4 ametochradin	122 behubutamid	T1C1-256.126->90.8	Contaming	1.1	2 255.128->64		ma2	12	255.126	90.845	155.125	None		Postive		
arrelyn	122 befullutarrid	TLC2 255.126->161	Conferring		2 255 125->64		m42		256.126	151,929	256.126	None		Positive		
amicarbacone Amicultorom	134 bendavid	T1 225175->148.00	TargetProj				ms2	1.0	325.175	146,000	226175	Name		Poster		
anipole anomidal	135 benalaql	T1C1 335.175->307	Confirming	• 1	1 325175->14		ma2		326.175	207.929	126175	None		Positive	- 1	
anlasine																- 1
aniotos2 aramite-NH4	Compound Details Pane															1 × 1
apon autor	Compound Name															
abaton																- 1
rhapine ahapine-deseth/	Jonization	Note	-													- 1
Attacine-deckopropyl	Chemical Formula			N	autral Mass   0											- 1
azaconazole azamethiphos	CAS No.				Category -											- 1
scinesulturon acimphosi ethyl	Compound Type	Analis		britem.	a Standard [											- 1
siprotyne	Compound Groups	wate														- 1
Azocyclotin azonystrobin	Compound Groups															- 1

Salact compounds

Choose from over 600 pesticides provided in the TraceFinder Compound Database to create a custom method, or choose one of the pre-tested, validated Master Methods for residue analysis.



LC-MS/MS method flexibility allows users to easily edit the pre-configured methods, including the SRM transitions and retention times, if required. Visualization features in the instrument software display dwell times and SRM density throughout the entire analytical run.

#### Acquire data



Data is acquired using optimal LC-MS/MS conditions for each multi-class, multi-residue method to deliver superb quantitative performance and sensitivity.



#### Review data



The color-flagging feature of TraceFinder software enables analysts to quickly review and confirm results.

#### Generate report



After data review, high-quality standard or custom reports can be used to rapidly generate useful results.



LC-MS/MS extracted ion chromatograms of white wine spiked with selected pesticides (overlay of injection #1 and injection # 300) demonstrate the robustness and reproducibility of the Vanquish Flex Binary UHPLC system and Accucore aQ column with the TSQ Quantis mass spectrometer.



The fast scan speed with polarity switching of the TSQ Quantis mass spectrometer provides accurate quantitation of residues in large pesticide panels. Over 500 pesticides were analyzed in a wine matrix over a 15-minute run. The inset shows the scans across the target pesticide, methabenzthiazuron, at 10 ng/g.

## thermo scientific

## The most comprehensive portfolio for pesticide analysis

In addition to the Pesticide Explorer Workflow, we offer complementary technologies for the analysis of pesticide residues. While each solution has specific advantages, collectively Thermo Scientific offers the most comprehensive pesticide analysis portfolio available on the market today.



## Orbitrap mass analyzer based for GC-MS, LC-MS and MS/MS

The Thermo Scientific<sup>™</sup> Orbitrap<sup>™</sup> mass spectrometers produce high-resolution accurate-mass (HRAM) data for confident trace-level screening and quantitation of pesticides, and identification of unknowns. Orbitrap-based instruments are renowned for high quality full-scan and MS/MS data with easy analytical set-up, efficient data processing, and retrospective data analysis.

#### Ion chromatography-MS/MS

For polar anionic pesticides that are challenging to analyze using conventional or hydrophilic interaction chromatography (HILIC) LC-MS/MS, the Thermo Scientific<sup>™</sup> Dionex<sup>™</sup> Integrion<sup>™</sup> HPIC system coupled with the Thermo Scientific<sup>™</sup> TSQ Altis<sup>™</sup> triple-stage quadrupole mass spectrometer provides a powerful solution. High-pressure ion chromatography (HPIC)-MS/MS with high-capacity ion exchange columns and post-column eluent suppression, combined with highsensitivity triple quadrupole mass spectrometer detection improves sensitivity and chromatographic separation for high-confidence analysis of polar anionic pesticides in complex food matrices.

#### Find out more at thermofisher.com/PesticideExplorer

For Research Use Only. Not for use in diagnostic procedures. © 2019 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. This information is presented as an example of the capabilities of Thermo Fisher Scientific Inc. products. It is not intended to encourage use of these products in any manner that might infringe the intellectual property rights of others. **BR73188-EN 0919M** 



## Chromatography Data System (CDS) software

Simplified data processing and extensive multi-vendor instrument control make the Thermo Scientific<sup>™</sup> Chromeleon<sup>™</sup> Chromatography Data System the choice for many laboratories. The software streamlines workflows from chromatography to routine quantitative MS analysis on IC, GC, LC, GC-MS/MS and LC-MS/MS systems by delivering superior instrument control, automation, data processing, and more. Run analyses in an enterprise environment—from method creation to final reporting—to boost productivity.

