


Non-Sterile Samco™ Capillary Transfer Pipette


INSTRUCTIONS FOR USE (thermofisher.com/capillarytransferpipettes)

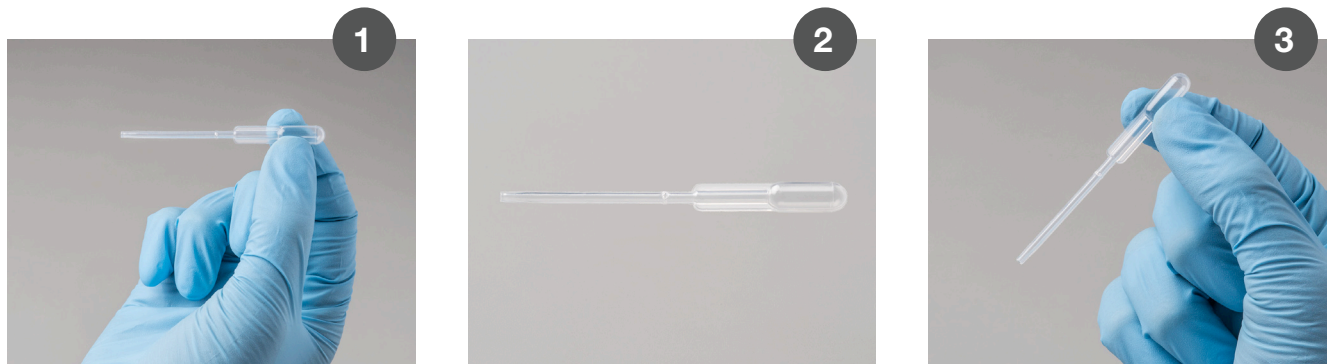

1. Hold horizontally, contact tip to droplet. **DO NOT SQUEEZE BULB, DRAW IS AUTOMATIC.** Angle the bulb down if slow draw.
2. Draw complete when draw stops and fluid reaches indicator band.
3. Squeeze bulb to dispense within 30 seconds of collection.
4. Discard used capillary transfer pipette according to your facilities disposal procedures.

CAUTION:

- This product is intended to be used for collection and transfer of liquids.
- Single use only.
- Recommend use with gloves to help reduce contamination risk.
- Do not clog vent hole.
- Do not rinse or soak in any solution.
- Not for long term sample storage.

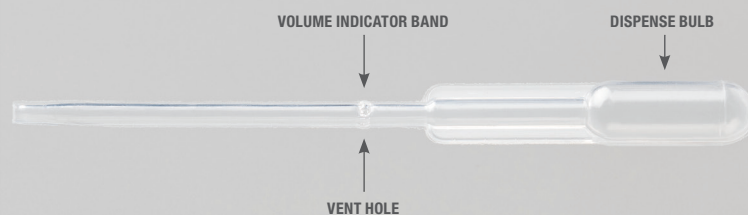

LABOMEX MBP, S. DE R.L. DE C.V.
Thermo Fisher Scientific
Braulio Maldona No. 6470
Tijuana, Baja California, CP 22626, Mexico

Made in Mexico

