

Thermo Scientific KingFisher Pure Kits



Pure performance
every time

Thermo
SCIENTIFIC

Thermo Scientific KingFisher Pure Kits

pure performance every time

We've expanded our unique nucleic acid purification workflow, providing optimized high-throughput methods and extreme flexibility!

- Flexible KingFisher system with eight optimized purification kits
- Dedicated kits for RNA or DNA in a wide variety of sample types
- Excellent yield, purity and reproducibility
- Walk-away solution for any throughput

Highly Reproducible Purification

The Thermo Scientific™ KingFisher™ system offers a rapid, automated and reproducible purification workflow, by using magnetic rods to move particles through the purification phases of binding, washing and elution. After sample lysis, nucleic acids bind efficiently with coated magnetic beads. Washing the nucleic acid/bead complex several times eliminates contaminants, producing highly pure DNA or RNA.

By eliminating laborious manual steps and purification problems such as filter clogging and risk of contamination, the KingFisher system consistently produces high purity yields.

Flexible Throughput

The KingFisher system can easily be scaled-up in terms of sample input volume and throughput – from 1 to 96 samples per run. Larger sample input volumes, for example up to 1 mL of blood, can be processed by the Thermo Scientific KingFisher Flex and KingFisher Duo instruments.

Minimized Hands-on Time

High sample purity and yields are obtained with minimal user input, and the maximized walk-away time ensures efficiency.

Safe and Easy

The easy-to-use KingFisher system and straightforward Thermo Scientific BindIt software protocols facilitate safe sample handling.

Optimized KingFisher Pure Kits and Protocols

Thermo Scientific KingFisher Pure Kits and protocols have been optimized with KingFisher systems for the rapid purification of the highest quality and yields of DNA or RNA, compatible with most common enzymatic downstream analyses and applications. There is no need for organic solvents.



Premium Quality Beads


The KingFisher Magnetic Beads are characterized by high surface area, uniform size distribution and excellent suspension stability, making them ideal for high throughput processing.

Wide Range of Samples

From routine use to advanced requirements, KingFisher Pure Kits are available for both DNA and RNA purification in a broad variety of sample types, including blood, serum, cells, tissues and plant material.

Excellent Performance


All eight customized KingFisher Pure Kits produce nucleic acids of high yield and quality. Kits are optimized for:

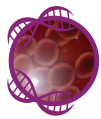

Blood DNA



Plant DNA

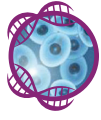


Tissue DNA



Plasmid DNA



Viral DNA/RNA

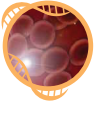

Blood RNA



Plant RNA



Tissue RNA

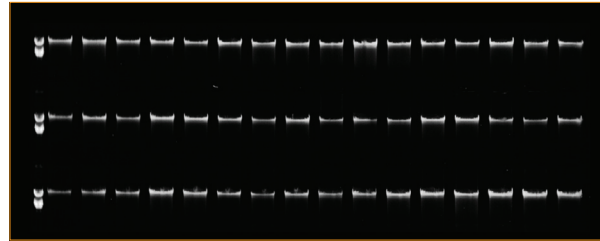


Figure 1. Genomic DNA purified from 200 µl of blood using the KingFisher Pure DNA Blood Kit shows excellent quality of gDNA on the agarose gel.

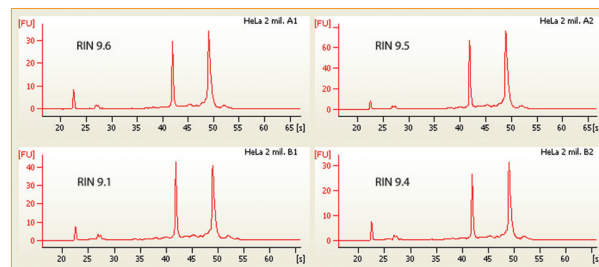


Figure 2. Total RNA purified from 2×10^6 HeLa cells using the KingFisher Pure RNA Tissue Kit shows intact RNA when analyzed on the Agilent Bioanalyzer 2100.

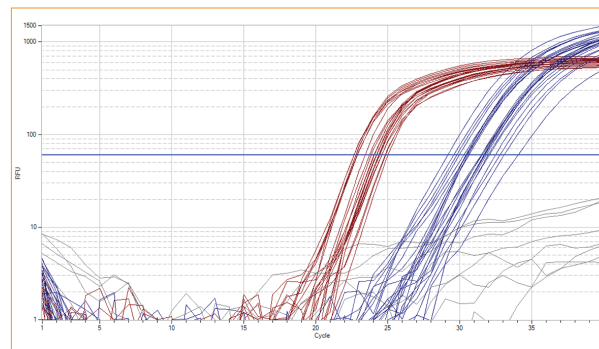


Figure 3. The RNA purified from 200 µl blood using KingFisher Pure RNA Blood Kit was used as a template in a gene expression study with Thermo Scientific PikoReal Real Time PCR System. Reference gene (GAPDH) in red and target gene in blue (PPIH).

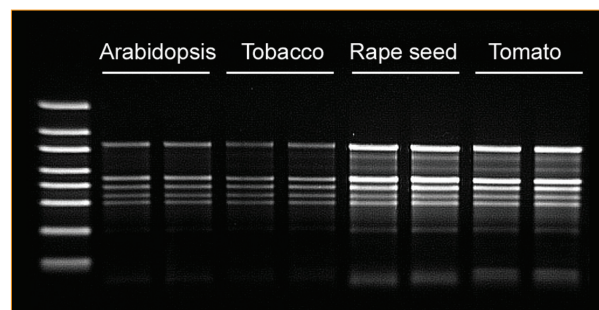


Figure 4. Total RNA purified from 15 mg plant leaves using the KingFisher Pure RNA Plant Kit shows intact RNA on the agarose gel.



Specification and Ordering Information

Product #	Description	Package size	Sample type	Sample input	Elution volume* (µL)	Run time (min)
98010196	KingFisher Pure DNA Blood Kit	96	Fresh or frozen whole blood treated with EDTA or citrate	200 µL or 1 mL**	200 µL: 100 µL (Duo) or 150 µL (Flex). 1 mL: 500 µL	40 (200 µL), 45 (1 mL)
98010496	KingFisher Pure DNA Blood Kit	384				
98020196	KingFisher Pure RNA Blood Kit	96	Fresh or frozen whole blood treated with EDTA or citrate, PaxGene and Tempus preserved blood	200 µL	100 µL (Duo) or 150 µL (Flex)	60
97030196	KingFisher Cell and Tissue DNA Kit	96	Tissues, cultured cells or bacteria	Up to 20 mg of tissue, 1 x 10 ⁷ cells, or 1 mL of overnight cultured bacteria	100 µL (Duo) or 150 µL (Flex)	25
97032496	KingFisher Cell and Tissue DNA Kit	24 x 96				
98040196	KingFisher Pure RNA Tissue Kit	96	Tissues or cultured cells	Up to 20 mg of tissue or 2 x 10 ⁶ cells	100 µL	50
98040496	KingFisher Pure RNA Tissue Kit	384				
98050196	KingFisher Pure DNA Plant Kit	96	Fresh or frozen plant tissues	Up to 50 mg of fresh plant tissue	100 µL (Duo) or 150 µL (Flex)	40
98050496	KingFisher Pure DNA Plant Kit	384				
98060196	KingFisher Pure RNA Plant Kit	96	Fresh or frozen plant tissues	Up to 50 mg of fresh plant tissue	100 µL	60
98060496	KingFisher Pure RNA Plant Kit	384				
98070196	KingFisher Pure Viral NA Kit	96	Cell-free body fluids, such as serum and plasma	200 µL	100 µL	40
98070496	KingFisher Pure Viral NA Kit	384				
98080196	KingFisher Pure Plasmid Kit	96	Bacterial culture	Up to 5 mL	100 µL	40
98080496	KingFisher Pure Plasmid Kit	384				

*Recommended elution volumes, but user adjustable

**With 1 mL samples reagents are sufficient for approx. 25 samples with 96 kit and for approx. 100 samples with 384 kit.

thermoscientific.com/kingfisherinfo

© 2013 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

North America:

USA/Canada +1 603 595 0505

USA toll free 800 345 0206

Europe:

Austria +43 1 801 40 0

Belgium +32 53 73 42 41

Finland +358 9 3291 0200

France +33 2 2803 2000

Germany national toll free 08001-536 376

Germany international +49 6184 90 6940

Italy +39 02 95059 552

Netherlands +31 76 571 4440

Nordic/Baltic countries +358 9 329 100

Russia/CIS +7 (495) 739 76 41

Spain/Portugal +34 93 223 09 18

Switzerland +41 44 454 12 12

UK/Ireland +44 870 609 9203

Asia:

Australia +613 9757 4474

China +86 21 6865 4588 or +86 10 8419 3588

China toll free 800-810-5118, 400-650-5118

India +91 22 6716 2200

Japan +81-3-5826-1616

Korea +82 11 796 7771

Other Asian countries +65 6872 9717

Countries not listed:

+49 6184 90 6940 or +33 2 2803 2000

Thermo
SCIENTIFIC

Part of Thermo Fisher Scientific