For a complete user manual and additional resources including installation videos for Connect Edge, visit www.thermofisher.com/connectedge

If additional assistance is needed, please contact Technical Support at +1 866 984 3766 (www.thermofisher.com/contactus)

Safety Considerations

IMPORTANT: Read carefully and understand the instructions and warnings contained in this document before installing or using the product. Keep this document for future reference. To lower the risk of personal injury, electric shock, fire, or damage to equipment, observe the instructions and warnings contained in this document. Failure to comply with the instructions and warnings contained in this document violates the standards of safety, design, manufacture, and intended use of the product. Thermo Fisher assumes no liability for any damage caused by failure to observe the instructions and warnings contained in this document. For concerns regarding the correct understanding of this document, contact Thermo Scientific technical services.

Device Maintenance

CAUTION: Do not attempt to disassemble the device. There are no user serviceable parts inside.

CAUTION: Do not misuse the device. Follow instructions on proper operation and only use as intended. Misuse could make the device inoperable, damage the device and/or other equipment or harm users.

CAUTION: Never try to dismantle the product yourself or push the objects of any kind into the product, as this may cause short circuits which could result in a fire, electric shock and injury up to and including death.

CAUTION: Do not apply excessive pressure or place unnecessary weight on the device. This could result in damage to the device or harm to users.

CAUTION: Do not use this device in explosive or hazardous environments.

CAUTION: Do not expose your device to any extreme environments such as those with high temperature or humidity. Such exposure could result in damage to the device or fire.

CAUTION: Do not expose the device to liquids. It is not waterproof. Exposure to liquids could result in damage to the device, electric shock or injury up to and including death.

CAUTION: Do not place the device alongside computer discs, credit or travel cards or other magnetic media. The information contained on discs or cards may be affected by the device.

CAUTION: Using accessories, such as antennas, that Thermo Fisher has not authorized or that are not compliant with Thermo Fisher accessory specifications may invalidate the warranty.

Caution: Changes or modifications to this unit not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

Safety Instructions

- Do not use this product for protection or as part of an automated emergency system or for any other application that involves protecting people and/or property.
- Customers and users of Thermo Fisher Scientific products are responsible for making sure that the product is fit for the intended usage. Do not open the product casing and do not disassemble or modify internal components in any manner.
- Thermo Fisher Scientific products do not contain any internal components that require user intervention or repair. If the device shows signs of improper operation, disconnect it immediately from its power source and contact Thermo Fisher Scientific technical support.

Electrical Warnings

- The Connect Edge gateway must be powered by a 24V DC power supply.
- The Connect Edge adapters must be powered by a 5V DC power supply or via the PoE Ethernet base.
- Do not open the Connect Edge gateway and adapters and do not dismantle internal components or modify them in any manner. The Connect Edge gateway and adapters do not contain any user-repairable parts. If the Connect Edge gateway and adapters show any sign of malfunction, unplug them immediately and contact Thermo Scientific for repair or replacement.
- Do not cause a short circuit with the electrical plug. Do not force either the AC or DC plug.
- Before removing the connector from the gateway or unplugging power cables, unplug the cable from the power outlet.
- Do not subject the gateway and adapters to physical shock, as it may cause serious malfunction or damage. Do not use or place the gateway and adapters in a wet or humid location. The Connect Edge gateway and adapters are not waterproof.

Maintaining your Connect Edge Hardware

Maintaining your gateway

Periodically inspect the product to verify its integrity and to ensure proper operation. To maintain the product, complete the following steps:

- Carefully read and understand the instructions contained in this document
- Safely remove the power supply.
- Verify the installation of the product.
- Clean the product.

How to safely remove the power supply

ELECTRIC SHOCK HAZARD

Failure to remove power correctly may create an electric shock hazard, which could result in personal injury or loss of life, and / or damage the equipment or other property.

To avoid injuries and safely remove power supply from the product, complete the following steps:

- Make sure your hands are dry.
- Turn OFF all the power supply sources.
- Disconnect all the cables.
- Make sure that all the circuits are discharged.

How to verify the installation of the product

To verify the installation of the product, complete the following steps:

- Verify that the product is clean and not damaged.
- Verify that the LED indicators are visible and not damaged.
- Verify that all the locking parts (for example: screws, bolts, nuts) are correctly fastened.
- Verify that the product is installed correctly.

How to clean the product

To clean the product, complete the following steps:

- Never use detergents, aerosol sprays, solvents, or abrasive sponges.
- To remove dust from the case of the product, use a dry, lint-free, cloth.
- To remove the dirt, use water-based, non-flammable, cleaner products.

CAUTION: Follow your facility safety guidelines and wear the required Personal Protective Equipment (PPE) during installation and/or troubleshooting.

Regulatory Information

Certifications and Compliance

FCC statements

This device complies with part 15 of the FCC rules. Operation is subject to the following two conditions:

1. This device may not cause harmful interference
2. This device must accept any interference received, including interference that may cause undesired operation

FCC RF Radiation Exposure Statement

This equipment complies with FCC radiation exposure limits set forth for an uncontrolled environment. To avoid the possibility of exceeding the FCC radio frequency exposure limits, human proximity to the antenna should not be less than 20 cm (8 inches) during normal operation. End users must follow the specific operating instructions for satisfying RF exposure compliance. This transmitter must not be co-located or operating in conjunction with any other antenna or transmitter.

FCC Class A Notice

This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause interference to radio communications. Operation of this equipment in a residential area may cause interference in which case the user will be required to correct the interference at their own expense.

IC Statements

This device complies with Industry Canada license-exempt RSS standard(s). Operation is subject to the following two conditions:

1. This device may not cause harmful interference
2. This device must accept any interference, including interference that may cause undesired operation of the device

Under Industry Canada regulations, this radio transmitter may only operate using an antenna of a type and maximum (or lesser) gain approved for the transmitter by Industry Canada. To reduce potential radio interference to other users, the antenna type and its gain should be so chosen that the equivalent isotropically radiated power (e.i.r.p.) is not more than that necessary for successful communication.

This Class A digital device complies with Canadian ICES-003.

CE - Conformity with European regulations

The product is in compliance with the essential requirements and other relevant provisions of directive 2014/53/EU (Radio Equipment Directive), 2014/30/EU (EMC Directive) and 2014/35/EU (Low Voltage Directive).

RoHS Statement

The wireless device complies with the EU Directive 2011/65/EU (Restriction of the Use of Certain Hazardous Substances in Electronic and Electrical Equipment) and the amendment of (EU) 2015/863. Do not dispose of this product with household trash. Thermo Scientific recycles this product under certain conditions. Contact us for more information.

CAUTION: Any changes or modifications not expressly approved by Thermo Scientific could void the user's authority to operate the equipment.

South Korea: (Applicable for the gateway)

Statement for class A equipment

Class A equipment (Broadcasting equipment for business) This equipment may cause radio interference when used in household environment certified for business use purpose.

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Statement for KN 32/35

This product may cause malfunction if it is used nearby wireless communication equipment such as a cell phone, a Wi-Fi or Bluetooth device etc.

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