

Large Volume Precipitation of Proteins with Ammonium Sulfate Using Thermo Scientific Fiberlite Carbon Fiber Rotors

Key Words

Ammonium Sulfate Precipitation, Superspeed Centrifuges, Carbon Fiber Rotors

Introduction

Ammonium sulfate precipitation is a method of protein purification by altering the solubility of protein. Ammonium sulfate is commonly used due to its high solubility that allows salt solutions with high ionic strength.

At sufficiently high ionic strength, the protein will be almost completely precipitated from the solution. Differential precipitation of proteins by ammonium sulfate is one of the most widely used preliminary purification procedures. It is based on proteins having differing solubility in ammonium sulfate solutions and can result in a two- to five-fold increase in specific activity.

Provided that appropriately buffered ammonium sulfate solutions are used to protect the desired activity, recoveries approaching 100% can be expected. A typical protocol consists of adding ammonium sulfate to give specific percentage saturation, followed by a period of time for proteins to precipitate and a centrifugation step to collect the precipitate. The adapted protocol¹ described below was carefully followed to precipitate protein removed by centrifugation using the Thermo Scientific Sorvall LYNX 6000 superspeed centrifuge and the large volume Thermo Scientific Fiberlite F12-6x500 LEX carbon fiber rotor. The total capacity of the Fiberlite® F12-6x500 LEX rotor is 3 liters and is rated for RCF up to 24,471 x g.

The ammonium sulfate concentration was increased to a value that will precipitate most of the protein of interest, while leaving the maximum amount of protein contaminants still in solution. The precipitated protein of interest was recovered by centrifugation and dissolved in fresh buffer for the next stage of purification.

PROTOCOL: Ammonium Sulfate Precipitation

1. Measure the volume of protein solution to obtain 1500 mL. If necessary add 1 M HEPES, pH 7.4, or 1 M Tris-HCl, pH 7.5, to a final concentration of 50 mM.



Figure 1: Thermo Scientific Fiberlite F12-6x500 LEX carbon fiber rotor.

2. While gently stirring, slowly add ammonium sulfate to 45% saturation (277 g ammonium sulfate/liter solution). Stir gently for a minimum of 4 hr and a maximum of 16 hr at 4 °C.

Note: Addition of ammonium sulfate acidifies the solution, therefore protein stability may be improved by buffering with at least 50 mM HEPES or Tris buffer. It is important that no foaming of the solution occurs during ammonium sulfate addition as this will cause denaturation of proteins. If foaming occurs, slow the addition of ammonium sulfate. Make sure that the ammonium sulfate dissolves nearly as quickly as it is added.

3. Transfer the ammonium sulfate precipitated solution to 500 mL Thermo Scientific Nalgene polycarbonate centrifuge bottles with screw caps (PN 3140-0500).

Perform centrifugation using the Fiberlite F12-6x500 LEX rotor and the Sovall® LYNX 6000 superspeed centrifuge with the following parameters: 25,000 x g for 20 min at 4 °C.

4. Decant the supernatant, being sure not to disturb the relatively loose protein-ammonium sulfate pellet.

Note: At this point, the pellet can be stored for several weeks at 4 °C without significant loss of protein stability. The supernatant can also be stored at 4 °C for later analysis.

5. Re-suspend the pellet in PBS, pH 7.2. The solution will remain turbid, but discreet flocculent is not visible.
6. Remove residual precipitate and collect the clarified proteins by performing centrifugation with the following parameters: 25,000 x g for 10 min at 4 °C.
7. Remove the clarified proteins to fresh tubes and store at 4 °C for up to one week.
8. Alternatively, if column purification is going to be performed immediately, dialyze protein against column binding buffer. Pass the solution through a 0.45 µm filter and apply to the appropriate column.

Conclusion

This protocol is useful to quickly remove large amounts of contaminant proteins, as done in the first step in many purification schemes. It is also often employed during the later stages of purification to concentrate protein from a dilute solution following procedures such as gel filtration.²

In addition to the rapid precipitation of protein contaminants, the Fiberlite F12-6x500 LEX rotor will harvest bacteria and yeast from cultures in shorter run times than rotors with similar total volumes.

References

1. Elder, G.H., Tovey, J. A. and D. A. Sheppard, (1983). Purification of uroporphyrinogen decarboxylase from human erythrocytes. Immunochemical evidence for a single protein with decarboxylase activity in human erythrocytes and liver. *Biochem. J.* 215: 45-55.
2. Koch, U., Choksi, S., Marcucci, L. and R. Korngold, (1998). A Synthetic CD4-CDR3 Peptide Analog Enhances Skin Allograft Survival Across a MHC Class II Barrier. *Journal of Immunology*, 161: 421-429.

thermoscientific.com/lynxsuperspeed

© 2012 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

Australia +61 39757 4300
Austria +43 1 801 40 0
Belgium +32 53 73 42 41
China +86 21 6865 4588 or
+86 10 8419 3588
France +33 2 2803 2180
Germany national toll free 0800 1 536 376
Germany international +49 6184 90 6000

India toll free 1800 22 8374
India +91 22 6716 2200
Italy +39 02 95059 554
Japan +81 45 453 9220
Netherlands +31 76 579 55 55
New Zealand +64 9 980 6700
Nordic/Baltic/CIS countries
+358 9 329 10200

Russia +7 812 703 42 15
Spain/Portugal +34 93 223 09 18
Switzerland +41 44 454 12 22
UK/Ireland +44 870 609 9203
USA/Canada +1 866 984 3766

Other Asian countries +852 2885 4613
Countries not listed +49 6184 90 6000

Thermo
SCIENTIFIC

Part of Thermo Fisher Scientific