



Aqua-Tec Water Preservation Cell

Protect your lab samples with confidence

Easily prevent microbial contamination from any CO₂ incubator or water bath reservoir without chemical disinfection solutions.

The Thermo Scientific™ Aqua-Tec™ Water Preservation Cell is a revolutionary way to protect your valuable samples from waterborne contaminants without the use of harsh chemicals. Simply place the 3,7-inch (~ 9,4 cm) cell into the water reservoir of any CO₂ incubator or water bath.* The Aqua-Tec cell prevents contamination from most types of bacteria, molds, and fungi for up to six months.

The Aqua-Tec Water Preservation Cell was developed specifically for laboratory applications across a range of temperatures. It's safe to handle, requires no special protocol for disposal, and enables virtually maintenance-free microbial protection. The Aqua-Tec cell automatically self-regulates to maintain proper disinfecting concentration and effectiveness, regardless of the water level.

Features and benefits

- **Designed and tested specifically for laboratory use** in CO₂ incubators and water baths
- **Proven effective for up to 6 months** against bacteria, molds, and fungi
- **Safe to handle**—requires no special disposal protocol
- **Self-regulates to maintain effectiveness**, regardless of water level
- **Easy to use** in equipment from any manufacturer

* Use only sterilized distilled water in CO₂ incubators.

Safe and effective for you and your research—helps protect your valuable samples

The Aqua-Tec Water Preservation Cell has been proven to inhibit the growth of a wide range of commonly encountered microbes for up to six months. Eliminate the need to handle dangerous chemicals or expose sensitive cells to the risk of potentially toxic offgassing from disinfectants. The Aqua-Tec cell allows your CO₂ incubators and water baths to operate at maximum efficiency and saves the significant time, handling, and exposure to chemicals that come with using traditional chemical disinfection methods.

Extra features assure protection

Each Aqua-Tec Water Preservation Cell comes with an optional suction cup for easy mounting on the reservoir wall if desired. Also included is a convenient sticker to remind you when to change to a new Aqua-Tec cell.



**Aqua-Tec Water Preservation Cell*
ordering information**

Description	Cat. No.
Aqua-Tec Water Preservation Cells value pack of 4 units	50109579

*Aqua-Tec Water Preservation Cell is intended for laboratory use only.

The Aqua-Tec Water Preservation Cell is proven effective against a variety of common microbial contaminants including bacteria, molds, and fungi.

thermo
scientific

Aqua-Tec
Water Preservation Cell

Year _____

JAN	FEB	MAR	APR	MAY	JUN	JUL	AUG	SEP	OCT	NOV	DEC
<input type="radio"/>	<input type="radio"/>	<input type="radio"/>	<input type="radio"/>	<input type="radio"/>	<input type="radio"/>	<input type="radio"/>	<input type="radio"/>	<input type="radio"/>	<input type="radio"/>	<input type="radio"/>	<input type="radio"/>

thermofisher.com/aquatec

Caution: Causes eye irritation and may be harmful if swallowed. Avoid contact with skin, eyes, or clothing, and wash thoroughly with soap and water after handling. This product is hazardous to fish.

EU BPR Registration CHZN2388

SilverSan AG, Rorschach Switzerland

Learn more at thermofisher.com/aquatec

thermo scientific

For Laboratory Use. It is the customer's responsibility to ensure that the performance of the product is suitable for customers' specific uses or applications. **This product is only available in selected countries.** Please consult your local sales representative for details. © 2019–2021 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. **COL016895 0821**