

Connection simplified

Technical features of the Centri-Vue App



The Thermo Scientific™ Centri-Vue™ Application provides real-time connection with select Thermo Scientific™ Superspeed and Large Capacity Centrifuges.

Check your run status or locate an open instrument on your mobile device – without leaving your office or lab space.

Key features


- Determine centrifuge status at-a-glance
- Establish secure connection for control
- Monitor your run status

The Centri-Vue application is available with the following Thermo Scientific centrifuges:

- Thermo Scientific™ Sorvall™ LYNX Superspeed Centrifuges
- Thermo Scientific™ Sorvall™ BP 8 and 16 Centrifuges
- Thermo Scientific™ Cryofuge™ 8 and 16 Centrifuges
- Thermo Scientific™ Sorvall™ BIOS A and 16 Centrifuges


Determine status of 1 to 100 centrifuges at-a-glance

Avoid hunting for an open instrument; instead find an idle centrifuge from the list

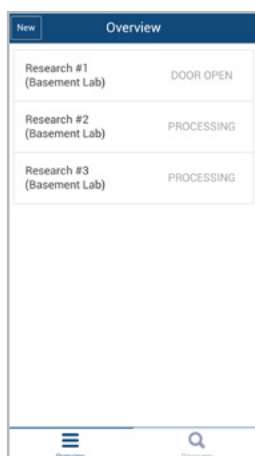


Establish secure centrifuge connection for start or stop control

Once you have set-up the run, start the instrument from the control panel OR your smart device

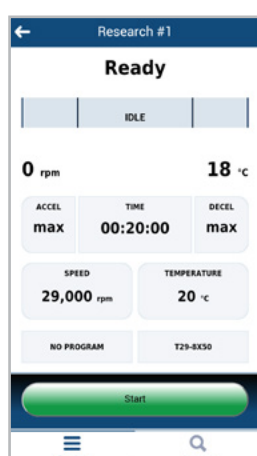


Replicate instrument main screen on your smart device


Monitor your run, with the option to stop the run from your smart device

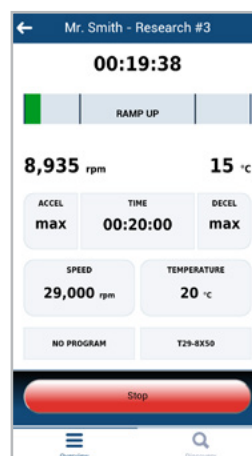

Check for diagnostic errors or if your run was stopped prior to completion

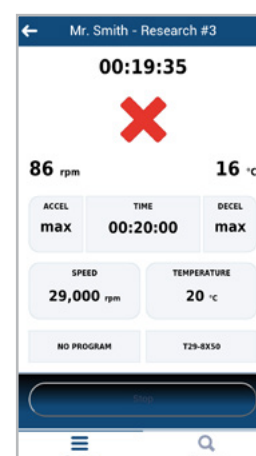

Application set-up instructions

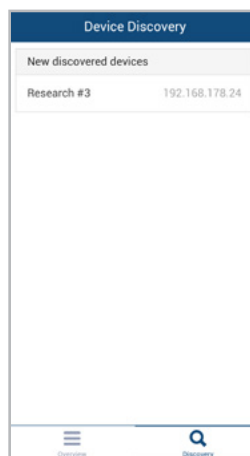

If connecting to a Sorvall LYNX centrifuge, make sure software version 604V7 (1.2.2) or above is installed. If not, schedule a service call to install the latest software update before proceeding.

Open the Configuration menu and select Device. On the Device screen, activate the DHCP check box.

Connect the centrifuge to a network using the onboard RJ-45 connector located behind the control panel or at the back of the centrifuge. Consult with your IT team to configure the IP address of the centrifuge on the network. Once configured, an IP address will be issued to the centrifuge by the network.

Refer to the Touchscreen User Interface manual for additional set-up instructions.

After installing the Centri-Vue app on a smart device, ensure the device is connected to the same network as the centrifuge. Then select Discover to locate you centrifuge and add it to your list of connected centrifuges. The IP address can also be added manually.

Optional: Edit the centrifuge information, providing a name for easy recognition.

Start using the Centri-Vue application to monitor and control your centrifuge!



Stay connected. Download the Centri-Vue application for instant remote monitoring and control of your Thermo Scientific centrifuges³, available for both iOS™ and Android™ devices.

1. Centri-Vue application requires iOS 7.0 or later or Android 5.1 or later.
2. User is responsible for network connection and set-up, as well as any smart device subscription charges.
3. Compatible with Sorvall LYNX centrifuges, Sorvall BIOS centrifuges, Sorvall BP 8 and 16 centrifuges as well as Cryofuge 8 and 16 centrifuges.

Learn more at thermofisher.com/floorcentrifuges

thermo scientific

For Laboratory Use. It is the customer's responsibility to ensure that the performance of the product is suitable for customers' specific uses or applications. © 2024 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. Apple and OS X are trademarks of Apple Inc., registered in the U.S. and other countries and region. Android is a trademark of Google LLC. Linux is the registered trademark of Linus Torvalds in the U.S. and other countries. **EXT6443 0224**