



Instructions for Using CryoMed 2.0 User Interface (UI) Emulator


- 1) Click on the “CryoMed 2.0 UI Emulator .zip folder”
 - a. The .zip folder will begin to download to your PC

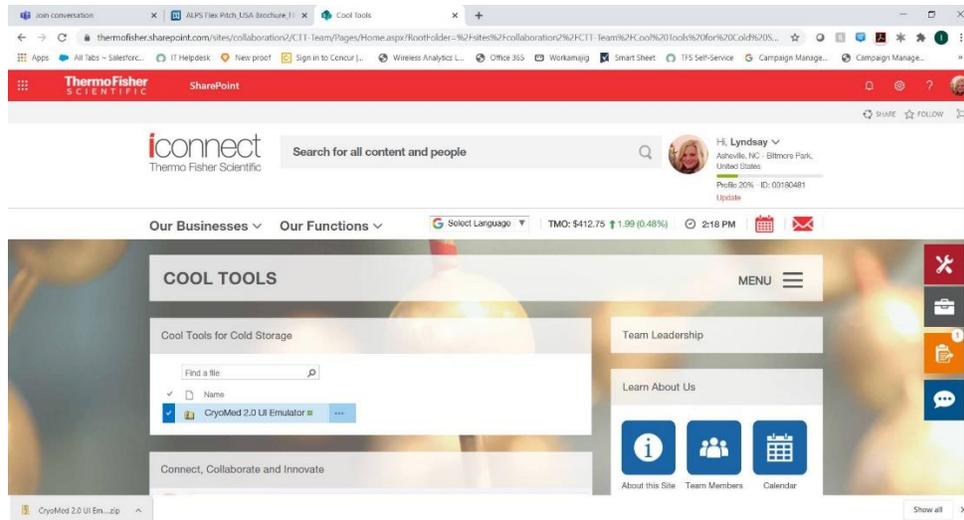

- 2) Once the folder is downloaded, it will appear at the bottom of your web browser, select the downloaded folder from the download bar.
- 3) After you clicked on the downloaded folder, the screen below will appear. Once this screen appears, select the “Extract all” options in the tool bar (or you can right-click on the .zip folder and select “Extract all”)

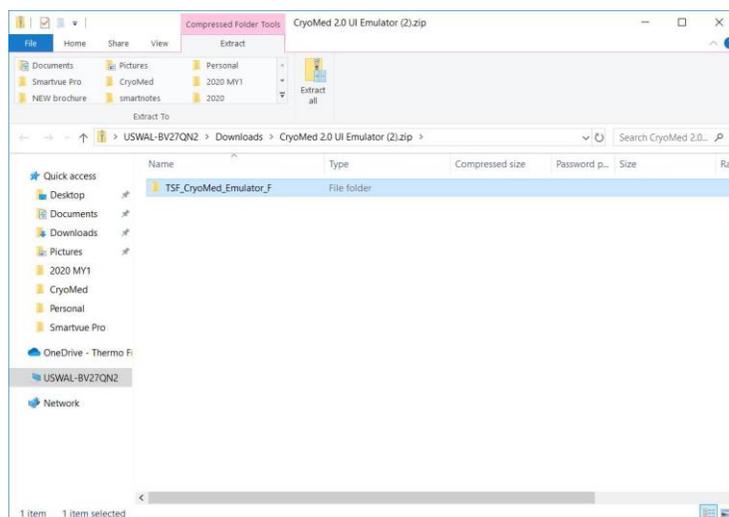

- 4) After selecting “Extract all,” the box below will appear on your screen. Select “Browse” and then save the contents to your “Desktop”

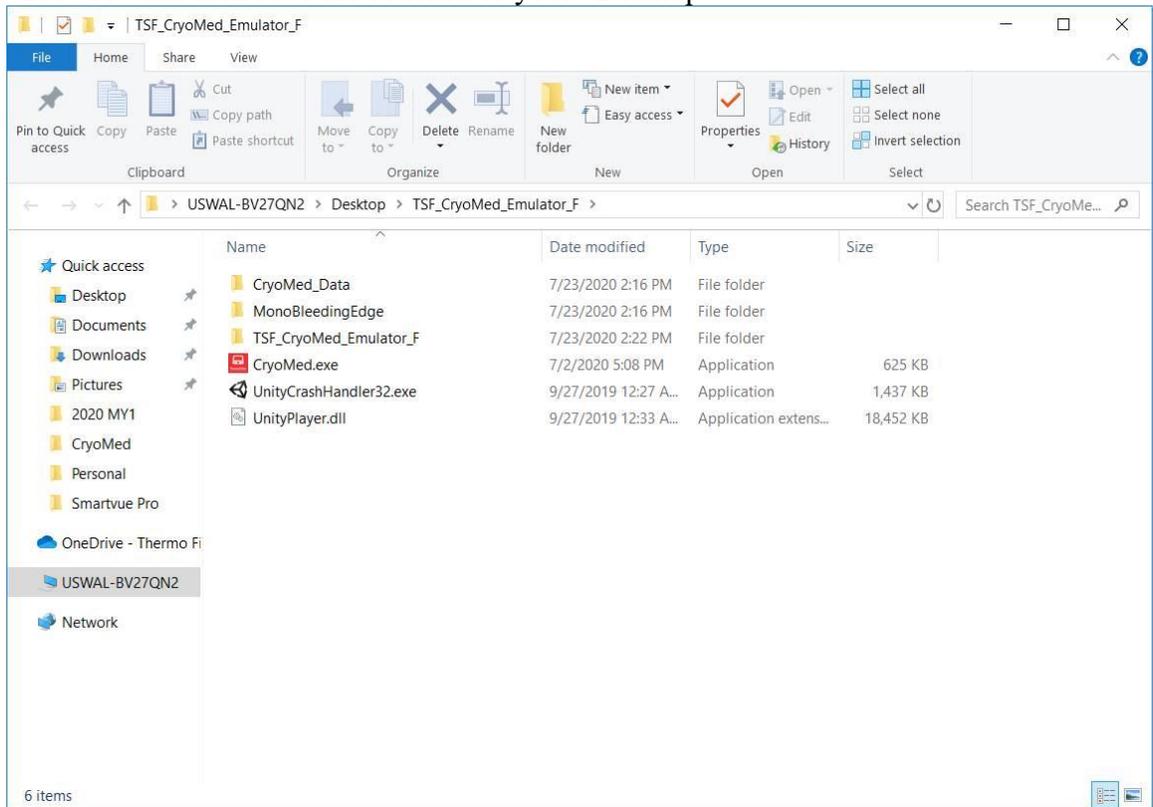


- 5) Once the folder is saved to your Desktop, select the  CryoMed icon.

- 6) After selecting the icon above, the following dialog box will appear. Select “Allow this file” in the dialog box



- 7) After you select “Allow this file” a second dialog box will appear. Select “OK” in this dialog box

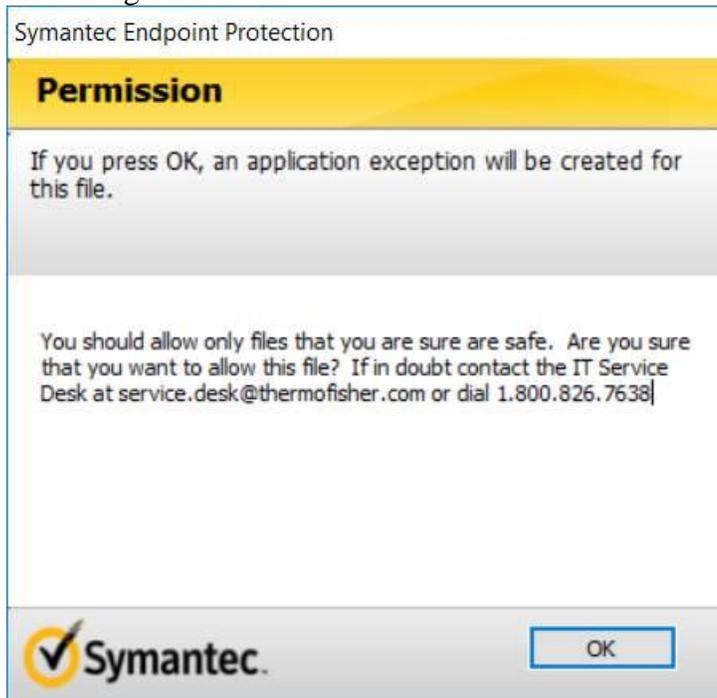

- 8) After selecting “OK” in the above dialog box, go back to your desktop folder where the Emulator files were stored and select the  CryoMed icon again.
- 9) After you select the CryoMed icon, the following screen appears. Click the “Start Setup” button. After you select the “Start Setup” button, the UI emulator will be active on your screen.

