

AccuSEQ[®] Software v2.0

Publication Part Number 4442195 Revision Date 25 January 2013 (Rev. B)

SUBJECT: Install AccuSEQ[®] Software v2.0

- Overview 1
- Computer requirements 2
- Before the installation 2
- Install AccuSEQ[®] Software v2.0..... 3
- Calibrate the instrument 4
- Uninstall AccuSEQ[®] Software v2.0 4

Overview

This user bulletin describes how to uninstall 7500 Software, SDS Software, or AccuSEQ[®] Software v1.0 and install AccuSEQ[®] Software v2.0 on an Applied Biosystems[®] 7500 Fast System.

The installation CD contains AccuSEQ[®] Software v2.0. The CD includes a registration code that you enter when you install the software.

IMPORTANT! You must use an Administrator account (unrestricted access) to install the software. After the software is installed, you can run AccuSEQ[®] Software v2.0 without using an Administrator account.

Computer requirements

Note: AccuSEQ® Software v2.0 has been tested using only Regional and Language options set to English (United States).

Component	Recommended Requirements
Computer	<ul style="list-style-type: none"> • 32-bit • Intel® CORE™ i7-2640M processor, 2.8 GHz • 4 GB RAM • Hard drive with 500 GB available • DVD-ROM drive • USB v2.0 • Ethernet network interface adapter (1000 BASE-T) • UL listed • CE marked • FCC labeled
Monitor	<ul style="list-style-type: none"> • 1280 × 1024 pixel resolution for full screen display • 17-inch • 32-bit color
Operating System	<ul style="list-style-type: none"> • Microsoft® Windows® 7, 32 bit, Service Pack 1 or later

Before the installation

Back up existing plates and templates

1. In Windows® Explorer, navigate to the location appropriate for your software:
 - **7500 Software:** x:\[directory]\7500\experiments
 - **SDS Software:** x:\[directory]\SDS Documents
 - **AccuSEQ® Software v1.0:** x:\[directory]\Applied Biosystems\AccuSEQ
 where *x* is the drive on which your software is installed.
2. Copy the data to a location outside the x:\Applied Biosystems\directory.

Uninstall the 7500 Software, SDS Software, or AccuSEQ® v1.0 Software

To uninstall the 7500 Software, SDS Software, or AccuSEQ® Software v1.0, use one of the following methods:

- Select **Start ▶ Settings ▶ Control Panel**, then select **Add or Remove Programs**. Select 7500 Software, SDS Software, or AccuSEQ® Software v1.0, then **Remove**, or
- Select **Start ▶ All Programs ▶ Applied Biosystems ▶ [Software] ▶ Uninstall [Software]** where *[Software]* is the name of the software that you want to uninstall.

Install AccuSEQ® Software v2.0

Log on requirements


To install AccuSEQ® Software v2.0, you must:

- Log on to the local computer (not a network domain).
- Use an Administrator account (unrestricted access).

Note: After the software is installed, you can run AccuSEQ® Software v2.0 without using an Administrator account.

Verify the installation user account

1. On the desktop, select **Start ▶ Control Panel**.
2. In the Control Panel window, click **User Account ▶ Manage User Accounts**.
3. In the Users tab, verify that the user account you will use to install the software belongs to the Administrators group and that the domain name is the same as the computer name.

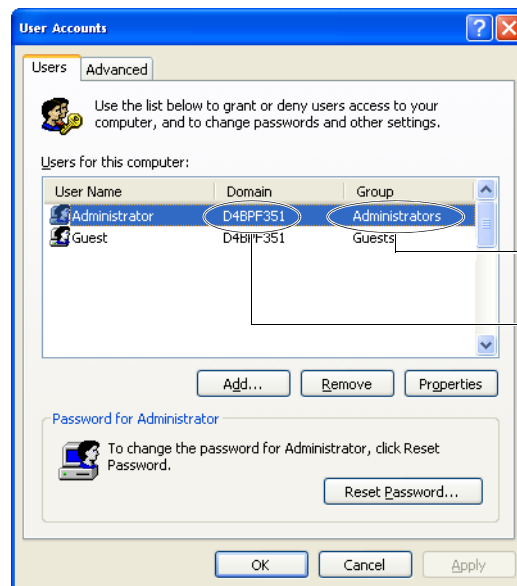

User account belongs to the Administrators group

Domain is the same as the computer name

To find the name of your computer:

- a. Click **Start ▶ Programs ▶ Accessories ▶ Command Prompt**.
- b. In the Command Prompt window, enter **hostname**.

Install the software

1. Insert the AccuSEQ® Software v2.0 CD in the CD-ROM or DVD-ROM drive, then follow the instructions to install.
2. When prompted, enter the registration code for the software (located on the CD packaging).



Calibrate the instrument

After you install the software, use the Instrument Maintenance Manager to perform ROI, background, optical, and pure dye calibrations. Then perform an RNase P verification run. For information, see the *Applied Biosystems® 7500/7500 Fast Real-Time PCR Systems Maintenance Guide* (Pub. no. 4387777).

Uninstall AccuSEQ® Software v2.0

If you need to uninstall AccuSEQ® Software v2.0, note the following:

- You must log on to the local computer with an Administrator account. See [page 3](#) for information.
- During the uninstallation, the software:
 - Creates a backup file containing the x:\Applied Biosystems\AccuSEQ\experiments folder
 - Names the file Backup_M-DD-YYYY_software version #_uninstall instance #, for example: Backup_6-15-2012_2.0
 - Stores the backup file x:\Applied Biosystems\AccuSEQ\Backup. This directory is not removed during the uninstallation.

For Research Use Only. Not for use in diagnostic procedures.

DISCLAIMER: LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF

NOTICE TO PURCHASER: PLEASE REFER TO THE ACCUSEQ SOFTWARE V2.0 MYCOPLASMA SEQ EXPERIMENTS GETTING STARTED GUIDE FOR LIMITED LABEL LICENSE OR DISCLAIMER INFORMATION.

© 2013 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation and/or its affiliate(s) or their respective owners. Intel is a registered trademark and Core is a trademark of Intel Corporation. Microsoft, Windows, and Internet Explorer are registered trademarks of Microsoft Corporation.

Headquarters

5791 Van Allen Way | Carlsbad, CA 92008 USA
Phone +1 760 603 7200 | Toll Free in USA 800 955 6288

For support visit www.lifetechnologies.com/support

www.lifetechnologies.com

