

MicroSEQ® *Listeria monocytogenes* Detection Kit

Fast, Accurate Food Pathogen Testing

- Easy-to-use, streamlined lyophilized format
- Fast time-to-results
- High sensitivity and specificity for confidence in results
- Optimal performance with wide range of food sample types
- Internal positive control helps eliminate false negatives



A Dangerous Foodborne Pathogen

Tasteless, odorless, and impossible to see with the naked eye, *Listeria monocytogenes* is responsible for listeriosis, a disease with a high fatality rate in susceptible populations, including newborns, immunocompromised individuals, and the elderly, causing the highest mortality rate among foodborne illnesses. Listeriosis is particularly dangerous for pregnant women, potentially causing miscarriage or stillbirth.

A Concern for the Food Industry

In addition to its impact on health, an outbreak of listeriosis can be devastating to food processors. If the outbreak is traced to a facility, the operator may be forced to recall and destroy contaminated products, shut down operations, and face lost revenue, damaged reputation, fines, and litigation. Containment is complicated by the fact that *Listeria* can grow at temperatures as low as 3°C, allowing it to multiply in refrigerated foods during storage, shipping, and retail display.

The Testing Challenge

Listeria monocytogenes is one of six species of the genus *Listeria* and the only identified species to cause listeriosis. *Listeria monocytogenes* bacteria grow slowly and are difficult to distinguish from other species of *Listeria* on agar culture media. This has prompted many food-testing laboratories to search for a test method that provides rapid, accurate results—with high sensitivity and specificity—that is easy to use and implement. The MicroSEQ® *Listeria monocytogenes* Detection Kit meets these criteria.

Benefits of Real-Time PCR

Real-time PCR is a proven method for pathogen detection and has been applied successfully to a wide range of foodborne pathogens, including *Salmonella*, *Listeria monocytogenes*, and *E. coli* O157:H7. Real-time PCR tests for food pathogens by amplifying and detecting a DNA target sequence that is specific to the organism under investigation. In this way, real-time

PCR can detect pathogens quickly, simply, and with outstanding sensitivity and specificity. And since the kit uses fast PCR chemistry, the PCR step takes only 40 minutes, compared with up to 2.5 hours using standard PCR.

Improved Time-to-Results

Real-time PCR offers faster time-to-results than other methods. *Listeria* testing using traditional culture-based methods can take up to 5 days. Immunoassay-based methods take up to 2 days. Time-to-results with the MicroSEQ® *Listeria monocytogenes* Detection Kit is less than 3 hours after preenrichment, and less than 27 hours total.

High Specificity and Sensitivity

Because the MicroSEQ® *Listeria monocytogenes* Detection Kit detects genetic material unique to the organism, it provides both high specificity and sensitivity. With culture-based methods, interpretation of the results is highly subjective, which may lead to missed or ambiguous identification of *Listeria monocytogenes*. Immunoassay methods, while less subjective than culture methods, are based on antibody-binding mechanisms that may be prone to interference and lead to false positives and/or false negatives.

The MicroSEQ® *Listeria monocytogenes* Detection Kit specifically detects the following *Listeria* serotypes: 1/2A, 1/2B, 1/2C, 3A, 3B, 3C, 4A, 4AB, 4B, 4C, 4D, 4E, and 7. The kit does not detect other pathogens.

In combination with the PrepSEQ™ sample preparation protocols, the kit can detect 1–3 colony forming units (cfu) per 25 grams of sample, with a limit of detection of 10⁴ cfu/mL.

Demonstrated Performance

Performance of the MicroSEQ® *Listeria monocytogenes* Detection Kit has been demonstrated on a variety of foods, including:

- Meat products
- Seafood products
- Fruits and juices
- Vegetable products
- Dairy products
- Infant formula
- Salad dressings



Figure 1. Reagents for the MicroSEQ® *Listeria monocytogenes* Detection Kit are provided in a ready-to-use lyophilized format for convenience and reproducibility.

Lyophilized for Efficiency and Ease of Use

For maximum ease of use, reliability, and consistency of results, the reagents used in the MicroSEQ® *Listeria monocytogenes* Detection Kit are lyophilized into preformatted assay beads. The beads hold the active enzyme, the target-specific primer and probe set, internal positive control (IPC), and other reagents for PCR. The IPC is provided to help eliminate false negatives by detecting the presence of materials that can inhibit target amplification.

Closed-Tube Integrity

The MicroSEQ® *Listeria monocytogenes* Detection Kit uses specially designed reaction tubes that remain closed throughout the assay process. Once the sample is added, the tubes are closed and remain that way until detection is complete, greatly reducing the chances of contamination.

No electrophoresis or post-PCR processing is required. All the operator has to do is to prepare the assay beads, add the samples and controls, and run the test. Sample handling is minimal, and every step is guided by the RapidFinder™ Express Software with on-screen instructions.

Optimized Sample Preparation

The MicroSEQ® *Listeria monocytogenes* Detection Kit is optimized for use with the PrepSEQ™ Sample Preparation Kits. The PrepSEQ™ kits remove PCR inhibitors, enhance assay performance, and enable the user to enrich and prepare high-quality



Figure 2. Closed-tube detection reduces chances of contamination.

samples from a broad range of food sample types for detection of 1–3 cfu per 25 grams of sample.

- For automated high-throughput applications, the PrepSEQ™ Nucleic Acid Extraction Kit prepares high-quality microbial DNA and RNA from broth cultures when used with the Applied Biosystems MagMAX™ Express-96 Sample Preparation System.
- For lower sample numbers, the PrepSEQ™ Rapid Spin Sample Preparation Kit provides a simple, cost-effective way to prepare DNA from broth cultures.

Fast, Actionable Answers

When the assay is complete, RapidFinder™ Express Software presents an easy-to-read screen that allows the user to view the results in each reaction location. Results are clearly displayed and can be labeled with flags, notifications, and prompts that enable

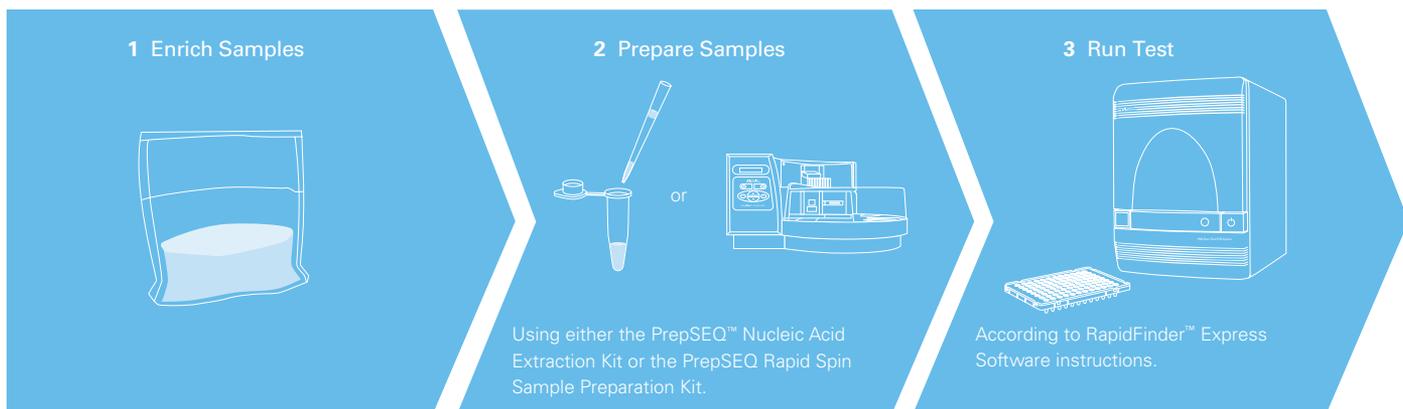


Figure 3. The MicroSEQ® *Listeria monocytogenes* Detection Kit streamlines the assay workflow with optimized sample preparation and assay procedures. Nontechnical personnel with minimal training or prior experience can perform the assay in 3 simple steps.

the operator to quickly interpret the data and take appropriate action.

A Complete Solution

The MicroSEQ® *Listeria monocytogenes* Detection Kit is part of a complete food-testing solution. The kit includes everything required to run 96 reactions. All components have been designed for rapid implementation in food-testing laboratories and performance-verified to make detection as fast, easy, and reliable as possible. Everything is provided ready to use.

- **Optimized for Sensitivity and Specificity:** The assay is designed to provide maximum sensitivity on the 7500 Fast Real-Time PCR System.
- **Ready to Use:** The active enzyme, reagents, primers and probes, and internal positive control are lyophilized into preformatted assay beads. No mixing is required.
- **Optimized Sample Preparation:** A choice of sample preparation kits helps ensure high-quality assay results.
- **Software-Guided:** Application-specific

RapidFinder™ Express Software guides the user through each step of the procedure—from run file setup to final results.

- **AOAC-Certified:** The full workflow including the MicroSEQ® *Listeria monocytogenes* Detection Kit, PrepSEQ™ Sample Preparation Kits, 7500 Fast instrument, and RapidFinder™ Express Software has earned the *Performance Tested Methods*™ certification from the AOAC Research Institute.

food pathogens. We also provide responsive, knowledgeable applications consulting, support, training, and technical service.

For more information about the MicroSEQ® *Listeria monocytogenes* Detection Kit and our other solutions for food pathogen testing, please contact your local Applied Biosystems sales representative or visit us at www.appliedbiosystems.com.

Resources for Food Safety

As the world leader in real-time PCR, Applied Biosystems is committed to providing the food industry with improved tools for pathogen detection. The MicroSEQ® *Listeria monocytogenes* Detection Kit is part of a growing family of fast and convenient food pathogen detection tools that utilize lyophilized reagents, application-specific software, optimized sample preparation, and fast real-time PCR instrumentation. Other solutions in our expanding portfolio include easy-to-use TaqMan® Pathogen Detection Kits for rapid detection of a broad menu of

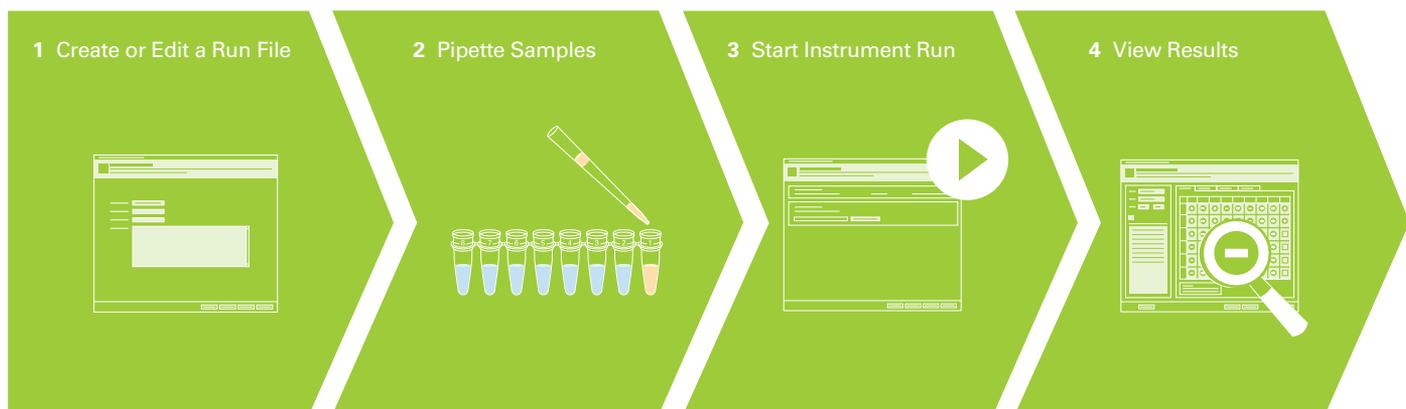


Figure 4. Software-guided procedure. After sample enrichment and preparation, RapidFinder™ Express Software guides the user through the entire workflow with on-screen instructions. Amplification, detection, data collection, and analysis are fully automated.

ORDERING INFORMATION

Description	Part Number
MicroSEQ® Pathogen Detection Kits	
MicroSEQ® <i>Listeria monocytogenes</i> Detection Kit	4403874
MicroSEQ® <i>Salmonella</i> spp. Detection Kit	4403930
TaqMan® Pathogen Detection Kits	
TaqMan® <i>Campylobacter jejuni</i> Detection Kit	4366098
TaqMan® <i>Cronobacter sakazakii</i> Detection Kit	4382492
TaqMan® <i>E. coli</i> O157:H7 Detection Kit	4366100
TaqMan® <i>Listeria monocytogenes</i> Detection Kit	4366102
TaqMan® <i>Pseudomonas aeruginosa</i> Detection Kit	4368604
TaqMan® <i>Salmonella enterica</i> Detection Kit	4366104
TaqMan® <i>Staphylococcus aureus</i> Detection Kit	4368606
Sample Preparation	
PrepSEQ™ Nucleic Acid Extraction Kit	4400799
PrepSEQ™ Rapid Spin Sample Preparation Kit with Proteinase K	4426714
PrepSEQ™ Rapid Spin Sample Preparation Kit – Extra Clean with Proteinase K	4426715
Instrumentation and Software	
Food Pathogen Detection System Package: Applied Biosystems 7500 Fast Real-Time PCR System with PC Tower and RapidFinder™ Express Software	4445785
Food Pathogen Detection System Package: Applied Biosystems 7500 Fast Real-Time PCR System with Notebook Computer and RapidFinder™ Express Software	4445787

For Research Use Only. Not intended for any animal or human therapeutic or diagnostic use.

For environmental (including soil, air and water), quality control, and food and finished products testing; for research use only. Not for use in diagnostic procedures (including, in the United States, epidemiology use with human or animal samples by public health officials). This product has not been cleared or otherwise approved by the United States Food and Drug Administration or by any other regulatory body in any country, or under the European IVD Directive, for diagnostic or any other clinical purpose. The purchaser or user of this product agrees not to use this product for diagnostic or other clinical purposes.

NOTICE TO PURCHASER: Limited Use Label License

The products shown in this bulletin may be covered by one or more Limited Use Label License(s). Please refer to the respective product documentation or the Applied Biosystems website under www.appliedbiosystems.com for the comprehensive license information. By use of these products, the purchaser accepts the terms and conditions of all applicable Limited Use Label Licenses. These products are sold for research use only, and are not intended for human or animal diagnostic or therapeutic uses unless otherwise specifically indicated in the applicable product documentation or the respective Limited Use Label License(s).

© 2010 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation or their respective owners. TaqMan is a registered trademark of Roche Molecular Systems, Inc.

Printed in the USA. 03/2010 Publication 116PB12-02 C011354 0410