

miRNA QC tool version 1.1.1 Release Notes

miRNA QC tool version 1.1.1

This software is also known as “AffyMir” version 1.1.1

Updates in version 1.1.1

1. An Array type check was added to verify the array type between the CEL and CDF files was the same.
2. The installer was updated to install both version 11 and version 15 library files.

Installation

1. Installs AffyMir version 1.1.1 executables
2. Installs library files for both v11 and v15 miRNA arrays
3. Compatible with Windows XP, Windows 7
4. Runs only on 32 bit systems.

Known Issues

Issue: Software exits when running analysis with mismatched library files.

Description:

When loading v2 CEL files with a v1 library files set (or vice versa), the user will receive an error message indicating that the CEL files and library files do not match.

Once the dialog is dismissed, the ‘Analysis’ menu option becomes enabled, as well as the ‘Run Analysis’ option. If you select the analysis with the incorrect library files, the application will exit and you will have to re-start it.

Ensure you have selected the correct library files before proceeding to run analysis.

Support

If you encounter issues which you are unable to resolve on your own, please contact your Affymetrix Support Center.

Affymetrix, Inc. (U.S.)
Tel. 888-DNA-CHIP (888-362-2447)
Fax. 408-731-5441
support@affymetrix.com
sales@affymetrix.com

Affymetrix UK Ltd.
UK and Others Tel. +44 (0) 1628 552550
France Tel. +33 800919505
Germany Tel. +49 1803001334
Fax. +44 (0) 1628 552598
supporteurope@affymetrix.com
saleseurope@affymetrix.com