

# Mouse IFN- $\gamma$ Antibody Pair

Quantity: 10 Plate Format

Catalog Number CMC4033

Lot #: 1721366

Pub No. MAN0014791 Rev. 1.0


**\*Note:** A letter at the end of the lot number signifies an additional packaging of this same lot.

## Product use and materials provided

The Antibody Pair for Mouse IFN- $\gamma$  contains components required to construct an enzyme-linked immunoassay for the specific and quantitative measurement of IFN- $\gamma$ . Sufficient quantities of all reagents are provided to yield 10 plates of 96 wells if the recommended assay procedure and recommended storage and handling of materials are followed as specified on this insert.

<b>Coating Antibody:</b>	<b>Mouse IFN-<math>\gamma</math> Coating Antibody (0.0625 mg/0.125 mL)</b>
Part Number:	AMC4739D
Lot Number:	1721367
Form:	Liquid, 4 vials, contains 0.1% sodium azide.
Storage:	Store at 2-8°C for 1 month. For longer periods, aliquot and store at $\leq -20^\circ\text{C}$ .
Recommended Dilution:	Dilute to 2.0 $\mu\text{g/mL}$ with Coating Buffer B (Cat. # CB01100, or see Recommended Buffers). For example, to make 10 mL (enough to coat 1 plate), add 40 $\mu\text{L}$ coating antibody to 9.960 mL of Coating Buffer B.
<b>Detection Antibody:</b>	<b>Mouse IFN-<math>\gamma</math> Detection Antibody (0.025 mg/0.125 mL)</b>
Part Number:	AMC4934D
Lot Number:	1721368
Form:	Liquid, 1 vial, contains 0.1% sodium azide.
Storage:	Store at 2-8°C for 1 month. For longer periods, aliquot and store at $\leq -20^\circ\text{C}$ .
Recommended Dilution:	Dilute to 0.06 $\mu\text{g/mL}$ with Assay Buffer <b>supplemented with 10% calf serum</b> (Cat. # DS98200, or see Recommended Buffers). For example, to make enough for 1 plate, add 1.8 $\mu\text{L}$ detection antibody to 5.998 mL Assay Buffer.
<b>Standard:</b>	<b>Mouse IFN-<math>\gamma</math> Standard</b>
Part Number:	SD0242 (inquire regarding additional vials)
Lot Number:	1736106
Form:	Lyophilized, 3 vials (single use).
Storage:	Store at 2-8°C.
Reconstituted Standard Concentration:	3,000 $\text{pg/mL}$ .
Reconstitution:	Reconstitute in Assay Buffer (Cat. # DS98200 or see Recommended Buffers) according to instructions on vial label. Allow standard to rehydrate for approximately 10 minutes before dilutions. If the standard stock is not being used immediately, please aliquot into polypropylene tubes and freeze at $-80^\circ\text{C}$ . <i>Do not store at room temperature or at <math>4^\circ\text{C}</math> for any extended time or subject to more than one freeze-thaw cycle.</i>
Recommended Starting Standard Curve:	Dilute standard stock to 1000 $\text{pg/mL}$ (a 1:3 dilution) followed by six 1:2 serial dilutions using at least 300 $\mu\text{L}$ of buffer. Mix thoroughly between dilutions. Avoid foaming. To an empty tube add 300 $\mu\text{L}$ of buffer and label as zero standard.
<b>Streptavidin-HRP:</b>	<b>0.1 mg/0.5 mL</b>
Part Number:	SNN4004W
Lot Number:	1721369
Form:	Liquid, 1 vial, contains animal serum and 50% glycerol in phosphate buffered saline with 0.05% thymol as a preservative.
Storage:	Store concentrate at 2-8°C for 1 month. For longer periods, aliquot and store at $\leq -20^\circ\text{C}$ . Diluted streptavidin-HRP should not be stored; discard remaining solution after use.
Recommended Dilution:	Dilute to 0.1 $\mu\text{g/mL}$ . For example, to make enough for 1 plate, add 5.0 $\mu\text{L}$ of streptavidin-HRP to 9.995 mL of Assay Buffer (Cat. # DS98200 or see Recommended Buffers).

Following the recommended assay procedure using the **Antibody Pair Buffer Kit (Cat. # CNB0011)**, the representative standard curve was generated.

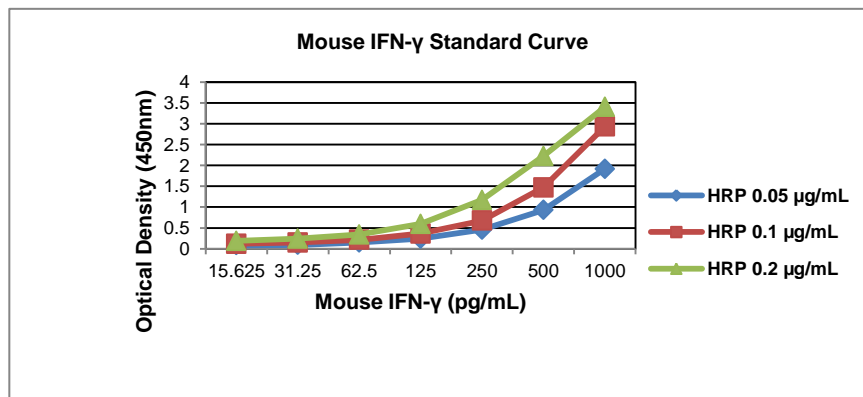

## Additional materials required

- 96 well NUNC™ MaxiSorp microplates; NUNC Cat. # 434797.
- Pipettes, plate covers or plate sealers, and timer.
- Microplate reader with a detector that can measure absorbance at 450 nm.
- 1 L graduated cylinder; plate washer or wash bottle.
- Polypropylene tubes for standards and sample dilutions, if needed.

## Recommended buffers and solutions

The Antibody Pair Buffer Set (Cat. # CNB0011) containing Coating Buffers A and B, Assay Buffer, Substrate Solution (TMB), Stop Solution, and Wash Buffer is recommended.

1. **Coating Buffer A:** *Coating Buffer A (Cat. # CB07100) from Invitrogen is recommended. Alternate buffer choice listed below.*  
8.0 g NaCl, 1.13 g Na<sub>2</sub>HPO<sub>4</sub>, 0.2 g KH<sub>2</sub>PO<sub>4</sub>, 0.2 g KCl, 0.1% ProClin™; q.s. to 1.0 L with distilled H<sub>2</sub>O, pH to 7.4.
2. **Coating Buffer B:** *Coating Buffer B (Cat. # CB01100) from Invitrogen is recommended. Alternate buffer choice listed below.*  
4.3 g NaHCO<sub>3</sub>, 5.3 g Na<sub>2</sub>CO<sub>3</sub>, 0.1% ProClin™; q.s. to 1.0 L with distilled H<sub>2</sub>O, pH to 9.4.
3. **Assay Buffer:** *Assay Buffer (Cat. # DS98200) from Invitrogen is recommended. Alternate buffer choice listed below.*  
8.0 g NaCl, 1.13 g Na<sub>2</sub>HPO<sub>4</sub>, 0.2 g KH<sub>2</sub>PO<sub>4</sub>, 0.2 g KCl, 5.0 g bovine serum albumin (fraction V), 1 mL Tween-20™ Detergent and 0.5% ProClin™ as a preservative; q.s. to 1.0 L with distilled H<sub>2</sub>O, pH to 7.4.
4. **Wash Buffer:** *Wash Buffer 25x (Cat. # WB01) from Invitrogen is recommended. Alternate buffer choice listed below.*  
0.2 g KH<sub>2</sub>PO<sub>4</sub>, 1.9 g K<sub>2</sub>HPO<sub>4</sub>·3H<sub>2</sub>O, 0.4 g EDTA, 0.5 mL Tween-20™ Detergent; q.s. to 1.0 L with distilled H<sub>2</sub>O, pH to 7.4.
5. **Substrate Solution:** *TMB (Cat. # SB01) from Invitrogen is recommended. Alternate solution choice listed below.*  
Tetramethylbenzidine (TMB) and Hydrogen Peroxide.
6. **Stop Solution:** *Stop Solution (Cat. # SS03100) from Invitrogen is recommended. Alternate solution choice listed below.* 0.4 N H<sub>2</sub>SO<sub>4</sub>.

## Assay optimization

Antibody Pair kits from Invitrogen™ are designed to be very flexible for your experiments. Consequently, the assay procedure contains only recommendations. The assay procedure has been optimized for use with tissue culture samples. However, serum and plasma samples may be used but may require that certain assay parameters be modified. Investigators are advised to determine optimal buffer formulations, concentrations and incubation times for individual applications.








## Recommended assay procedure

1. Prepare coating solution by diluting the coating antibody. See “Coating Antibody” section for the recommended coating antibody dilution.
2. Coat plates with 100 µL per well of the coating solution. Cover plates and incubate overnight (12-18 hr.) at 4°C.
3. Aspirate wells and wash 1 time with > 400 µL of Wash Buffer per well. Following wash, invert and tap on absorbent paper to remove excess liquid.
4. Block plate with 300 µL per well of Assay Buffer for 1 hour at room temperature.
5. Aspirate, invert, and tap on absorbent paper to remove excess liquid.
6. Prepare standards and sample dilutions in Assay Buffer (or in a diluent that most closely matches the matrix of your sample).
7. Pipette 100 µL of standards (in duplicate), samples and controls into designated wells.
8. Immediately following step 7, add 50 µL of the working detection antibody into each well. For recommended dilutions, see “Detection Antibody” section. Gently tap the plate on the side 10 times to mix. *Cover plate and incubate for 2 hours at room temperature.*
9. Aspirate and wash 5 times using the method in step 3.
10. Add 100 µL of the working streptavidin-HRP solution into each well. For recommended dilutions, see the “Streptavidin-HRP” section. *Cover plate and incubate for 30 minutes at room temperature.*
11. Aspirate and wash 5 times using the method in step 3.
12. Add 100 µL of the TMB substrate to each well. *Incubate plate without a plate cover for 30 minutes in the dark at room temperature.*
13. Add 100 µL of Stop Solution to each well.
14. Measure absorbance at 450 nm (reference absorbance: 650 nm) within 30 minutes of adding Stop Solution. Calculate results using a log-log or 4-parameter curve fit.

## Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at [www.lifetechnologies.com/termsandconditions](http://www.lifetechnologies.com/termsandconditions). If you have any questions, please contact Life Technologies at [www.lifetechnologies.com/support](http://www.lifetechnologies.com/support).

## Product label explanation of symbols and warnings

 REF	Catalog Number	 LOT	Batch code	 Temperature limitation	 Use by	 Manufacturer	 Consult instructions for use	 Caution, consult accompanying documents
--	----------------	---	------------	--	--	--	--	---

**DISCLAIMER:** LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

Corporate entity: Life Technologies | Carlsbad, CA 92008 USA | Toll Free in USA 1.800.955.6288

© 2015 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

**For research use only. Not for use in diagnostic procedures.**

Manufacturing Site \* 7335 Executive Way \* Frederick \* MD 21704 \* Tel: 800.955.6288 \* E-mail: [techsupport@lifetech.com](mailto:techsupport@lifetech.com)

[thermofisher.com](http://thermofisher.com)

MAN0014791 Rev: 1.0 Effective: 13 Jan 2016 (Page 2 of 2)