


SureTect *Listeria* species PCR Assay Workflow

NF VALIDATION ISO 16140: Relative Level of Detection

Ana-Maria Leonte and Jonathan Cloke, Thermo Fisher Scientific, Wade Road, Basingstoke, Hampshire, RG24 8PW, UK

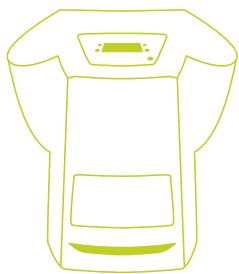

SureTect™ PikoReal™
Real-Time PCR
Instrument

Summary

As part of the NF VALIDATION™ ISO 16140 validation of the Thermo Scientific™ SureTect™ *Listeria* species PCR Assay workflow (alternative method), a relative level of detection (RLOD) study was conducted by ADRIA Développement, Quimper, France. The following is a summary of that study.

Methodology

Choice of strains: Five *Listeria* spp isolates were selected from the culture collection at ADRIA Développement and were spiked into all the matrices analyzed during the NF VALIDATION ISO 16140 relative level of detection study.

Protocol: For each matrix/strain combination, four levels of contamination were prepared and for each level of contamination, six replicates were analyzed. The target contamination levels were: 0 CFU/g or ml, the contamination level required to obtain 0% to 50% positive samples, the contamination level required to obtain 50% to 75% positive samples and the contamination level required to get 75% to 100% positive samples. The samples were analyzed using the reference method detailed in ISO 11290-1:1996, including Amendment 1:2004 prior to inoculation in order to verify the absence of *Listeria* spp. After inoculation, samples were tested using the ISO reference method and the alternative method.

Alternative method: Twenty-five gram samples were homogenized with 225 ml of supplemented Thermo Scientific™ Oxoid™ 24 *Listeria* Enrichment Broth (24 LEB). The sample was enriched by incubating for 22 to 26 hours at 37 ± 1 °C. Ten microlitres of SureTect Proteinase K Reagent and 10 µl of SureTect Lysis Reagent 2 were added to each of the required SureTect Lysis Tubes (supplied prefilled with Lysis Reagent 1). Ten microlitres of each of the required number of enriched samples were added to the Lysis Tubes, which were then heated at 37 ± 1 °C for 10 minutes, followed by 95 ± 1 °C for 5 minutes. The tubes were cooled for at least 2 minutes at room temperature prior to transferring 20 µl aliquots of the lysates to SureTect *Listeria* species PCR Tubes containing SureTect *Listeria* species PCR tablets. The PCR Tubes were then immediately sealed and transferred to the Thermo Scientific™ SureTect™ PikoReal™ Real-Time PCR Instrument for processing.

ISO method: Twenty-five gram samples were analyzed according to the reference method detailed in ISO 11290-1:1996, including Amendment 1:2004. Each sample was enriched by incubating at 30±1 °C for 22 to 26 hours in 225 ml of Half Fraser Broth. After primary enrichment, 10 µl of the Half Fraser Broth enrichment was inoculated onto both Thermo Scientific™ Oxoid™ Chromogenic Listeria Agar (OCLA) (ISO) and PALCAM Agar and incubated at 37±1 °C, for 22 to 26 hours. A further 100 µl of the Half Fraser Broth enrichment was inoculated into 10 ml Fraser Broth and incubated for 46 to 50 hours at 37±1 °C, from which OCLA (ISO) and PALCAM Agar plates were streaked and incubated for 22 to 26 hours at 37±1 °C. Presumptive positive growth from up to five colonies was confirmed by Gram stain, haemolysis, catalase, CAMP test and biochemical identification.

Results: The levels of detection for the alternative method and ISO reference method were determined according to the Spearman-Kärber (LOD₅₀) method (Table 1). The aim was to obtain the relative levels of detection for each of the “strain-matrix” combinations analyzed during the AFNOR Certification validation study.

Table 1: Relative detection level results for the alternative method and ISO reference method according to the Spearman-Kärber test

Strain / matrix pairs	ISO reference method (CFU/25 g)	Alternative method (CFU/25 g)
Rillettes / <i>L. monocytogenes</i> Ad 669	0.658 [0.339-1.277]	0.405 [0.229-0.714]
Smoked salmon / <i>L. innocua</i> 1	0.745 [0.492-1.154]	0.587 [0.366-0.942]
Raw milk / <i>L. ivanovii</i> Ad 991	0.705 [0.381-1.305]	0.792 [0.408-1.536]
Raw vegetables / <i>L. seeligeri</i> Ad 1293	0.603 [0.339-1.071]	0.603 [0.356-1.021]
Process water / <i>L. welshimeri</i> Ad 1252	0.828 [0.541-1.268]	0.828 [0.541-1.268]

Conclusions

The relative level of detection study conducted as part of the NF VALIDATION study demonstrated that the SureTect Listeria species PCR Assay workflow showed similar detection levels for the matrices analyzed compared to the ISO reference method, detailed in ISO 11290-1:1996, including Amendment 1:2004. For the ISO reference method, the level of detection range for all food categories studied during the validation was 0.3-1.3 CFU/25 g, compared to a range of 0.2-0.7 CFU/25 g for the alternative method, demonstrating that the SureTect Listeria species PCR assay workflow is an accurate alternative method for the detection of *Listeria* spp for the matrices tested as part of the NF VALIDATION study. The NF VALIDATION certificate and the validation report summary for this study are available from <http://nf-validation.afnor.org/en/>.

www.thermofisher.com/SureTect

© 2016 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. ATCC is a trademark of the American Type Culture Collection. NF VALIDATION is a trademark of Association Française de Normalisation (AFNOR). This information is presented as an example of the capabilities of Thermo Fisher Scientific Inc. products. It is not intended to encourage use of these products in any manners that might infringe the intellectual property rights of others.

Contact Information:

microbiology@thermofisher.com
USA +1 800 255 6730
International +44 (0) 1256 841144

993-220
LT2238A
March 2016

Thermo
SCIENTIFIC

A Thermo Fisher Scientific Brand