

Software Installation Guide

Thermo Scientific™ Oxoid™ Microbact™ Computer Aided Identification System

System Requirements

The Microbact Computer Aided Identification System should be compatible with any standard PC capable of running Windows® 7, 10, or 11.

Microsoft® and Windows® are registered trademarks of Microsoft Corporation.

**CAUTION:**

Microbact Computer Aided Identification System may not install or function correctly if this guide is not followed.

**CAUTION:**


If you have an existing earlier version of Microbact Computer Aided Identification System, close it before proceeding with the upgrade.

Uninstalling Previous Versions


If re-installing the application on an existing system, the previous version will be uninstalled automatically.

Installing Microbact Computer Aided Identification System Desktop Software


1. If the Microbact Computer Aided Identification System desktop software is open, close it.
2. Download the zipped software package MB1244A.zip.
3. Extract all files from MB1244A.zip to a temporary location on the PC's hard drive.

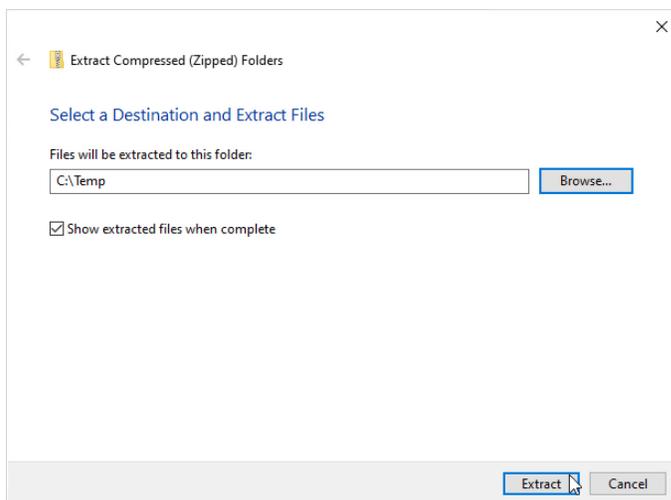


4. With Windows Explorer navigate to the folder from step 3, right-click the file “setup.exe”, and select “Run as administrator” from the pop-up menu.

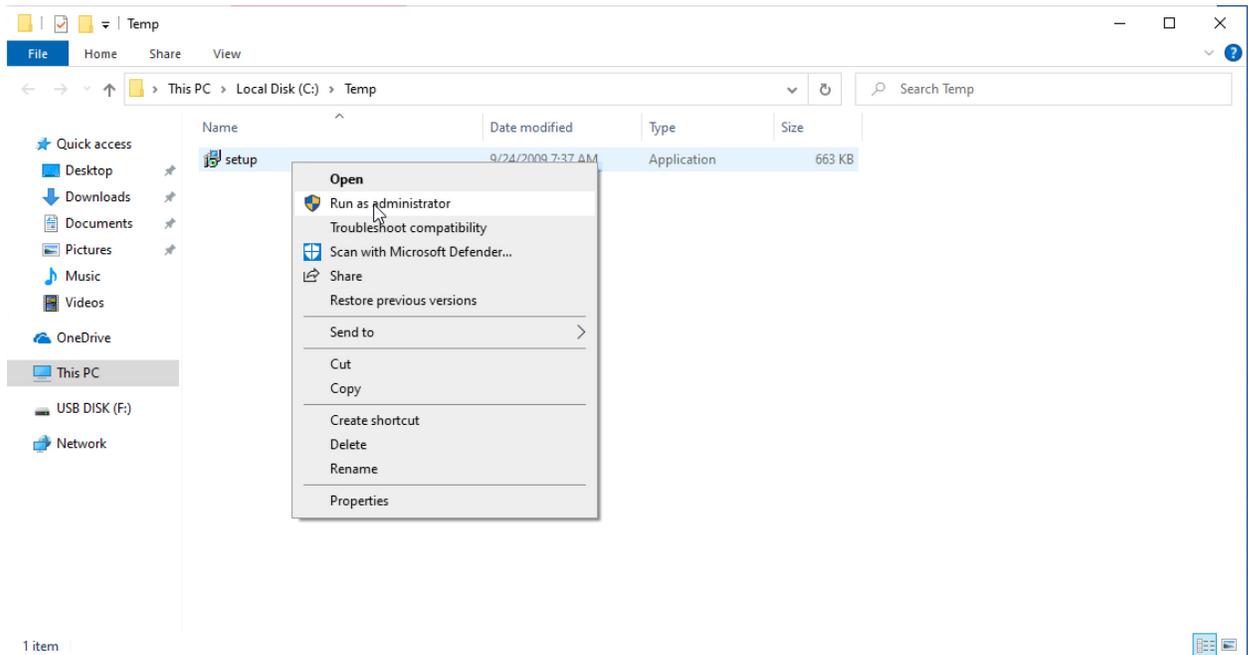


5. Acknowledge the User Account Control permission (enter the administrator account password if prompted).

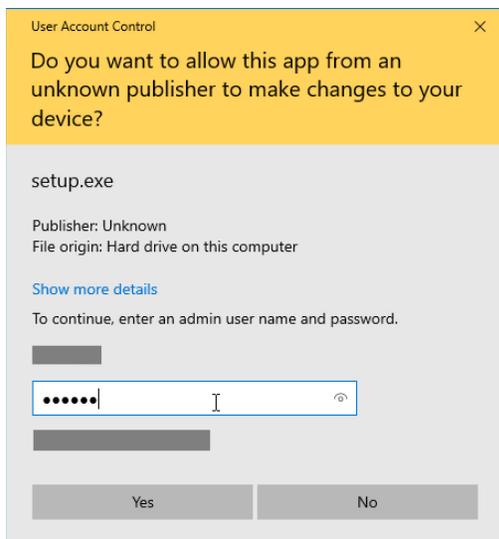


- Click “Next” when the “Microbact 2009 Setup” window opens.

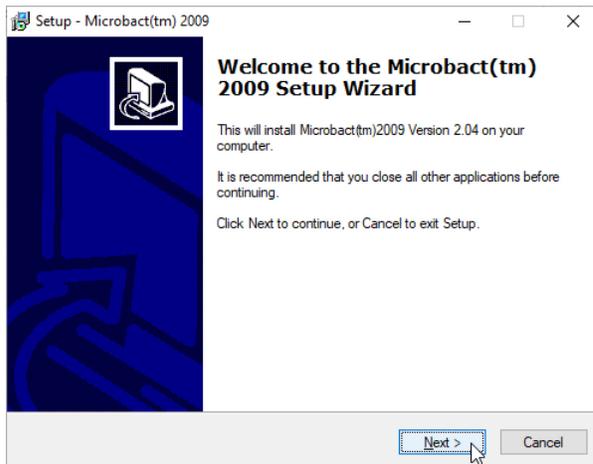


- Accept the agreement and click “Next”.

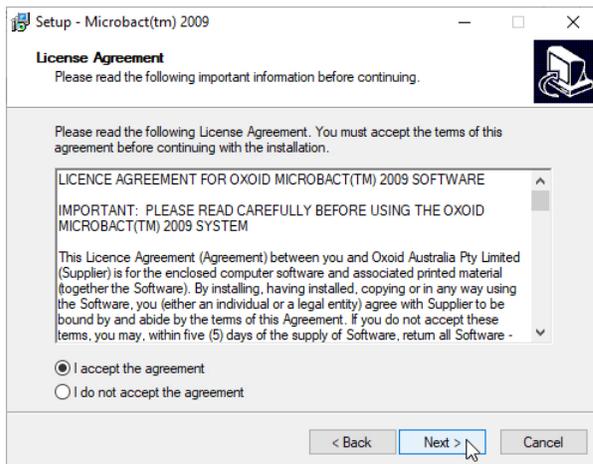


- Choose a destination folder and click “Next”.

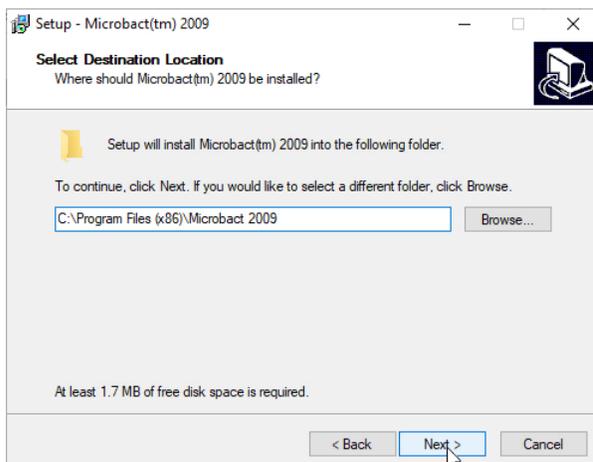


9. Select the Start Menu folder and click “Next”.

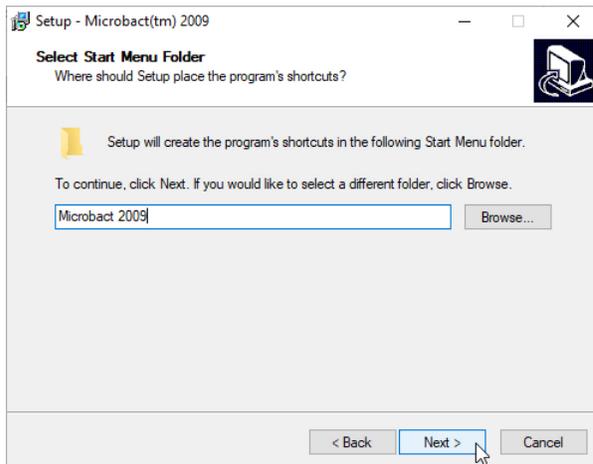


10. Agree to create a desktop icon if desired. Click “Next”.

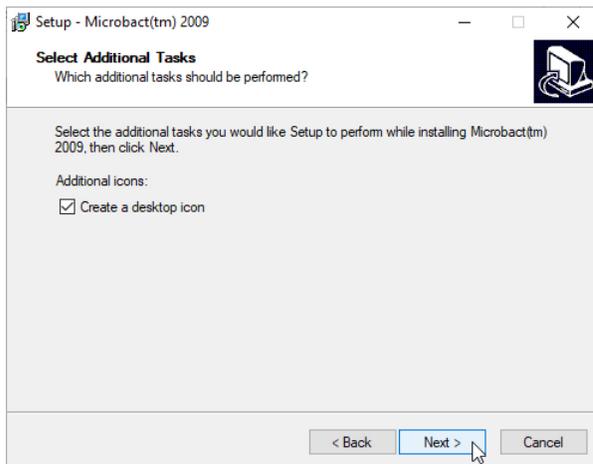

11. Review installation settings and click “Install” when ready.

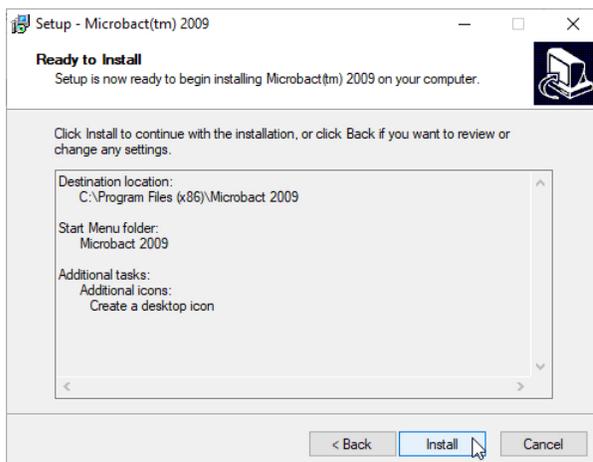

12. When setup is complete, clear the “Launch Microbact 2009” checkbox and click “Finish”.

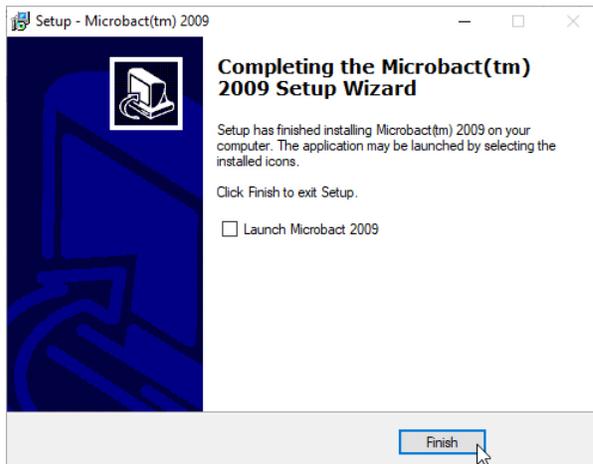


Enabling Use for Windows Users Without Administrator Privileges

If the software will be run from an account without administrator privileges, an administrator of the PC must ensure that user accounts are allowed write access to the main Microbact folder. If this is not done, users may see a “File access denied” and/or a “Cannot create file” message. See examples below:




Allow Write Access to Microbact 2009 Folder


1. Right-click the “Microbact 2009” folder (by default located under C:\Program Files (x86)) and select “Properties”.

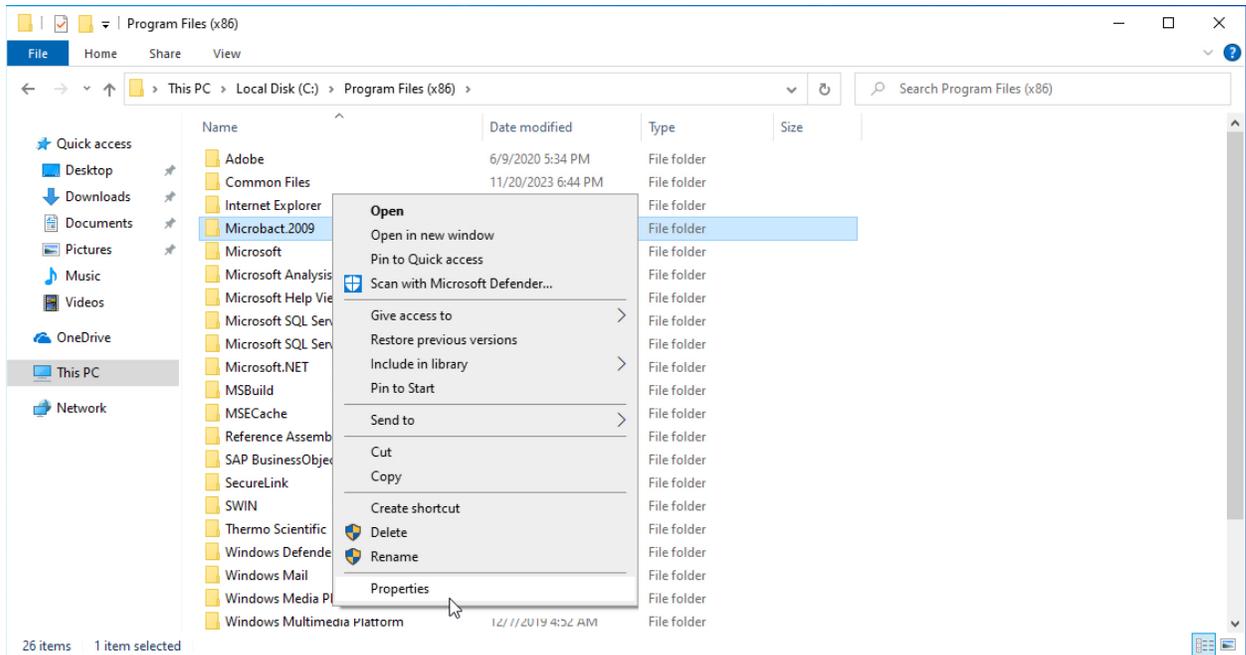


2. Click on the “Security” tab, locate the “Users” group for the PC, and click “Edit”.

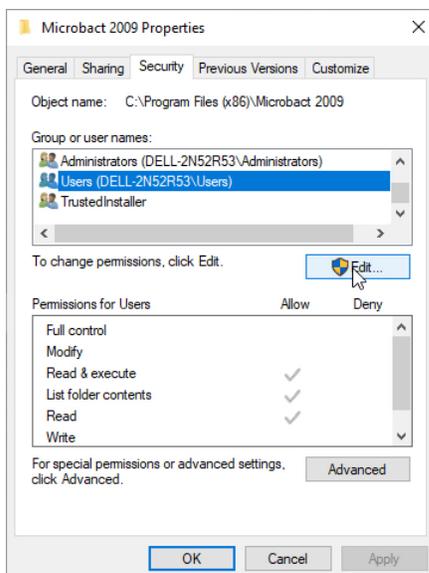

3. Check the box to Allow Write permissions for Users, then click “OK”.

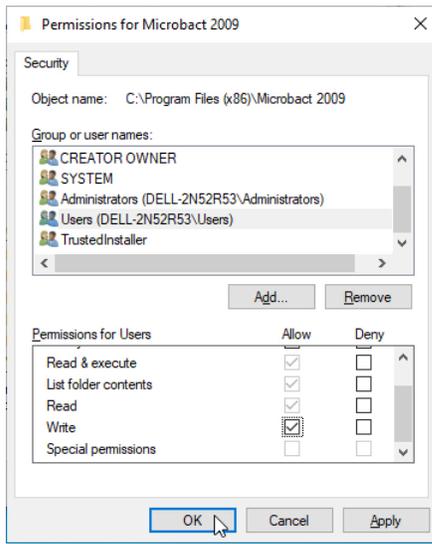

4. Confirm any changes by clicking “OK”.

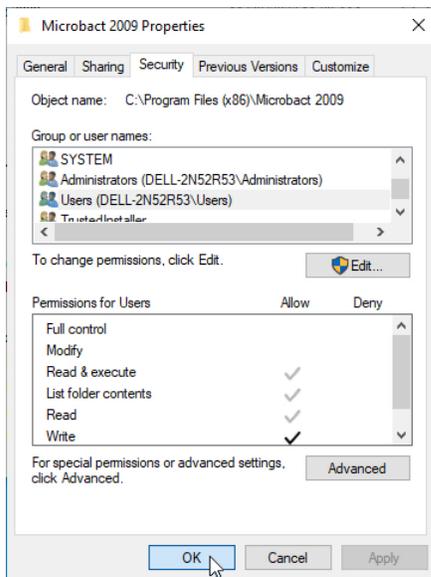

5. It should now be possible for standard users to run Microbact successfully.

For more information and for installation support, please contact your technical support team. Contact information for technical support is provided on the next page.

Technical Service Contact Information

Country	Email Address	Phone Number
Austria	microbiology.techsupport.de@thermofisher.com	+49 (0)281 152 266
Belgium	microbiology.techsupport.benelux@thermofisher.com	+31 629 550 379
Czech Republic	mikrobiologie.tech.podpora.cz@thermofisher.com	+42 0800 101 038
Denmark	microbiology.techsupport.nordic@thermofisher.com	+45 449 797 35
Finland	mikrobiologia.tekninentuki.fi@thermofisher.com	+358 (0)10 329 2496
France	microbiologie.techsupport.fr@thermofisher.com	+33 (0)472 523 372
Germany	microbiology.techsupport.de@thermofisher.com	+49 (0)281 152 266
Italy	microbiologia.supportotecnico.it@thermofisher.com	+39 (0)295 059 350
Netherlands	microbiology.techsupport.Benelux@thermofisher.com	+31 629 550 379
Norway	microbiology.techsupport.nordic@thermofisher.com	+47 230 396 90
Portugal	microbiologia.apoiocliente.pt@thermofisher.com	+35 121 423 5355
Slovakia	mikrobiologie.tech.podpora.cz@thermofisher.com	+42 0800 101 930
Spain	microbiologia.soporte.es@thermofisher.com	+34 91 382 20 23
Sweden	microbiology.techsupport.nordic@thermofisher.com	+46 (0)86 266 050
Switzerland	microbiology.techsupport.de@thermofisher.com	+49 (0)281 152 266
United Kingdom	microbiology.techsupport.uk@thermofisher.com	+44 (0)1256 694 238
United States	USCLEtechsupport@thermofisher.com	1-800-642-7029 (option 2)
European Instrument Services	microbiology.instrumentservice.europe@thermofisher.com	-